



# **TF/TFC & SPS**

## **Laboratory Dispensers and Solution Makers**

### **G2 Operation Manual**

Rev 2.0  
17<sup>th</sup> June 2011

## TF G2 Operation Manual

Edition: XXXXXX 2009 (1st edition)  
Related to TFC v 2.0.0.1b2 version of software

Document number: TFCG2OPMN02\_EN.DOC

© Copyright 2009 Datacolor. All Rights Reserved. This material may not be reproduced or copied, in whole or in part, without the written permission of Datacolor.

Whilst every effort has been made to ensure the accuracy of this manual, Datacolor can accept no responsibility for any omissions or errors it may contain. Datacolor reserves the right to make improvements and/or changes in the product(s) and/or program(s) described in this manual at any time.

MS-DOS™ and Windows 98/ME/2000/XP/Vista/7 and SQL Server are the trademarks of the Microsoft Corporation.

### PC specifications:

- 2 GHz Intel Quad Core 2 32-bit (x86) or 64-bit (x64) processor.
- 2 GB of system memory.
- 256 MB of graphics memory.
- Support for DirectX 9 graphics.
- 320GB of hard drive capacity with 15 GB free space.
- DVD-ROM Drive.
- Audio output capability.
- Internet access capability.
- 1 serial port
- 2 USB Ports
- Monitor - 17" - 19" LCD
- Windows XP Professional ( 32 or 64 bits) SP3 or above
- Windows 7 ( 32 or 64 bits) SP1 or above

### Corporate Headquarters - United States

Datacolor Corporate Headquarters  
5 Princess Road  
Lawrenceville, NJ 0864

## CONTENT

<b>DOCUMENT VERSION UPDATES .....</b>	<b>6</b>
<b>ABOUT THIS MANUAL .....</b>	<b>8</b>
<b>WHO SHOULD READ THIS MANUAL? .....</b>	<b>9</b>
<b>HOW TO USE THIS MANUAL.....</b>	<b>9</b>
<b>CHAPTER 1: GETTING STARTED .....</b>	<b>10</b>
<b>1-1 About your TFC system.....</b>	<b>11</b>
<b>1-2 System overview .....</b>	<b>11</b>
<b>1-3 Features and benefits .....</b>	<b>13</b>
<b>1-4 Starting up the System .....</b>	<b>15</b>
<b>1-5 Shutting Down the System.....</b>	<b>16</b>
<b>1-6 Using the ALAB Folder .....</b>	<b>17</b>
<b>CHAPTER 2: TF DP PROGRAM OVERVIEW .....</b>	<b>21</b>
<b>2-1 TF DP Program – Configuration.....</b>	<b>24</b>
2-1-1 TF DP PROGRAM – CONFIGURATION - MACHINE .....	25
2-1-2 TF DP PROGRAM – CONFIGURATION – DATABASE SETUP .....	27
2-1-3 TF DP PROGRAM – CONFIGURATION – USER ACCOUNT .....	28
2-1-4 TF DP PROGRAM – CONFIGURATION – AUTHORITY .....	29
2-1-5 TF DP PROGRAM – CONFIGURATION – SELECT LANGUAGE .....	30
<b>2-2 TF DP Program – Parameters.....</b>	<b>31</b>
2-2-1 TF DP PROGRAM – PARAMETERS - PERSONAL .....	32
2-2-2 TF DP PROGRAM – PARAMETERS - DILUTION.....	34
2-2-3 TF DP PROGRAM – PARAMETERS – VALVE .....	35
2-2-4 TF DP PROGRAM – PARAMETERS – PASTE .....	36
2-2-5 TF DP PROGRAM – PARAMETERS – IMPORT PARAMETER.....	37
2-2-6 TF DP PROGRAM – PARAMETERS – EXPORT PARAMETER.....	41
2-2-7 TF DP PROGRAM – PARAMETERS – BATCH/RECIPE PARAMETER.....	44
2-2-8 TF DP PROGRAM – PARAMETERS – PRINT PARAMETER .....	46
<b>2-3 TF DP Program – Utility .....</b>	<b>48</b>
2-3-1 TF DP PROGRAM – UTILITY – BACKUP.....	49
2-3-2 TF DP PROGRAM – UTILITY – RESTORE .....	50
2-3-3 TF DP PROGRAM – UTILITY – INJECTOR ANALYSIS .....	51
2-3-4 TF DP PROGRAM – UTILITY – RECIPE ANALYSIS.....	52
<b>2-4 TF DP Program – Database .....</b>	<b>53</b>
2-4-1 TF DP PROGRAM – DATABASE – PRODUCT .....	54

2-4-2 TF DP PROGRAM – DATABASE – PROCESS .....	56
2-4-3 TF DP PROGRAM – DATABASE – SOLUTION .....	58
2-4-4 TF DP PROGRAM – DATABASE – DYEPOD TRAY.....	61
2-5 TF DP PROGRAM – BATCH/RECIPE MANAGEMENT .....	62
2-5-1 TF DP PROGRAM – RECIPE MANAGEMENT – RECIPE HEADER.....	63
2-5-2 TF DP PROGRAM – RECIPE MANAGEMENT – RECIPE DETAIL .....	66
2-5-4 TF DP PROGRAM – BATCH MANAGEMENT .....	70
<b>2-6 TF DP Program – Solution Stock .....</b>	<b>78</b>
<b>2-7 TF DP Program – History Report .....</b>	<b>81</b>
<b>2-8 TF DP Program – File.....</b>	<b>83</b>
<b>CHAPTER 3: TF/TFCCTRL PROGRAM .....</b>	<b>84</b>
<b>3-1 TF/TFC Ctrl program – Overview .....</b>	<b>84</b>
TF/TFC CTRL PROGRAM – OVERVIEW – MANU BAR .....	85
TF/TFC CTRL PROGRAM – OVERVIEW – STATUS BAR.....	87
<b>3-2 TF/TFC Ctrl program – Dispense .....</b>	<b>88</b>
<b>3-3 TF/TFC Ctrl program – Machine.....</b>	<b>90</b>
3-3-1 TF/TFC CTRL PROGRAM – DIAGNOSTIC .....	90
3-3-2 TF/TFC CTRL PROGRAM – SPECIFIC WT. CALIBRATION .....	110
3-3-3 TF/TFC CTRL PROGRAM – CYLINDER CALIBRATION .....	111
<b>3-4 TF/TFC Ctrl program – Parameter.....</b>	<b>114</b>
3-4-1 TF/TFC CTRL PROGRAM – PARAMETER – SETUP.....	114
3-4-2 TF/TFC CTRL PROGRAM – PARAMETER – AUX SETUP .....	121
3-4-3 TF/TFC CTRL PROGRAM – PARAMETER – PASTE SETUP.....	125
3-4-4 TF/TFC CTRL PROGRAM – PARAMETER – PRINT SETUP .....	127
3-4-5 TF/TFC CTRL PROGRAM – PARAMETER – AGITATION SETUP.....	129
<b>3-5 Summary of features of TFC-P system .....</b>	<b>130</b>
<b>PART 4: AUTOLAB SPS PROGRAM .....</b>	<b>132</b>
<b>4-1 Autolab SPS Program – Overview .....</b>	<b>132</b>
<b>4-2 Autolab SPS Program – Dispense .....</b>	<b>135</b>
4-2-1 AUTOLAB SPS PROGRAM – DISPENSE – MAKING SOLUTION .....	135
4-2-2 AUTOLAB SPS PROGRAM – DISPENSE – MANUAL DISPENSE WATER .....	137
4-2-3 AUTOLAB SPS PROGRAM – DISPENSE – SOLUTION LIST.....	138
<b>4-3 Autolab SPS Program – Database .....</b>	<b>139</b>
4-3-1 AUTOLAB SPS PROGRAM – DATABASE – SOLUTION DATA.....	139
4-3-2 AUTOLAB SPS PROGRAM – DATABASE – PROCESS DATA .....	141

4-3-3 AUTOLAB SPS PROGRAM – DATABASE – HISTORY DATA.....	142
<b>4-4 Autolab SPS Program – Parameter Setup.....</b>	<b>143</b>
<b>4-5 Autolab SPS Program – Machine.....</b>	<b>148</b>
4-5-1 AUTOLAB SPS PROGRAM – MACHINE – DIAGNOSTIC.....	148
4-5-2 AUTOLAB SPS PROGRAM – MACHINE – VALVE CALIBRATION .....	155

## Document version updates

---

Revision 2 has the following updates made:

- 1) Administrative rights - Stock solution can be updated even no such rights given by supervisor (bug) fixed on 2<sup>nd</sup> march
- 2) Dilution of multiple bottles – not updating date and time ( bug - corrected)
- 3) Recipe calculation unit – (bug - corrected)
- 4) The restore option is password protected – fixed on 2<sup>nd</sup> march  
if customer is chief group user or engineer then can skip and will show another passwords form just key in our engineer password “alabt” or chief user's password or engineer's password set by customer and restore program will start.
- 5) The water volume should appear on print or print preview in DP program Batch /recipe menu. – Feature development - fixed on 2<sup>nd</sup> march
- 6) The Control was updated for V5 injector - updated 2<sup>nd</sup> March 2011
- 7) SPS was updated for some default value - updated 2<sup>nd</sup> March 2011
- 8) UDR was update for fieldname of table of database , because last time i changed the some field name , if user print the product , it will get an error messages, and can't select the new field name on Edit Mode - updated 2<sup>nd</sup> March 2011
- 9) Importing file from ITMprocess/dcprocess - now it is possible to change the file name for import. Updated 2<sup>nd</sup> march 2011.
- 10) The restore path should be same as backup. We cannot restore the files located elsewhere than backup path and we have to copy the files to backup directory only.
- 11) Dilution amount in TF (request from Portugal) – dilution amount is changeable but there is no indication how to do it. Now we have added new button “Change amount” which will help user to change the amount as per his requirement ( V2.0.2 May 2011)
- 12) New Import and Export option under Product Menu
- 13) Batch /Recipe Management screen – modified the Batch / Recipe management screen to make it more simple. Modifications are as follows : ( V2.0.2 May 2011)
  - a) “Reset Position” won't be there always , it will appear only if the product input rows are more and out of screen window
  - b) Chemical recipe – There was no indication on how to add “chemical recipe” in existing dye recipe, now there is option to ‘Insert” or user can insert by double clicking on selected chemical recipe from the list.
  - c) New “Insert” and “Delete” buttons for recipe product add windows. – Two buttons are added for user to easily add or delete row as per old G1 software
  - d) “Import Now” button – Removed “Import Now” button on main screen and incorporated it in “Transfer Log” tab. In Transfer log tab, you will get “ “ ( like refresh on Input tab) and also with caption of “Import”.
  - e) New option for Dyepot Capacity – as per request from European and American customers (Rudolph/ Tam), there is option to select the Dyepot Capacity with each dyepot Tray. This is again optional and will active only if selected. In Dye Pot Tray menu – while you define dye pot

tray, you will get this option and system will not save recipe if the total volume is exceed the dye pot capacity.

### Control Program

- 1) Out Conveyor timing to be manually configurable – for customer – XinShen in china
- 2) V5 injector setup is out of engineer setting and now there is graphic option for customer /engineer to understand how to put the V5 injectors in system.

**Addendum:** Important safety note before operating or using the Datacolor Autolab Dispensing System: In light of the prevalent use of smartphones and smart devices in the daily operations of our customers, it is crucial to underscore the significance of adhering to safety protocols, ensuring ongoing training, and regularly updating documentation related to product usage. It is strongly recommended that all personnel refrain from carrying smartphones, smart devices, or magnetic items in close proximity to the Autolab Dispensing System. In rare instances, we have observed that these devices may cause interference, potentially compromising the dispenser's intended operation and posing a safety risk.



## About this Manual

---

This is the *TF Laboratory Dispenser Operations Manual* covering the following TF systems:

- SPS solution maker
- TF dispenser
- TFC dispenser with conveyor
- TFC-P colorant/paste dispenser



## WHO SHOULD READ THIS MANUAL?

---

This Manual is designed for users of TF Laboratory dispenser and solution maker systems, who need to know how to begin using the system. Once you are familiar with how the system works, this Manual will provide a reference to help you carry out specific tasks using the system.

## HOW TO USE THIS MANUAL

---

This manual is divided into four main parts:

- Part 1, *Getting Started*, gives you the basic information you need to start using the system: to get the best out of your system, we recommend that you read **all** of this section before you start
- Part 2, How to use TF DP program
- Part 3, How to use TF Control program
- Part 4, How to use SPS program

### To find the information you are looking for:

Consult the Contents List at the beginning of the Manual.

## CHAPTER 1: GETTING STARTED

---

Part 1, *Getting Started*, is essential reading for anyone using a TF system for the first time. To get the best out of your system, it gives you the basic information you need, without going into detail about the individual options. It is recommended that you read all of this section before you start using the system. A summary of each chapter is given below:

- **About your TFC system**  
The aim of this chapter is to familiarize you with your TF system.
  - Gives an overview of the TF system
  - Describes the key features and benefits of the system
- **Starting up the system**  
this chapter describes how to start your TF system, and what you see on the screen while the system is starting up.
- **Shutting down the system**  
This chapter describes how to shut down your TF system.
- **Using the TF file Folder**  
This chapter explains the main features of the TF file folder, including how to use the menus and screens, and housekeeping tasks such as browsing and deleting information.

## 1-1 About your TF system

---

The aim of this chapter is to familiarize you with your TF system. This chapter:

- Gives an overview of the TF system
- Describes the key features and benefits of the system

## 1-2 System overview

---

The TF/TFC Laboratory Dispenser is an automatic dyestuff solution dispenser, designed to produce accurate and repeatable recipes in an easy and efficient manner. The system is available in different sizes and configurations. All of the TF Laboratory Dispensers include:

- Robot works as transport and dispensing media.
- Stock solution area, where solution bottles are there.
- Bottles for the various dyestuff solutions used
- Bottles for the auxiliary solutions used (Only systems with conveyor, TFC)
- Individual injector for each solution bottle for transferring solution to dye pot.
- A system for automatically refilling the auxiliary bottles (Only systems with conveyor)
- Pressurised paste tank, for storing paste to be dispensed. (TF-P systems only)
- Conveyor system, which moves dye pots into dispensing area (Only systems with conveyor, TFC)
- A manual operating panel
- An internal dispenser control unit
- A dispensing area, where the recipes are dispensed
- A computer for controlling the system

The solutions are transferred to dispensing area, and dispensed into recipe pots, via individual injector, by using robot. Dyestuffs are dispensed gravimetrically (the injectors close when a specified weight of solution has been weighed into the beaker). Water and auxiliaries are usually dispensed volumetrically, based on time interval (valves close when a certain amount of solution has passed thorough the valve). However water and auxiliaries can be dispensed gravimetrically, if required. Paste is dispensed gravimetrically (the paste dispensing valve is closed, when a specified weight of solution has been weighed into the beaker).

### Dyestuff store

The number of bottles on your system will depend on the size and model of your system. Magnetic stirrers under each bottle prevent sedimentation.

Each bottle has individual injector.

### Auxiliary store

There are 7 auxiliary and 1 water bottles. A water filter is used to ensure the

water supply is free from impurities.

### Auxiliary refill

There are 7 auxiliary refill bottles.

### Paste store

The number of paste dispensing modules on your system will depend on the order of your TF-P system, up to a maximum of 3 paste tanks is allowed. It is also possible to add additional paste modules after the TF-P system has been installed, but cannot exceed 3 modules in total.

## Operating panels

The AGITATION speed control regulator is used to control the stirring devices under the solution bottles. There are 4 regulators maximum, for every 40 bottles has one regulator. The POWER button switches the power on and off. When this switch is in the "0" position, all electricity will be turned off **except** night-time agitation. To turn off all power, use the main switch on the side of the system.

As a safety feature, full power is automatically turned off if you open the electric cabinet door of the system. This feature is called "interlock". It is possible to disable the interlock by an engineer.

**WARNING: If the INTERLOCK DISABLE button is lit, the power will stay on when you open the side doors.**

The large red emergency stop button "**EMG. STOP**" will automatically cut all power to the system. To restart the system if the emergency stop has been activated, the button should be twisted to disable the cut off of power.

## Main control cabinet

The control cabinet at the front of the system consists of:

- LA50B CPU Module
- LA50B SERVO module
- LA50B D.I.O. Module

This is for controlling the movement of robot and dispensing.

## 1-3 Features and benefits

---

### Accuracy

Using the gravimetric dispensing method, an accuracy of  $\pm 0.02\text{g}$  is achievable, enhancing reproducibility in dyeing in the laboratory and, consequently, in production.

### Speed

A batch of 12 recipes can be dispensed in about 12-20 minutes (assuming each recipe has a volume of 180 ml and contains three dyestuffs and one to three auxiliaries).

### Multi-tasking

The TF computer software is multi-tasking — while the system is dispensing recipes, you can continue using the computer, e.g. input new recipes ready for dispensing when the first set is completed.

### Monitoring stock solutions

The system can detect and warn you when a solution is at a low level. You can also specify how long a solution will remain usable (i.e. specify its shelf life). The system will warn you when a solution's lifetime has expired.

### Anti-sedimentation

A magnetic stirring device prevents sedimentation: you can adjust the speed of stirring.

### Measurement units

You can choose grams per litre, or percentage of the weight of fabric or weight of solution, or a mixture of units

### Automatic selection of the best concentration

If you have a dye on the system in several different concentrations, the system can calculate and automatically dispense the optimum concentration for a recipe.

### Automatic water addition

You can set up the system so that water is automatically added to a recipe in order to achieve the correct total volumes and liquor ratios.

## Printing results and reports

The results of all the recipes you dispense are automatically printed. Also other reports, such as a list of the recipes or solutions specified on the system, can be printed.

## Alarms

The system can detect if a dye pot tray is not placed on the scale correctly, and will sound an alarm and give a warning message on the screen.

In the same way, the system indicates when dispensing has finished, and if all the safety doors are closed properly.

## Operational flexibility

The system can allow for products to be added manually after the recipe has been dispensed. You can create groups of dyes and auxiliaries to simplify recipe input.

## Linkage

Linking with the “DCI Match” is quick and error free – manual keying-in is an error prone operation.

## 1-4 Starting up the System

---

This chapter describes how to start your TF system.

⇒ **To start your TF system:**

1. Turn on the main power switch on the side of the system
2. Ensure the emergency stop button is not depressed.
3. Push the dispenser ON/OFF switch to the ON position.
4. Turn the AGITATION speed tuner on the dispenser to the correct speed.
5. Turn ON the printer; check that the printer is on-line and paper is loaded. If you require more information consult the manual supplied with your printer.
6. Turn ON the computer. (The ON/OFF switch is usually on the front of your computer).
7. The TFC software is loaded automatically. What you see when the software has loaded depends on how your system has been set up, and which system you are running.

**If you see a “*PLC Module Not Connected*” message, this indicates that the dispenser is turned off. Please turn the dispenser ON.**

## 1-5 Shutting Down the System

---

This chapter describes how to shut down your system. If you wish to shut down the system, e.g. at weekends or holidays, follow the procedure below.

⇒ **To shut down your system:**

1. Turn the printer OFF. Check your printer manual if you are not sure where the ON/OFF switch is.
2. Quit all of the application programs that are running.
3. Use Windows standard procedure to shut down the computer.
4. Turn OFF the power to the computer.
6. Turn the dispenser ON/OFF switch to the OFF position.
7. If night time agitation is not needed, you can also switch off the power to the dispenser.

It is suggested that if the system is to be left unused for a long period of time, all solutions are removed from the system and all tubes are cleaned with water.



## 1-6 Using the ALAB Folder

---

This chapter describes how to shut do

This chapter explains:

- Running programs from the ALAB File folder and menu
- How navigate the menus and select functions
- Using the Toolbar buttons
- Using the Buttons
- Using the Message Box
- Special keys you can use

### ALAB File Folder and Menu

The software can be accessed in a number of different ways. Firstly, all programs will be automatically started up when turning on the system.

Additional programs can be accessed from the ALAB file Folder, which can be found on the main Window desktop. Clicking on the ALAB folder will open the ALAB Window, which will then display all the available programs, as icons. Double clicking on an icon will execute that program.

Programs can also be run from the ALAB Menu. This can be accessed by selecting the Start button on the Taskbar on the bottom left hand side of the screen. Then select Programs and then ALAB from the Menu options shown. Within the ALAB Menu will be displayed all the available programs, and selecting an option will execute that program.

The software is split into 3 main programs:

1. TF or TFC DP for the inputting of solution, dyestuff and recipes information, and management of the dispense batch.
2. TF or TFC dispense control program, for dispensing of recipes and maintenance of the dispenser.
3. All 2 programs are run at the same time, and you can swap between programs at any time. This is called Multi-Tasking. To move between programs, you have to select the required program from the Taskbar. This is not normally visible, and will only be displayed once the mouse pointer is moved to the bottom of the screen. Once the Taskbar is displayed, the different programs that are presently running will be displayed as buttons on the Taskbar. To move from one program to another, just click on the button required. Alternatively, pressing ALT+⇐⇒ (the TAB button) will swap from one program to another, in a cycle.

## Using the menu function

A menu is a list of alternative options from which to choose from. Each option allows you to carry out a particular task. Within a main menu option, there may be additional sub-options available, and these will be shown as a sub-menu underneath. When an option or sub-option is shown in light grey colour instead of the normal dark gray colour, it means that you cannot access that function.

⇒ **To select an option from a menu:**

1. Use the mouse to click on a menu option on the top main menu bar, then select the function required, using one of the following methods:
  - Using cursor keys only:
    - Use the ← and → cursor keys to move from left to right
    - Use the ↑ and ↓ cursor key to move up and down a sub menu
    - Press ENTER key to select the option highlighted.
  - Using a combination of cursor keys and highlighted letter keys:
    - Use the ← and → cursor keys to move from left to right
    - Type the key of the letter underlined within the option required.
  - Using the mouse only:
    - Move the mouse left and right and up and down the menus and sub-menus, then click on the function required, using the right mouse button.

## Using the Toolbar

Toolbar allows you to quickly access regularly used functions.

## Using the Message box

When using the software, the system will sometimes prompt you with the message box. There are two kinds of message box, which are as follows:

- Warning message box:
  - This is a warning message only. Use the mouse to acknowledge the condition and continue, by clicking on the **'OK'** button displayed, or by pressing **ENTER**.
- Confirm message box:
  - When a Confirm message box is displayed, a description of the event which has occurred, or which will occur, will be shown, along with two action buttons, normally **'OK'** or **'CANCEL'**, or **'YES'** and **'NO'**. Select **'OK'** or **'YES'** to accept the action, **'CANCEL'** or **'NO'** to stop the action and return to the previous step.

## Typing in information

When inputting information, for example adding a recipe to the system, you use a special screen called a form screen, with boxes or fields that need to be filled in with information.

⇒ **To fill in form screens:**

- When inputting information into a form screen with more than one field input box, the position where data is to be presently inputted will be denoted by a flashing or coloured cursor. New information can be directly typed in. Once completed presses **ENTER** to accept the input and move to the next field.
- If data already exists in the field selected, then this will be highlighted in blue. To delete the existing data from the field, and add new data, start typing in the new information directly, and this will automatically clear the existing data. If the existing data is to be modified, then use the cursor keys to move to the position in the field where the modification is to be made, and type in the new information. Use the **DELETE** key to remove data not required. Once completed, presses **ENTER** to accept the input.
- The mouse can also be used to move from field to field, and clicking within a box will allow modification to that field.
- The **TAB** key, ⇔, can also be used to move from field to field.

## Special keys

Certain keys on your keyboard have special actions:

Use this key:	To:
ESC	To close the current screen
F1	Insert a line in Recipe Add/Modify form
F12	Delete a line in Recipe Add/Modify form
BACKSPACE	Delete one character to the left of the cursor.
← (Cursor key)	Move the cursor left one character.
→ (Cursor key)	Move the cursor right one character.

## The Number keypad

The block of keys on the far right of your keyboard is the Number keypad. You can use this keypad in two ways:

- If the **NUM LOCK** light is lit, you can use the pad to type in numbers
- If the **NUM LOCK** light is not lit, you can use the cursor keys.

To change between the two modes, press the **NUM LOCK** key.


## CHAPTER 2: TF DP program overview

---

This program is designed to allow input, modification and set up of all information relative to the products, processes, solutions, recipes and batches used in the laboratory and on the dispenser.


Before starting to enter the data into the computer, you need to collect all of the information regarding the products used. e.g., how the different dyestuff solutions should be made up, what are the different dyestuffs that are used, what different concentrations of solutions are required, which solutions will be dispensed on the system, and so on.

There is an icon , **ALTFDp** is for TF, and **ALTFCDP** is for TF systems, on the desktop of computer. Double click the icon to launch DP program, and you will see the screen like below:

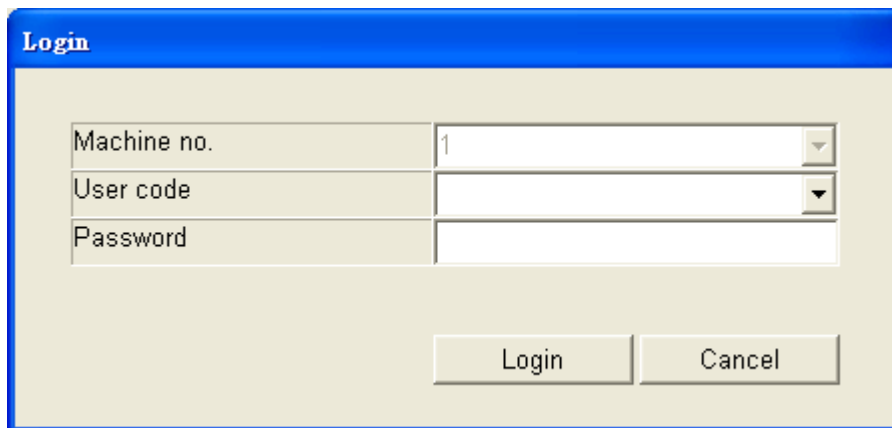


The image shows a 'Login' dialog box with a blue title bar. It contains three input fields: 'Machine no.' with the value '1', 'User code', and 'Password'. Each field has a dropdown arrow on its right side. At the bottom right, there are two buttons: 'Login' and 'Cancel'.

Firstly, engineers have to give the dispenser a machine number to let DP software create the databases. Machine number is a very important configuration value for identifying different machines. Different machines have different machine numbers.

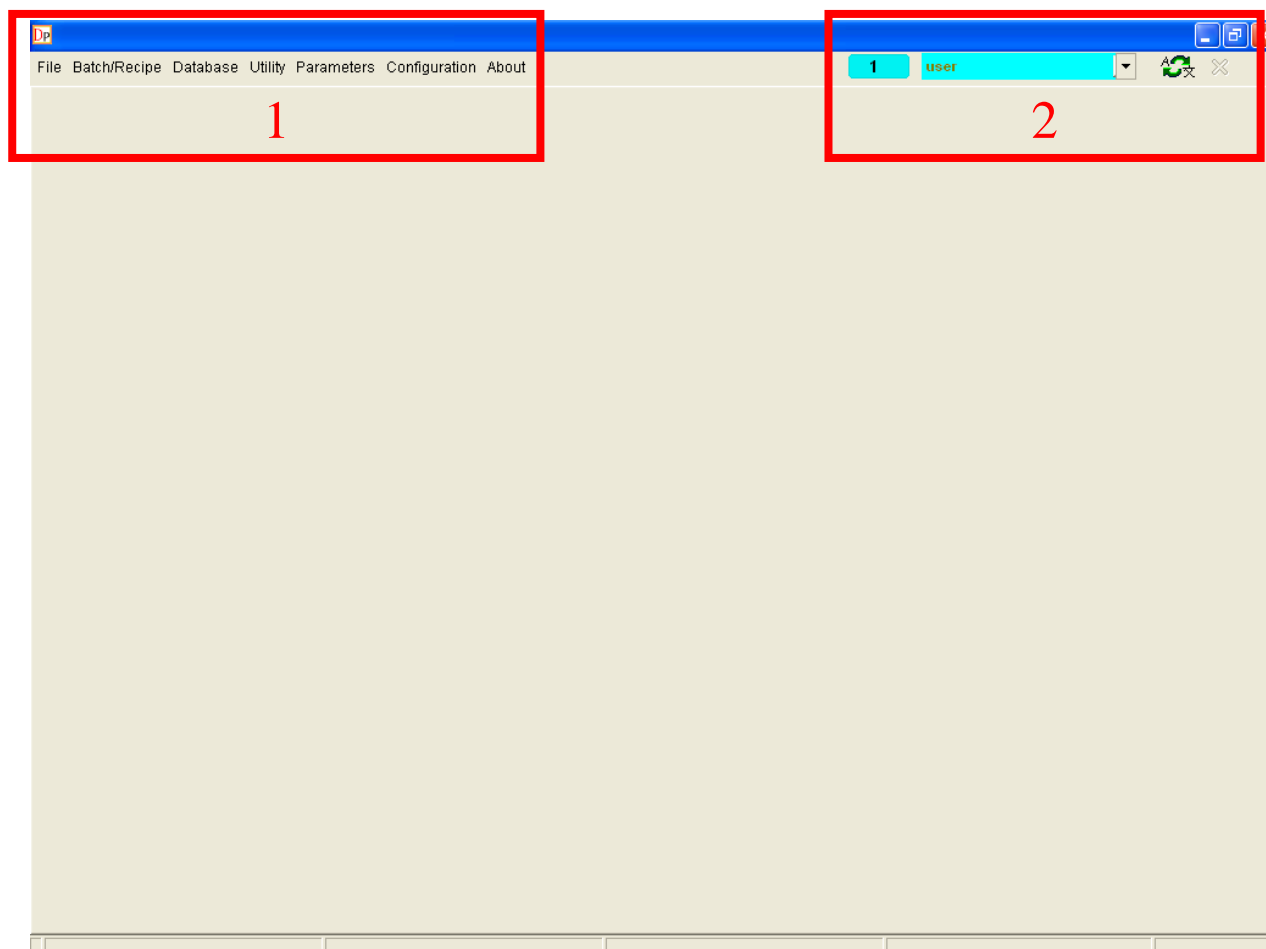


# 1. Menu Bar :

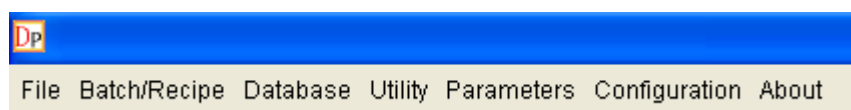

# 2. Tool Bars




<p>1</p>	<p>The machine number of the dispenser which physically link with the computer.</p>
<p>Remote 1</p>	<p>The machine number of the database current logging in. If the machine logging in is not for the actual dispenser, it will show Remote in front of machine number.</p>
<p>Employee ID 001</p>	<p>The user who is using the software. It is allows to quick switch users who are in the same authority group by selecting in the drop down list.</p>
<p>A 文</p>	<p>Select Language, start the language selection form</p>

	Close the opened function tab.
---	--------------------------------

## 2-1 TF DP Program – Configuration

---

There are 5 items in the Configuration, and they are password protected. Permission of changing the configurations is depending on the authority of login user. They are divided into 5 sections as follows:

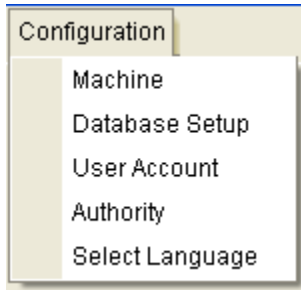

- **Machine**
- **Database Setup**
- **User account**
- **Authority**
- **Select Language**



## 2-1-1 TF DP Program – Configuration - Machine

The screen below is “**Machine Configuration**”

Machine model	LAG01-160
Physical dispense No.	1
Enable sample weight scale	No
Enable paste function	No
Global injector dispense accuracy (g)	0.02
Global chem/water valve accuracy (g)	0.10
Global paste accuracy (g)	1.00
Batch dispensing order	Injector dispensing first

- **Machine Model**

Show the machine model of the database which is logging in.

- **Physical Dispenser No.**

Show the machine number which is physically connected with the computer.

- **Enable sample weight scale**

If you have fabric scale is connect, select “Yes”, the options of weighing and zeroing fabric scale in “Recipe Add/Modify” form will be enabled automatically. And the Parameters for fabric scale will be also enabled in the “Fabric Weigh Scale”.

After enable the option, the fabric scale icon  will appear beside the parameter. Click the icon to set the fabric scale parameter.

Balance Data		Weight	
Scale Setting			
Scale type	METTLER		
<input checked="" type="checkbox"/> Define tare command	Z		
Other Scale Type Setting			
Scale start command	SIR		
<input type="checkbox"/> Define command pre code	27		
RS232 Setting			
COM port	2		
Baudrate	9600		
Save		Start	
		Tare	

Suggest setting the correct “COM Port” and Scale Type only, and leaving the other settings as default.

- **Enable Paste function**

If your system is TF-P system, select “Yes” to enable paste function. The reserved bottle position numbers 231 to 233 in “paste valve” in solution database will then become available, which should be used to define paste product data.

- **Global injector dispense accuracy (g)**

Define the global injector dispense accuracy for all injectors. Minimum is 0.02g.

- **Global chemi/water dispense accuracy (g)**

Define the global chemical or water valves dispense accuracy. Minimum is 0.1g.

- **Global paste dispense accuracy (g)**

Define the global paste valves dispense accuracy. Minimum is 1g.

- **Batch dispensing order**

There are 3 options, “injector first”, “Valve first” and “Paste first”.

If choose the option “Injector first” the dispensing order will be Injector-valve-paste.

If choose the option “Valve first”, the dispensing order will be Valve (Gravimetrically)-injector-paste.

If choose the option “Paste first”, the dispensing order will be paste (Gravimetrically)-injector-valve.

## 2-1-2 TF DP Program – Configuration – Database Setup

---

DP software is a centralized database structured program. Database Setup is the setting of accessing databases, which stored in a server computer.

The screen below is “**Database Setup**” setting form.

- **Database server**

This is the server name of the database stored. Server name is combined by 2 parts, “PC name\SQLEXPRESS”. In front of “\” is the computer name and after “\” is a fixed word “SQLEXPRESS”. The default of the server name will be identified by local computer name. Engineers can type it or by select in the drop down list. If the computer is the server, the setting can keep default.

- **Authentication**

There are 2 types of authentication to log in G2 software databases. One is Windows authentication and another one is SQL Server authentication. SQL Server authentication is the default option.

- **Database login name**

This field is only for SQL Server authentication to put SQL Server account.

- **Password**

This field is only for SQL Server authentication to put SQL Server login password.

## 2-1-3 TF DP Program – Configuration – User Account

---

The screen below is “**User Account**” screen.

In G2 software, different users can have their own personal parameters to adapt their work habit. Also it is for supervisor give different authority for the users in different user groups.

User code [dropdown] New

User group [dropdown]

User password [masked]

Total : 2 ☒ Keep previous

User code	User group
TEMP	CHIEF

Save Delete

- **User Code**  
Set user code
- **User Group**  
Set the user group for the user. Different user group has different authority to access functions.
- **User Password**  
Set user password to log in database.

## 2-1-4 TF DP Program – Configuration – Authority

The screen below is “**Authority**” screen.

There are 3 default groups, “User”, “Power User” and “Chief”. The users in different User Groups have different authorities. Depend on the needs, it allows to define “Enable”, “Disable” or “Browse” authority to the functions for the user group.

- **Database**  
It is the authority to use database function.
- **Utility**  
It is the authority to use utility function.
- **Parameter**  
It is the authority to use parameter function.
- **Configuration**  
It is the authority to use configuration function.
- **Function**  
It is the authority to some specific functions.

## 2-1-5 TF DP Program – Configuration – Select Language

---

The screen below is “**Language**” screen.




There are 20 languages for selecting, select the language, and click “OK” to close the screen. And then all wordings showed will be the language you selected.

## 2-2 TF DP Program – Parameters

---

There are many parameters in the Parameter. Permission of changing the parameters is depending on authority of the user group which the login user assigned to.

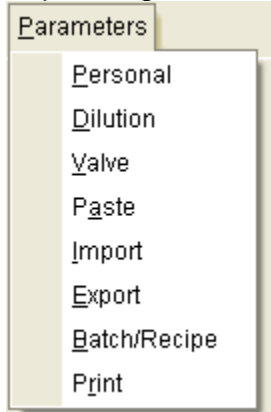

- **Personal**
- **Dilution**
- **Valve**
- **Paste**
- **Import**
- **Export**
- **Batch/Recipe**
- **Print**

## 2-2-1 TF DP Program – Parameters - Personal

Personal parameter is the parameters for the login user only. Users can set the parameters by personal working habits or assigned missions. After quick switch users or new user login, the personal parameters will be loaded for the new login/switch user.

The screen below is “**Personal**” screen.

- User Code**  
 The user code for the current user. It is browse only.
- User Group**  
 The user group which the user assigned to. It is browse only.
- User batch serial no.**  
 The latest batch serial no. of the user. When you set a new user, the serial No. will begin in 001 and be added in sequence. Press “Reset” button will reset the serial No.
- Sample weight**  
 Set the default sample weight for exhaust dyeing recipes. This setting will be automatically applied when creating recipes.
- Liquor ratio**  
 Set the default liquor ratio for exhaust dyeing recipes. This setting will be automatically applied when creating recipes.



- **Total volume**

Set the default total volume for exhaust dyeing recipes. This setting will be automatically applied when creating recipes.

- **Dispense water**


It is the parameter to set if dispensing water into the recipes. If choose "No", the dispenser will not dispense water into the beaker.

## 2-2-2 TF DP Program – Parameters - Dilution

---

Dilution parameter is the parameters for the function “Dilution in TF system”. The software will follow the parameter to filter the stronger solution as dilution source.

The screen below is “**Dilution**” screen.

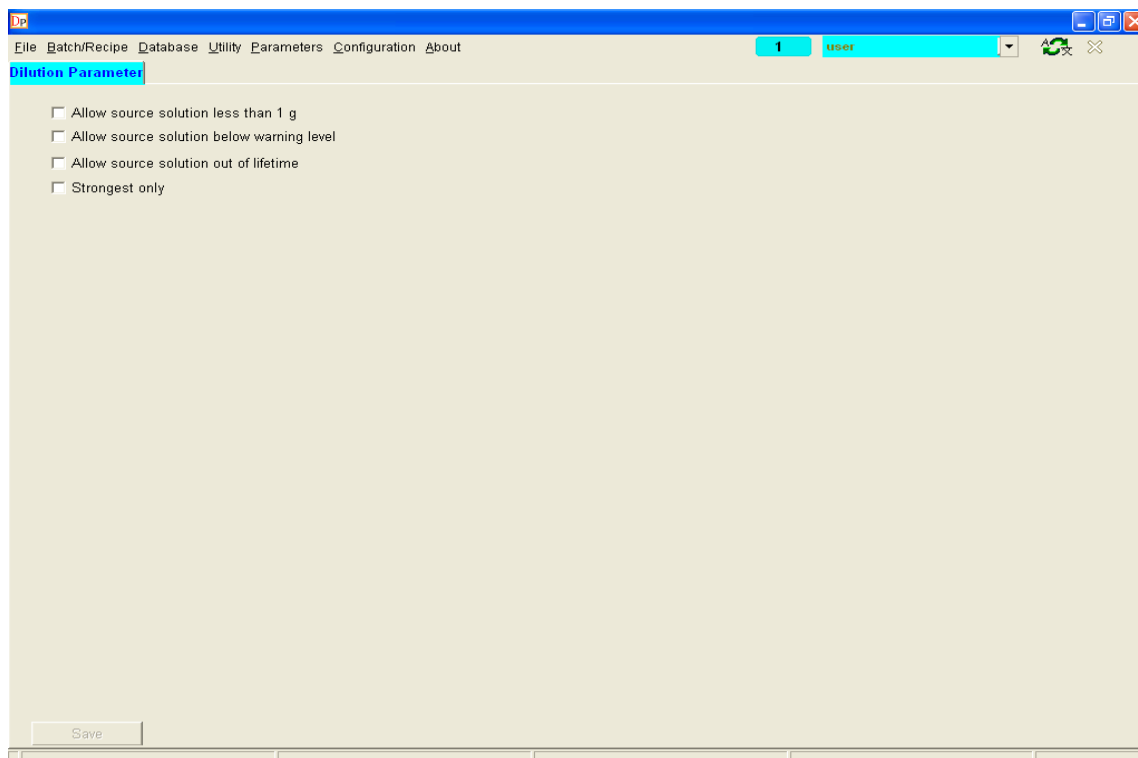

- **Allow source solution less than 1 g**

When enable this parameter, the software will also list the dilution source solution, which the dispense target weight of the solution is less than 1g.

- **Allow source solution below warning level**

When enable this parameter, the software will also list the dilution source solution, which the stock is below warning level.

- **Allow source solution out of life time**

When enable this parameter, the software will also list the dilution source solution, which the life time is out of date.

- **Strongest only**

When enable this parameter, the software will list only the dilution source solution, which the concentration is the strongest.

## 2-2-3 TF DP Program – Parameters – Valve

---

The screen below is “**Chemical**” setting form.

Bottle no.	Gravimetric dispense	Auto change to gravimetric if <g	Enable timeout	Hours
161	<input checked="" type="checkbox"/>	0	<input checked="" type="checkbox"/>	0

Save

With this option, valve volumetric dispense will become gravimetric dispense while match the settings.

If Check weight is enabled, and valve target dispense weight below the weight setting, valve will do gravimetric dispense.

If Check timeout is enabled, and valve has not been dispense exceed the hours setting, valve will go gravimetric dispense.

## 2-2-4 TF DP Program – Parameters – Paste

---

The screen below is “**Paste**” setting form.

Bottle no.	Product code	Enable timeout	Hours	Refresh(sec)
231	Product code	<input type="checkbox"/>	24	
232	Product code	<input type="checkbox"/>	24	
233	Product code	<input type="checkbox"/>	24	

After creating the database for paste, paste auto-refresh time-out settings will appear in the “Paste” page in parameter setup.

If Enable timeout was enabled, when the specific paste has not been dispense exceed the hours setting, the control s/w will refresh the paste pipe according to the Refresh (sec) settings.

## 2-2-5 TF DP Program – Parameters – Import parameter

---

The screen below is “**Import Parameter**” setting form.

This software has the capability of importing recipes from recipe management systems to allow automatic transfer of matching results to the Autolab dispenser, ready for dispensing. This ensures that no errors occur when typing in recipes by hand, as well as ensuring the time involved in the process of matching and dispensing is reduced.

All recipes sent by recipe management system, such as DCI Match and ITM Process are automatically imported to the TF system database. The TF software **automatically checks** the network according to the time interval set, and if recipes are present, then they will be first checked for validity, and if OK, imported into the TF database, available for immediate dispensing.

Any non-valid recipes, for example, containing a product code that does not exist in the TF product database, or a recipe whose concentrations are outside the permitted range, will be logged in the Transfer error Log. All valid recipes will be stored in the TF recipe database, and will be logged in the Transfer log.

These logs can then be used to check which recipes have been transferred, which are invalid, and also allows easy transferral of the valid recipes to a batch.

Recipes transferred from DCI Match and ITM Process will be stored in the TF database using the following condition rules:

TF will check the highest version number existing for that recipe. If no recipe exists, then it will be stored under version number 0. If the recipe exists and the highest version number have been dispensed, and then the recipe sent will be stored under the next version number. If the highest version number has not been dispensed, then it will be overwritten with the recipe sent.

In G2 DP software, it allows to import recipes from more than one source at the same time.

DP

File Batch/Recipe Database Utility Parameters Configuration About

1 user

**Import Parameter**

Import Code  ☐ Activate

Import settings

Import database type

Recipe import path  ...

File name

Sample weight unit ☒ g ☐ Kg

Total volume/weight unit ☒ c.c/g ☐ liter/Kg

☒ Auto import scan time  minutes

☐ Merge blend recipes into one recipe code

☐ Use 24 characters recipe name field in transfer file as recipe code in AutoLabTF

☒ Use recipe trial number

☒ Keep previous

Activate	Import code	File name	Auto import	Minutes	Use tab	Merge blend	Use 24 characters	Use recipe trial number
<input type="checkbox"/>								

Save Delete Load Defaults

- **Import Code**

The code of the recipe import sources. Enable the check box “Activate” to enable the import source and DP software will follow the settings to import recipes.

- **Import database type**

Set the import database type for the import code.

The option “TXT (RCP)” is the format for importing the recipes from Datacolor match or ITM process system.

- **Recipe import path**

Set the path of the imported file for the import code

- **File name**

The file name is fixed with import database type. When the import database type is “TXT(RCP)”, the imported file name is “fm\_ds5.rcp”. You can now change the file name as per your system.

- **Sample weight unit**

The default setting is “g”, if the unit is set to “Kg”, the sample weight inside transfer file will be multiple by 1000.

- **Total volume/weight unit**

The default setting is “cc”, if the unit is set to “Litre”, the total volume inside transfer file will be multiple by 1000.

- **Auto import scan time**

For setting the time interval of scanning imported file.

- **Use 24 characters recipe name field in transfer files as recipe Code**

DCIMatch stores in the transfer file the recipe code used in DCIMatch in two formats. It stores the recipe code in the "recipe code" field, but only up to 12 characters. This limitation of 12 characters is due to the fact the transfer file protocol used is the same protocol as was used for DS5, which had a limit of 12 characters for its recipe codes.

DCIMatch also stores a copy of its recipe code in the "recipe name" field in the transfer file, but up to 24 characters. The DP program allows the use of this 24-character field to be used to create the recipe code in Autolab. Therefore if a 24-character recipe code is required in TF, then the tick box should be enabled.

For example, If the recipe name transferred from Datacolor Match is "12345678901234/188/1" (the original recipe name was "12345678901234/188", "/1" at the end denotes it is first side of a blend recipe), if this is imported with "Use 24 characters recipe name as recipe code in TF" enabled and "Merge blend recipes into one recipe code" enabled, the recipe code stored will become "12345678901234/188".

If the same recipe is then imported with "Use 24 characters recipe name as recipe code in TF" enabled but "Merge blend recipes into one recipe code" disabled, then the recipe code will still be "12345678901234/188/1" (Uses the original recipe name, but then adds "/1" at the end because it is the first part of a blend recipe, which is identified with the suffix /1 in the code because "Merge blend recipes into one recipe code" is disabled.)

In this kind situation, the second imported recipe will overwrite the previous recipe, if the previous recipe has not yet been dispensed. If the previous recipe had been dispensed, the second imported recipe will be imported as same recipe code but will be second dispense shot.

- **Use recipe trial number**

In DCIMatch it is possible to create more than one recipe, for the same recipe code, which is called "trials". This is not the same as "version", because versions are recipes using the same dyes, but with modified concentrations, for example after doing a correction. A trial is a complete different recipe, with different dye combinations. Typically customer may generate 2 or 3 trials at the same time, and want to send them to the dispenser at the same time, and dispense them at the same time in different dye pots. They are alternative recipes to dye to the same color.

The recipe ID is a number, generated within the matching program, which consists of 2 parts. The first part is the recipe ID number, e.g. 279-2, and the last part is the trial version, .1, .2. So a complete recipe ID would be for example 279-2.1, 279-2.2, and 234-1.1 etc. The number, 279-2, or 234-1 is a unique number automatically created by DCIMatch, and has no significance, and is not generated by the customer.

So 279-2.1 is the first trial for this recipe shade, and 279-2.2 is the second trial for the same recipe shade. This recipe ID, including the trial number is now transferred in the transfer file in the recipe code field. The recipe name in DCIMatch is transferred into recipe name field in the transfer file, but the trial number is now added to the end of the name. For customers who want to create more than one trial for the same recipe, we

have to use this .1, .2 coding, otherwise, if they send over 3 trials of the same recipe at the same time, just with the code 279-2, without the .1, .2 and .3, what will happen is that as TFC DP software is importing the recipes, it will overwrite the first trial with the second, because the code would be the same, and then the third will overwrite the second trial. So after importing you would end up with just one recipe, the third formula, under the code 279-2. So the recipe code in TF DP program needs to include this .1, .2 etc, to make the coding unique.

However, for customers who do not want to transfer more than one trial for the same recipe, they may not want to use this trial number. They will probably want to import the recipe using just the recipe name or code, without the trial number. The import function in TF DP program is able to configure it to use or not use the trial number



## 2-2-6 TF DP Program – Parameters – Export parameter

---

The screen below is “**Export Parameter**” setting form.

Export parameter is the function to set the format of the recipes which exported from TF software to customer recipe management software.

In G2 software, it allows to export recipes in different formats to multiple destinations.

Export code: [dropdown] New ☐ Active

Export data type: Standard format

Export path: [text] ...

☐ Extra export path: [text] ...

Recipe select:

- ☐ All Recipe
- ☐ Dispensed (Successful and failure)
- ☐ Dispensed (Successful only)

Re-sort/print:

- ☐ Re-sort shot no
- ☐ Auto print out

Recipe delete:

- ☐ All Shots
- ☐ Not Selected
- ☐ Delete None
- ☐ Rename recipe code as exported

Total : 0 ☒ Keep previous

Active	Export code	Export data type	Export path	Export filename	Extra export path	Extra export filename
--------	-------------	------------------	-------------	-----------------	-------------------	-----------------------

Save Delete Load Defaults Refresh

### • Export Code

The identity code of the recipe export format. Enable the check box “Activate” to enable exporting recipes to the destination.

### • Export data type

Set the export data type for the export code. There are 2 options, “Emulate G1 format” and “Stand format”. The exported recipe format will follow the data type set.

### • Recipe Select

Only the recipes match the requirement will be exported:

- **All Recipe:** Export all the recipes
- **Dispensed (all):** Export all the dispensed recipe, whatever the dispensing result is.
- **Dispensed (No Error):** Only export the dispensed recipes with no error in dispensing.
- **Auto Export:** Enable this function, the dispensed result will be exported automatically, as soon as a recipe is dispensed.

### • Recipe Delete

Delete the recipes from the database, after it has been exported.

- **All Recipe:** Delete all the recipes
- **Not Selected:** Delete the not selected shots.
- **Delete None:** Do not delete any recipe

### • Products to Include

For selecting products will be included in the export recipe, dyestuff, chemical and water.

### • Recipe Sort/Print

It means the system will arrange the Shot No. in sequence after you exported the recipe and delete the non-selected Shot No.

- **Recipe Data to Include**

For selecting data will be included in the export recipe.

## 2-2-7 TF DP Program – Parameters – Batch/Recipe parameter

The screen below is “**Batch/Recipe Parameter**” setting form.

It is the parameters to set the rules for creating recipes and batches.

Priority	From Weight	To Weight
1	0.00	10.00
2	10.00	30.00
3	1.00	3.00
4	0.00	1.00
5	30.00	9999.00

### Recipe Parameter

- **Choose bottle setting**

The 4 dispensing weights setting decides the priority of the bottles to choose when there is the same product solution but in different concentrations. The table on the right side shows the priority of the available bottles. For example, if there are 2 same products in different concentrations in the solution database, when creating recipes, the higher priority bottle will be chosen to dispense.

- **Underflow weight (g):** The underflow limit of dispense target weight.
- **Target dispensing weight low limit (g):** The low limit of dispense target weight.
- **Best target dispense weight (g):** The best target weight to dispense. If there are 2 available bottles with same priority, the software will choose the one closer to this number.
- **Target dispensing weight up limit (g):** The up limit of dispense target weight.

- **Allow save recipe if over total volume**

Enable the option to allow saving a recipe with all products are over recipe total volume.

- **Allow save recipe if underflow**

Enable the option to allow saving a recipe with a product is under “Underflow weight”.

- **Add shot number automatically**

Enable the option; the shot number will be automatically increased by 1 when recalling a recipe for modifying. If disable the option, a button “Add New Shot” will appear beside recipe shot number field, operators have to add shot number by clicking the button.

- **Allow create new shot number when this recipe not dispense yet**

Enable the option, the shot number can be increased by 1 manually or automatically, even the recipe has not been dispensed yet.

- **Allow product selection by product name**


Enable the option, the drop down combo box in the Product NAME field will list/sort the products for selecting the product required and the position of Product Name field and Product Code will be switched in recipe creating screen.

- **Use Quality Code**

Enable the option, the focus will move to Quality field after inputting recipe code in recipe management screen to make operators input the quality. If disable the option, the focus will pass by the Quality field and goes to next field.

- **Default price unit**

There are four units can be choice:

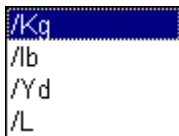

### **Batch Parameter**

- **Show message if have swap or manual bottle**

Enable the option; the system will give users a warning message when batches are transferred from pre-batch list to dispensing queue, if the batches contain swap or manual dispensing bottles.

- **Allow dispense if solution out of lifetime**

Enable the option, the system will allow the batch which contain out of life time solution, be transferred from pre-batch list to dispensing queue.

- **Allow dispense if solution is below low level (100c.c)**

Enable the option, the system will allow the batch which contain low stock solution, be transferred from pre-batch list to dispensing queue.

- **Individual butch list**

The user can only watch his own pre-butches or can watch all user’s pre-butches.


## 2-2-8 TF DP Program – Parameters – Print parameter

The screen below is “**Print Parameter**” setting form.

In G2 software, there is plug-in software called “Autolab UDR”. This software allows operators to create their own report format. DP software will call this Autolab UDR to print out reports.

This parameter is for setting up the templates for different reports. After AutolabUDR software is installed, default templates are ready to use.

The detail of AutolabUDR describe in the AutolabUDR operation manual.

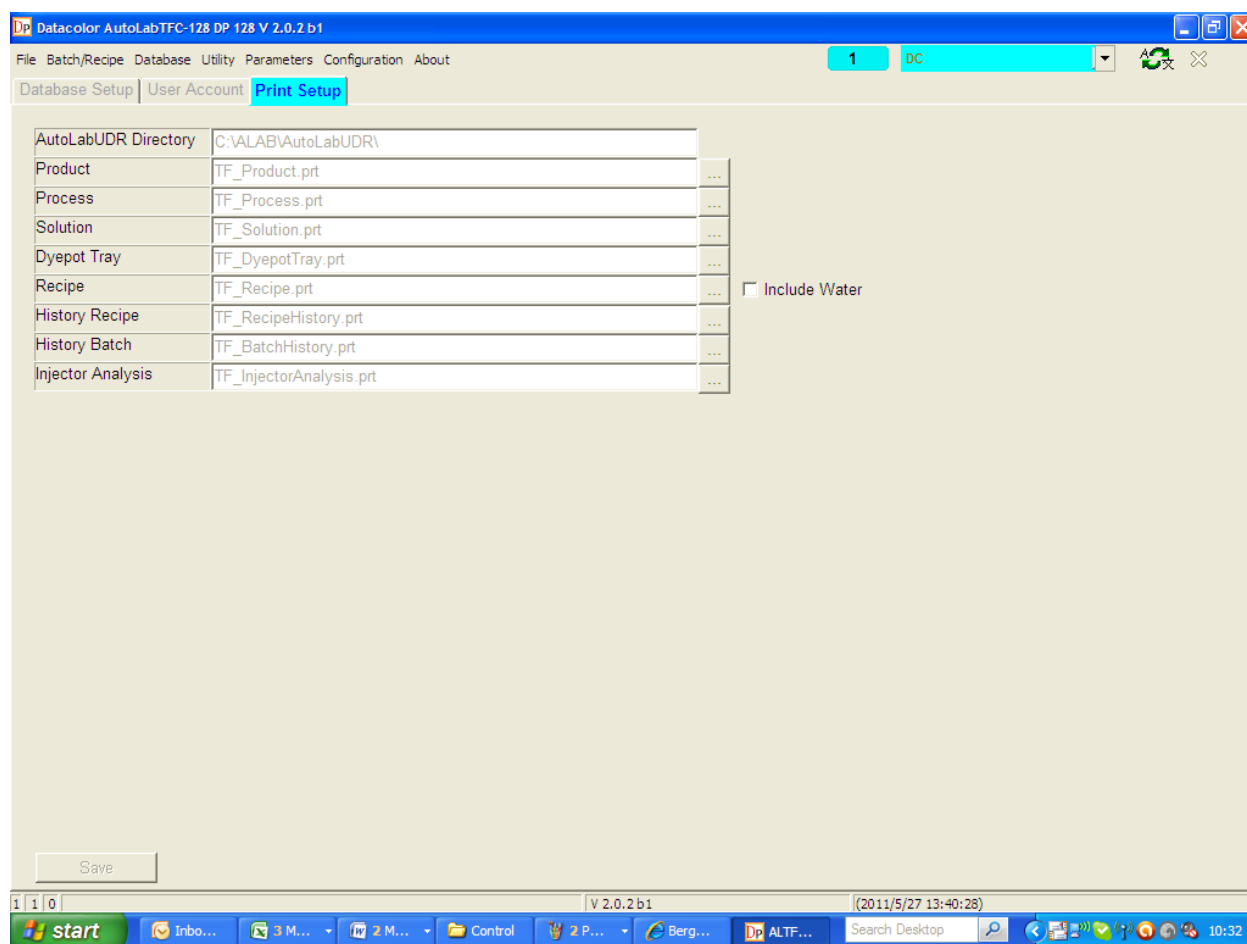

- **AutolabUDR Directory**

DP software will detect the folder C:\ALAB\AutolabUDR\ automatically when launching DP software. AutolabUDR software has to put in this fixed directory.

- **Product**

The name of product report. The default template is TF\_Product.prt.

- **Process**

The name of process report. The default template is TF\_Process.prt.

- **Solution**

The name of solution report. The default template is TF\_Solution.prt.

- **Dyepot tray**

The name of dyepot tray report. The default template is TF\_DyepotTray.prt.

- **Recipe**

The name of recipe report. The default template is TF\_Recipe.prt.

- **Include Water :**

If you want to have water amount displayed in Preview and print from then click this option. In earlier case it was not possible to get water amount on in preview and print in DP program. The water amount would appear only after dispensing is finished and it was possible to get through control program only.

- **History Recipe**

The name of recipe history report. The default template is TF\_RecipeHistory.prt.

- **History Batch**

The name of batch history report. The default template is TF\_BatchHistory.prt.

- **Injector Analysis**


The name of injector analysis report. The default template is TF\_InjectorAnalysis.prt.

## 2-3 TF DP Program – Utility

---

Utility function is for database maintenance and engineering analysis. There are 4 functions in Utility.

- **Backup**
- **Restore**
- **Injector Analysis**
- **Recipe Analysis**

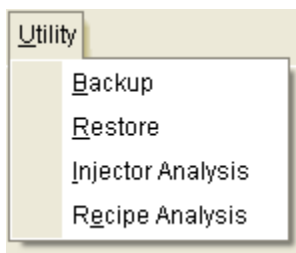




## 2-3-1 TF DP Program – Utility – Backup

---


The screen below is “Utility-Backup” form.

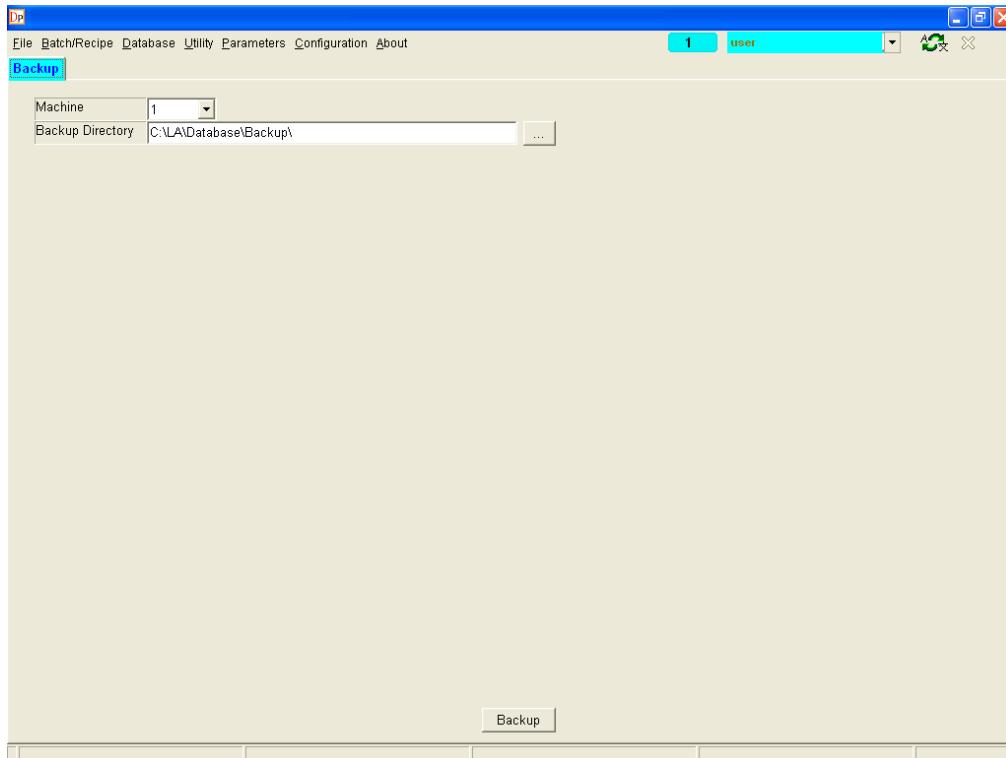

The backup function is to backup database so that it can be protected when data loss happens due to hard disk crash or pc failure.

You can select a machine database to backup, **It is important that the backup data can only be restored to the same machine number.**

**Before restore database, it is important to make sure all users have logged off the machine number.**

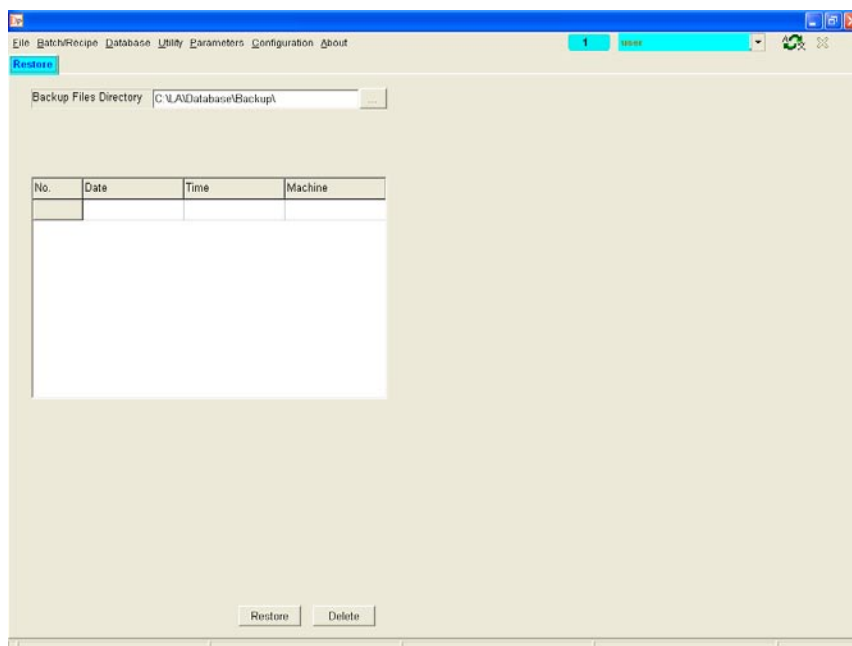

The backup records will be listed in the table. There are information of database backup date & time & machine number.

**It is important that the software only allows restoring the database which backup from the same machine number.**

if you don't have this password then click cancel and system will ask another password. You can input the chief or engineer password and proceed further. In case you don't have chief password then inform Datacolor engineer and he will help to do this task.

## 2-3-3 TF DP Program – Utility – Injector Analysis

The screen below is “Utility-Injector Analysis” form.

This function is for engineers to diagnostic the accuracy of dispensing. Engineers can base on the result of the analysis to find injectors need to be replaced.

After selecting the factor of date range, dispensing target weight and tolerance, click “Select” button.

Injector	Total	Success	Failure	Percent %
1	1	1	0	100
2	1	1	0	100
33	1	1	0	100
121	1	1	0	100

The software will list the times of successful or failure dispensing by injector numbers, also the rate of accuracy dispensing.

The summary of the analysis will be showed on the right down corner to present the accuracy dispensing ratio by the condition selected.

## 2-3-4 TF DP Program – Utility – Recipe Analysis

The screen below is “Utility-Recipe Analysis” form.

This function is for supervisors to find the numbers of recipe, which contain the products, dispensed less than a defined limit. The smaller dispensing target weight will cause inaccuracy dyeing result.

After selecting the factor of date range and dispensing target weight lower limit, click “Select” button.

Target Weight	Record Count	Percentage
0.01 ~ 1.00 g	1	100.00 %

The software will list the quantities of recipes satisfied with the condition defined also the percentage to all recipes.

The summary of the analysis will be showed on the right down corner to present the result.

## 2-4 TF DP Program – Database

---

The option is designed to allow input, modification and set up of all information relative to the products, and solutions used in the laboratory, product, process, etc...



- **Product**
- **Process**
- **Solution**
- **Dyepot Tray**

## 2-4-1 TF DP Program – Database – Product

The screen below is “**Product**” screen.

Input information about all the products, including dyes, chemicals and pastes that are used in the Laboratory.

The screenshot shows the 'Product' screen of the Datacolor AutoLabTF-80 DP 80 V 2.0.2 b1 software. The top menu bar includes 'File', 'Batch/Recipe', 'Database', 'Utility', 'Parameters', 'Configuration', and 'About'. The 'Database' menu is open, showing 'Authority', 'Recipe Analysis', and 'Product' (which is selected). The 'Product' form contains the following fields:

- Product code: CBRFBBS
- Product name: (empty)
- Product type: Dyestuff
- Calculate unit: %
- Min. solution conc.: 0.0
- Max. solution conc.: 100
- Price: 100
- Supplier: ciba
- Color: (red box)
- Modify time: (empty)

Below the form, there is a table of products. The table has columns: Product code, Product name, Product type, Calculate, Price, Modify time, Supplier, Color, and Min. soluti. The table contains 15 rows of data, including products like SOLOL TN, TBK SRL, TBRBU3RL, TBRBUBGE, TEL Blue, TEL Red, TEL Yell, test1234, TOR 2RL, TRD 5G, TRD R, and TVT BL.

At the bottom of the screen, there are buttons for 'Save', 'Delete', 'Load Defaults', 'Export', 'Import', 'Refresh', 'Preview', and 'Print'. The status bar at the bottom shows 'v 2.0.2 b1' and the date/time '(2011/5/27 13:40:28)'.

- **Product code**

Type a code to identify the new dyestuff or auxiliary to input (Maximum 32 characters. It is suggested that if the system is linked to a Color Match system, that the code length is limited to 8 characters.) This code will be used when you are inputting recipes.

- **Product name**

Type a name to identify the dyestuff or auxiliary or paste (max. 48 characters).

- **Product type**

Select a product type. There are 3 types, dyestuff, chemical and paste.

- **calculate unit**

You can select the different recipe calculation unit for the product like %, gpl, g/kg etc

- **Min. solution conc.**

Set the minimum solution concentration. The software will not allow operators to create a solution, which the concentration is smaller than the setting.

- **Max. solution conc.**

Set the maximum solution concentration. The software will not allow operators to create a solution, which the concentration is bigger than the setting.

- **Price**

For setting the unit price of products.

- **Supplier**

For inputting product supplier.

- **Color**

Set the color of the product.

- **Modification date**

For giving you idea about when this product was last modified.

- **Import**

If customer has changed pc and exported the product data from old pc then he can import the data using this function.


- **Export**

Customer can now export product data for future use. The default export file would be in \*.dat but you can change the extension a per your requirement. The export would be in normal ASCII file and can be opened using notepad.

## 2-4-2 TF DP Program – Database – Process

The screen below is “**Process**” screen.

This is the function to define the solution making process.

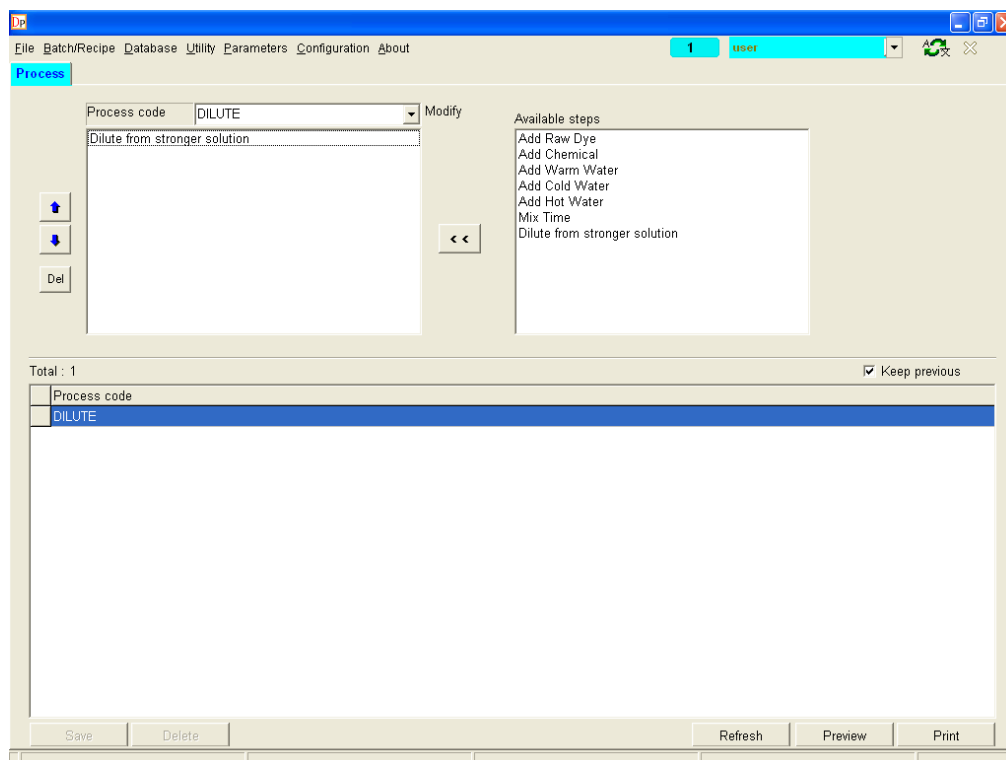

- **Process code**

For selecting a process code or input a new process code.



: For selecting a process code or input a new process code.



: For moving the selected step up.



: For moving the selected step down.



: For deleting the selected step.




The step “Add Raw Dye” can be separated into several steps to fulfill the application of making a high concentration solution. It is possible to add part of the dye, some water, then mix the solution, and then add the remaining required amount of dye required and more water. This can make dissolving of the dye easier.


The way to separate adding the raw dye quantity in more than one step is similar to the way of adding water. After selecting the step “Add Raw Dye” when creating process, a dialogue will pop up and users can key in the percentage of raw dye which they intend to add at this step.

At the end of defining the process, the total percentage of added raw dye must be equal to 100%; otherwise it is not possible save the process.

When a solution process was defined with the sum of the “add water” steps in the process being 100%, and there was a mix step at the end of the process after the last “add water” step, it would complete the making of the solution in the last “add water” step as defined by the operator, before the last mix step, not after the mix step.

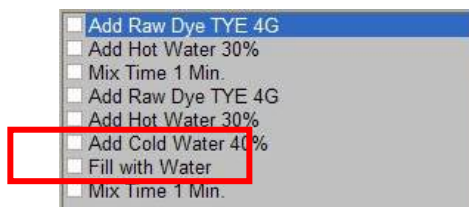

After mixing finished, operators need to click the mixer position on SPS main screen and the software will show a message to confirm the solution making is finished.

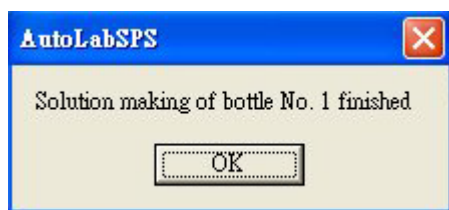

Note that if the total of the steps in a process do not add up to 100%, then automatically at the end of the process, the system will make up to the total amount, in an additional step automatically created by the system, after whatever the last was defined by the operator.

In this G2 software it is possible to make multiple “dilute” processes and also modify the dilute process to add chemical or water as per customer’s need.

### 2-4-3 TF DP Program – Database – Solution

The screen below is “**Solution**” screen.

This is the function to create solutions information for the dispenser.

- **Bottle Group**

In G2 software, bottles are divided into 4 groups, Main Bottle, Valve, Valve\_Paste and External Bottle.

- **Main Bottle:** The solutions in the bottles are dispensed by injectors.
- **Valve:** The solutions in the bottles are dispensed by solenoid valves
- **Valve Paste:** The solutions in the bottles are dispensed by paste valve. Only the product, which the product type is paste can be created in this group.
- **Extend Bottle:** The reserved bottles. The solutions are rare to use can be created in this group. When need it, operators can use quick-switch function to switch the bottle into Main bottle group.

On the right side, it shows the bottle range for the group selected.

- **Bottle No.**

For selecting bottle number.

- **Product code**

Key in the product code which you created in "Product".

- **Product name**

Show the name of the product.

- **Solution Conc.**

Specify the concentration of the solution in this bottle. The concentration has to be within the product minimum and maximum concentration defined in product database.

- **Solution Unit**

You can set four different units of dye solution. It includes

- **%: The quantity of dyestuff in 100 gram dye solution**
- **gpl: The grams of product in 1 litre solution**
- **g/K: The grams of product in 1 kilogram solution**
- **r%: Divide the quantity of dyestuff into the weight of water**

The limits for the input of concentration are as follows:

0.0001 to 100 for %

0.01 to 999 for gpl

0.01 to 999 for g/Kg

- **Default process**

Select a default solution making process for the solution to be made in SPS system.

- **Dilute process**

Select a solution dilute process for the solution to be made in SPS system.

- **Dilute from**

Select a stronger solution as the dilute source for the solution to be made in SPS system.

- **Specific gravity**

Set the specific gravity of the solution. If the solution unit is gpl, then the specific gravity will be calculated when making solution. Normally the value always set to 1.0.

- **Strength**

The manufacturer's standard strength of the dyestuff should be inputted here. Note that this value is used in the calculation of the amount of solution to add based on the standard strength of that product, which is considered to be a relative strength of 100%.

- **High priority**

The system will dispense the solution which set as priority first.

- **Rinse**

Set this parameter to let TFC clean the injector by squeezing it before sucking the solution. Avoiding any sedimentation inside the dip tube of injector.

- **Pull speed**

Set the injector pull speed when the system rinses injectors to match up with the solution

viscosity being dispensed. High viscosity solution needs slower rinse speed. “Normal” is the fast speed; “Low” is middle speed and “Very low” is the slow speed. (Standard default setting set to “Normal”)

- **Life time (hr)**

Set the expiring time of the dye solution. The unit is “hour”. If the solution is out of life time, warning message will show when adding recipes into dispensing queue.

- **Check lifetime**

Set the parameter for the single solution. Choose “Yes” to let software check solution life time when adding recipes into dispensing queue.

- **Amount to make**

Set the solution amount when making the solution in SPS system.

- **Warning level**

For setting the Low Stock Warning Level of the solution.

- **Check level**

Set the parameter for the single solution. Choose “Yes” to let software check solution stock when adding recipes into dispensing queue.

- **Product form**

Set the product form for the product assigned for the bottle number. In Main Bottle, Valve and Valve Paste groups, the product form has to be liquid. In External Bottle group, the product form can be solid, such as salt, which may be added manually.

- **Fill up use hot water**

If set “Yes” at the last step of making solution to fill water up, the hot water pipe will also opened to increase speed.

- **Same solution**

The drop list shows all the bottle numbers where same concentration product assigned.

- **Dispense method**

Define the dispense method of the solution. It can be gravimetric dispense or volumetric dispense.

- **Gravimetric dispense tolerance**

There are 2 ways to identify dispensing tolerance. One is actual gram and another one is percentage of target weight.

When choose actual gram, no matter how big the dispense target weight is, the dispense tolerance will be fixed in the given value.

When choose percentage, the tolerance will be varied by dispense target weight in the percentage given.

- **Approach accuracy**

When enable this function, the software will approach to the accuracy even the tolerance already reached.

## 2-4-4 TF DP Program – Database – Dyepot Tray

The screen below is “**Dye pot Tray**” screen.

This is the function to define the Dyepot tray used in the lab.

Dyepot tray code: ahiba 6 Modify

Dyepot tray name: test

No. of dyepots: 6

Check dyepot capacity: No

Total : 2 ☒ Keep previous

Dyepot tray code	Dyepot tray name	No. of dyepots
test1		6
ahiba 6	test	6

Save Delete Load Defaults Refresh Preview Print

V 2.0.2 b1 (2011/5/27 13:40:28)

- **Dyepot tray code**  
Type a code to identify the dyepot tray.
- **Dyepot tray name**  
Type a name to identify the dyepot tray.
- **No. of dyepots**  
Select the dyepot numbers of the dyepot tray. There are 4 options, 6, 4, 2 and 1 dyepot.
- **Check Dyepot Capacity**  
If you select this option then you can define dyepot capacity and system will check the total volume while recipe calculation and also display message if total volume is more than dyepot capacity.

## 2-5 TF DP Program – Batch/Recipe Management

The screen below is “**Batch/Recipe Management**” screen.

In G2 software, recipe management function and batch management are integrated into one function. Operators can easily manage recipes and batches in the same screen.

The up side on the screen is recipe management and the down side on the screen is batch management.

**Recipe Management Section:**

Recipe code: test14      Quality: [Dropdown]      Shot no.: 1      New Shot

Recipe name: [Empty]      Industry: Exhaust dyeing      Quality Weight (g/Yd): [Empty]

Sample weight(g): 5.00      Weigh      Zero      Chem. recipe: No      Price: 0.00 /Kg

Recipe status: New      Memo 1: [Empty]      Memo 2: [Empty]

Modify/Complete: [Empty]      [Empty]

SIDE 1      (F1 = Insert one row)      (F3 = Insert chemical recipe)      (F12 = Delete one row)

Liquor ratio	No	Product code	Recipe target c	Calc. uni	Bottle N	Target Wt	Actual Wt	Product name
Total Volume	50.00	1 ABL5GL	1	%	* 666	5.00		Astrazon Blue 5GL 200
✓ Dispense water	45.00	2 [Dropdown]						
	3							
	4							
	5							
	6							
	7							

Save      ☐ Keep previous

**Batch Management Section:**

Input Log      Recipe Browse      Batch Browse      Export      Transfer Log      Transfer error Log

Recipe List      Batch List      Dispense Queue

Recipe: [Empty]      Shot: [Empty]      Qua...: [Empty]      Count: [Empty]

Batch List: DC\_110122\_003

Dispense Queue: A [Empty]      B [Empty]

User Code: DC      Dyeport tray: test1      6

0 i 0      V 2.0.2 b1      (2011/5/27 13:40:28)

## 2-5-1 TF DP Program – Recipe Management – Recipe Header

The screen below is “**Recipe Management-Recipe Header**” screen.

TFC Recipe Management allows the creation of exhaust dyeing, continuous dyeing and textile printing recipes.

Recipe code		Quality		Shot no.	1	New	New Shot
Recipe name		Industry	Exhaust dyeing	Quality Weight (g/Yd)			
Sample weight(g)	5.00	Chem. recipe	No	Price	0.00	/Kg	
Recipe status	New	Memo 1		Memo 2			
Modify/Complete							

- **Recipe code**

This code is used to identify the recipe. The code used can be up to 34 characters long.

- **Quality**

It means the quality of the recipe. If users don't use quality, they can disable the setting “Use Quality Code” in batch/parameter function to make cursor bypass quality box when using keyboard creating recipes.

There are 2 ways to input quality code. One is select qualities by drop down list. The qualities in drop down list are pre-defined in database.

The other one is to key in a quality code, which is not existed in quality database, in the box manually. The keyed-in quality code will not be stored in quality database. The quality showed is only a text note.

- **Shot no.**

The shot number of the recipe. There are 2 methods to add shots depend on check item “**add shot number automatically**” on parameter-batch/recipe page checked or not. If the item is enabled, the button “**New Shot**” beside will disappear. If the item is not enabled, the button “**New Shot**” will appear. Users have to click “**New shot**” button to increase shot number.

There is another parameter item setting “Allow to add new shot if the recipe is not complete”. If enable this function and add shot method is “Automatically”, the recipe will add shot number automatically even the recipe is not dispensed after recalling the recipe back. If enable this function and add shot method is “Manually”, the shot number of the recipe will not increase after recalling the recipe back even the recipe is not dispensed yet. But users can click “New shot” button to add new shot for the recipe. Increased shot number must be in sequence.

The number beside “Shot No.” box is dispensing count for this time. It helps users identify how many times they dispense the same recipe at the same shot number. In the same shot number, the dispensing count will increase automatically in sequence. If shot number is changed, dispensing count will be reset.

Situation 1: After recalling recipe “AAA” back, shot number remains “1”. Users can click “New shot” button to add shot number and reset dispensing count. Or use shot number “1” to dispense again and dispensing count will increase 1.

Shot no means the recipe modify times to approach target dyeing result.

Dispense count means how many dispense times of this shot.

- **Recipe name**

The description of the recipe, up to 48 characters in length.

- **Sample weight**

This is the weight of the fabric sample to be dyed. This field is important for calculating the recipe. You can input the weight using the keyboard or press the Weigh button to automatically weight the sample on the Fabric Weighing Scale.

This input box will only appear in exhaust dyeing method.

- **Recipe status**

Show the status of the recipe. This status is for whole recipe status. There are 8 statuses:

1. not dispensed
2. dispensing
3. success dispensed
4. error dispensed
5. exported and not dispensed
6. exported and dispensing
7. exported and success dispensed
8. exported and error dispensed

- **Industry**

There are 3 different industries, exhaust dyeing, continuous dyeing and printing. The default when creating a new recipe will depend on personal parameter setting. Different industry will affect the required information.

Exhaust dyeing: The required must-be inputs are "sample weight", "liquor ratio", "total volume". The allowed solution calculation units are "%", "gpl", and "gm".

When solution calculation unit is "%", use sample weight as calculation factor.

When solution calculation unit is "gpl", use total volume as calculation factor.

When solution calculation unit is "gm", use input number as calculation factor.

Continuous dyeing: The required must-be inputs are "box factor", "pick up ratio", "total length" and "total volume". The allowed solution calculation units are 2, which are "gpl" and "gm".

When solution calculation unit is "gpl", use total volume as calculation factor.

When solution calculation unit is "gm", use input number as calculation factor.

Printing: The required must-be input is "total weight". The allowed solution calculation units are 3, which are "gpk" and "gm".

When solution calculation unit is "%", use total weight as calculation factor.

When solution calculation unit is "gpk", use total weight as calculation factor.

When solution calculation unit is "gm", use input number as calculation factor.

- **Chemical recipe**

You can add some recipes with auxiliaries which you usually use. When you need to add a new recipe with these auxiliaries you can apply the recipes you pre-created and so on



you can save your time. You can use F3 function key to add chemical recipe to existing recipes or you can also select the chemical recipe from list and right click on then select Apply to add chemical recipe to existing dyestuff recipe.

- **Modify/Dispensed**

“Modify” field shows the date and time which the recipe created or modified.

“Dispensed” field shows the date and time which the recipe physically dispensed. No matter success or fail.

- **Price**

The cost of the recipe per **unit** of fabric or per pound in exhaust dyeing. For continuous dyeing, it shows the cost of the recipe per liter of total volume. For textile printing, it shows the cost of the recipe per total weight.

- **Memo1/Memo 2.**

These 2 boxes are for operator input some notes. The title “memo1” and “memo2” can be changed manually. Double click the tile and a message will appear. Operators can change the description of the tile and put a new name for the title. Then users can put required text information in the box, for example: such as “customer name” and “order number”.

## 2-5-2 TF DP Program – Recipe Management – Recipe Detail

The screen below is “Recipe Management-Recipe Detail” screen.

SIDE 1		( F1 = Insert one row ) ( F3 = Insert chemical recipe ) ( F12 = Delete one row )							
Liquor ratio	No	Product code	Recipe target c	Calc. uni	Bottle N	Target Wt	Actual Wt	Product name	
Total Volume	50.00	1 ABL5GL	1	%	* 666	5.00		Astrazon Blue 5GL 200	
<input checked="" type="checkbox"/> Dispense water	45.00	2							
		3							
		4							
		5							
		6							
		7							

Save ☐ Keep previ

New function keys ( F1, F3 and F12 ) visible once you go to product selection code

### Exhaust dyeing

Liquor ratio	10.00
Total Volume	50.00
<input checked="" type="checkbox"/> Dispense water	

- **Liquor Ratio**

This is the required Liquor Ratio of the recipe. This input box will only appear in exhaust dyeing method.

- **Total Volume**

This shows the calculated total volume of the recipe to be dispensed. Total Volume = Sample Weight \* Liquid Ratio. In continuous dyeing, the total volume of the recipe required should be inputted.

- **Dispense Water**

This field indicate how much water will be dispensed in this recipe.

### Continuous dyeing

Total Volume	0.00
<input checked="" type="checkbox"/> Dispense water	

- **Total Volume**

The total volume of the recipe, operators can input the number directly.

- **Dispense Water**

This field indicate how much water will be dispensed in this recipe.

## Textile printing

Total Weight	0.00
<input checked="" type="checkbox"/> Dispense water	

- **Total Weight**

The total volume of the recipe, Operators can input the number directly.

- **Dispense Water**

This field indicate how much water will be dispensed in this recipe.

## Filling in the screen for RECIPE DETAILS

No	Product code	Recipe target c	Calc. unit	Bottle N	Target Wt	Actual Wt	Product name
1	NAOH38	20.00000	%	75	2.00		NaOH 38 Caustic soda
2	RNY GG	1.69609	%	49	0.17		Remazol Navy GG gran.
3	SOD-SIL-	95.00000	%	623	9.50		SOD-SIL-
4	RRD RB	0.64681	%	38	0.06		Remazol Red RB gran. 133%
5	RGYE RNL	0.23829	%	27	0.02		Remazol Golden Yellow RNL gran. 150%
6	LEO SRP	1.00000	%	21	0.10		LEONIL SRP
7							

- No**

The sequence number of product.

- Product code**

Select a Product Code from the combo box. The products shown are those as defined "Database" solutions.

- Target Conc.**

Input the required concentration of the product.

- Calc. Unit**

### Exhaust Dyeing

Dyes: 2 options, % and gm.

Chemicals: 3 options, %, gpl and gm.

### Continuous Dyeing

Dyes: 2 options, gpl and gm.

Chemicals: 2 options, gpl and gm.

### Textile Printing

Dyes: 3 options, %, gpk and gm.

Chemicals: 3 options, %, gpk and gm.

- Bottle No.**

This field is protected. The most suitable bottle number will be automatically selected, once the recipe is saved. Or operators can specify a bottle number for the product. For manual select bottles, a "\*" mark will appear in front of Bottle No. field.

- Target Wt.**

This field is protected. The amount of solution will be calculated and displayed, once the recipe has been saved. If the target weight is under target weight low or high limit, the field will be highlighted in red shade.

- Actual Wt.**


This field is protected. After the recipe is actual dispensed, the actual dispensed weight will be fill in this field automatically.

- Product name**

This field is protected. The Product Name will be displayed, as defined in Product Data.

## 2-5-4 TF DP Program – Batch Management

The screen below is “Recipe Management-Batch Management” screen.

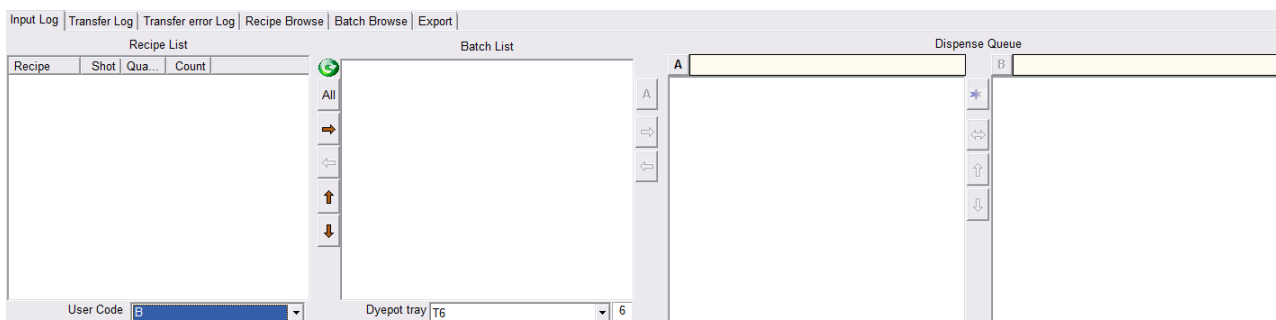


- **Input Log**

The manual created recipes in input log are machine dependent. The created recipes can only be seen by user code selection.

In this log, the user can select multi recipes by using standard Windows multi-select method and transfer these recipes into batches.

To delete the recipes in this log, select the recipes and right click mouse. A command “delete” will appear, click delete command to delete recipes.

- **Transfer Log**

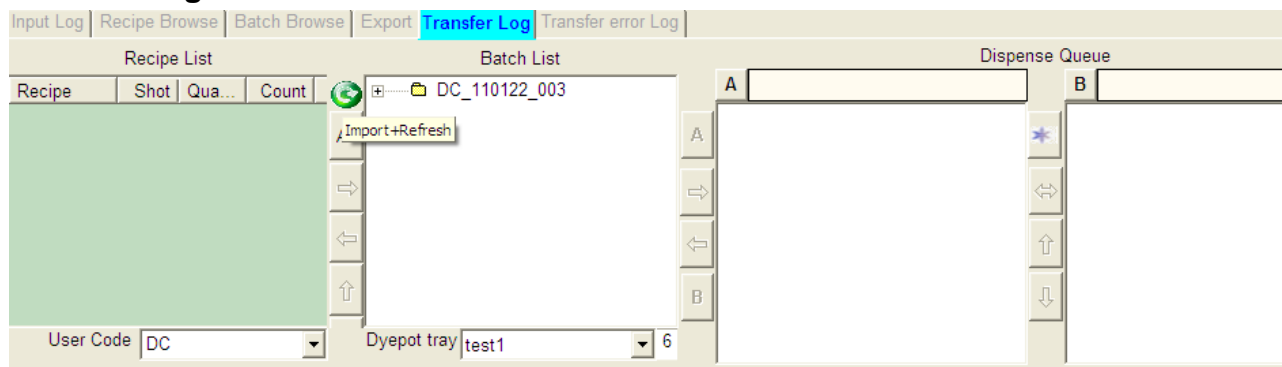

List the recipes which imported from “Datacolor Match” and “Datacolor Process” system. The recipes in transfer log are machine dependent. The recipes in one machine import log can be seen for any authorized users who log in the machine. After modify the imported recipes, the recipe will be moved from Transfer log to Input Log

To delete the recipes in this log, select the recipes and right click mouse. A command “delete” will appear, click delete command to delete recipes.



This button can be used to manual / immediate import or refresh the import.

## • Transfer error Log

Input Log			Transfer Log			Transfer error Log			Recipe Browse			Batch Browse			Export					
Recipe code			Shot			Sd			Recipe code			Error code			Error content			Product code		
1124/3 X.1/1			1			1			1 1124/3 X.1/1			10020			Product code not found in AutoLab solution			test		
1124/3 X.1/2			1			1			2 1124/3 X.1/1			10016			Product code not found in AutoLab product			test		
									3 1124/3 X.1/1			10017			Recipe underflow. Calculated weight of solu			test		
									4											
									5											
									6											
									7											

List the imported recipes which have error information. The software will list the error reason of the recipes. After modify the imported error recipes, the recipes will be also put in Transfer log.

Error code	Condition Description
10011	There are more than 20 products within the recipe.
10012	Unknown Record ID (Invalid recipe code)
10013	The sample weight is out of range. The default sample weight will be stored.
10014	Invalid total volume.
10015	The sum of the concentrations of the products does not match the check sum.
10016	Product code not found in Autolab product database.
10017	Recipe Underflow. Calculated weight of solution to dispense is less than the minimum dispenses weight, as set in Parameter Manager.
10018	Recipe Overflow. Total of the amount of each solution to dispense is greater than the Total volume. The liquor ratio for this recipe will be incorrect.
10019	Recipe to transfer was being edited at the time of the transfer, and the recipe could not be transferred.

## • Recipe Browse

Input Log

Recipe Browse

Batch Browse

Export

Transfer Log

Transfer error Log

User code

Recipe option

Date filter

Recipe code

DC

All recipe

06/17/2011

06/18/2011

Delete

Recipe code	Quality code	Shot no	Dispense count	Modify time	Finish time
test1			1	10/15/2010 07:57:36	
test1			1	10/15/2010 07:57:36	
• 200211-Q3990303-054.1			1	11/23/2010 11:51:19	
• 200211-Q3990303-054.1			2	12/15/2010 23:34:07	
200211-Q3990303-054.1			3	10/05/2010 13:59:14	
▶ 200211-Q3990303-054.1			4	10/05/2010 14:05:02	
200211-Q39903			2	10/05/2010 13:57:21	

Transfer to Batch

Delete

List all saved recipes in the database. There are 3 filters, by user code, by recipe option and by created date. If the operator needs to find a specific recipe, he can input the recipe code in recipe code box to find it.

Users can delete recipes in this screen. Multi-selecting recipes by Windows standard multi-select way is available. Also users can pick up recipes in recipe browse to dispensing directly. After selecting recipes, right click mouse to choose the command "Transfer to batch".

- Batch Browse**

Batch code	Finish time	Recipe code	Quality code	Recipe code	Quality code	Recipe code	Quality code	Recipe code	Quality code	Recipe code	Quality code	Recipe code	Quality code
DC_101015_001		test1											
DC_101123_002	11/23/20	test1											
DC_101123_005		test1											
DC_101210_001	12/10/20	abc											
DC_101213_001	12/13/20	testsm1											
DC_101230_002	12/30/20	test2344											
DC_101230_003	12/30/20	test2344											

List all dispensed batches in the database. There are 3 filters, by user code, by batch option and by created date. If the operator needs to find a specific batch, he can input the recipe code in recipe code to find it.

Users can delete batches in this screen. Multi-selecting batches by Windows standard multi-select way is available. Also users can pick up batches in batch browse to dispensing directly. After selecting batches, right click mouse to choose the command "Redispense". The selected batches will be transfer to dispensing queue.

- Export**

You can export the recipe using this function. Select the recipe code(s) and then select "Export now" and system will export the recipes in export file as per parameters set.

- Dyepot tray**

It needs to select the dye pot tray first before creating a batch. If the dye pots of the selected recipes are not fit in the selected dye pot tray, show warranting message and stop the process. Recipe will remain in the input log.

- Batch List**

Created batches will be put in this box. A batch code is generated by following rule.

Batch code = "user code"\_"date (6 digitals yymmdd)"\_ serial number (3 digitals)




The batches have tree view function, like Windows File Management function. After exploring a batch, the recipes in the batch will be listed. Users can click the recipes. View the recipe content or modify the recipe.

In this box, it will list all users batches by default, if you need to display personal batches in the box, please enable “individual batch list” in the Batch/Recipe management parameter settings.



If there are positions available for inserting recipes, after selecting recipes in Input Log or Transfer Log, just drag the selected recipes into the batch then the recipes will be inserted into the batch.

Users can change the start up time of a “Timer start up” batch by right click mouse after selection batches. A timer will appear and operators can set the time to dispense the batch.

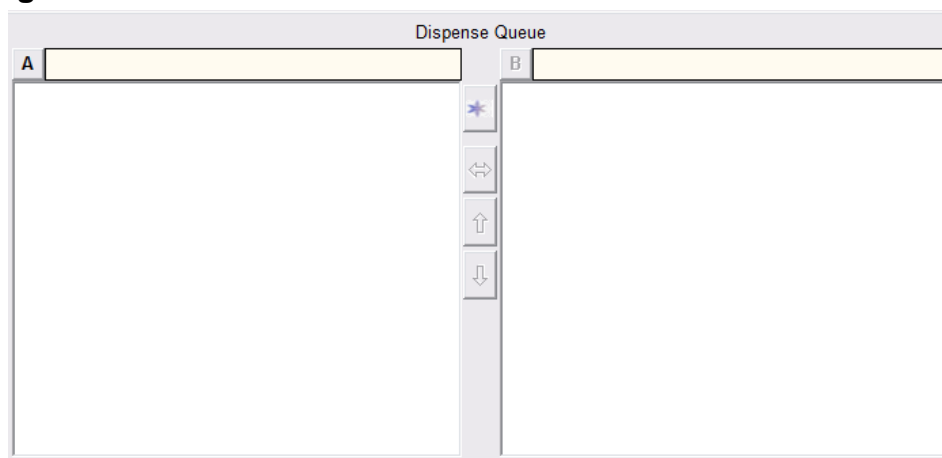- **Dispensing Queue**



There is only 1 dispensing queue in all model TFC and TF40 system. In TF system except TF40, there are 2 dispensing queues for A scale side and B scale side. The box beside “A” and “B” shows the current dispensed batch code.

There is a button “Freeze” . When users need to adjust batches in the batch dispensing queue, in the freeze status, the current batch will still dispensing, but next batch will not dispensing until the user de-freeze the system by pressing  again.

Batches can be moved from one queue to another queue by using the 2 direction arrow button.

2 buttons, “up arrow button” and “down arrow button” will allow user to move batch up and down.

If you have more than 1 bottle with the same dye solution and concentration, when transferring batches from “batch list” box to “Dispense queue” box, the software will identify if there are bottles can switch when the original is under low level or out of life time. bottle priority rules which are listed below to transfer batches and swap same solution bottles automatically if needed.


The searching slave bottle rules are, in order of priority.

1. The system will use the bottle which is in injector dispensing area.
2. The system will use the bottle which has shortest remaining life time.
3. The system will use the bottle where the solution quantity is above the safety stock (200g default).
4. The system will use the bottle where the solution quantity is between safety stock (200g) and minimum stock (100g).
5. If no bottle exists in the dispense area that fulfills the above criteria, then the system will then use a bottle from the extended bottle area, and follow the rule point 2, 3 and 4 to identify the extended bottle priority and selection of the bottle.


The priority of the slave bottles can be changed anytime due to the solution life time and stock levels continue to change.

The identifying slave bottle priority rules are presented in flow chart below. The software will identify the priority of the bottles in slave bottle list then use the highest priority bottle when the master bottle can't pass solution checking.

The flow chart below is the complete algorithm of transferring batches from Pre-batch box to Work-batch box. This flow chart will base on the settings in "Batch settings" to identify if the batches can be transferred to work batch box. If the solution in a batch has same solution slave bottles for replacement and it needs to be replaced, the software will follow the flow chart in next page to identify the priority of slave bottles then use the highest priority bottle to replace the master bottle.

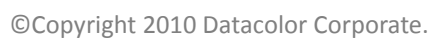

**The algorithm of swapping to slave bottle**

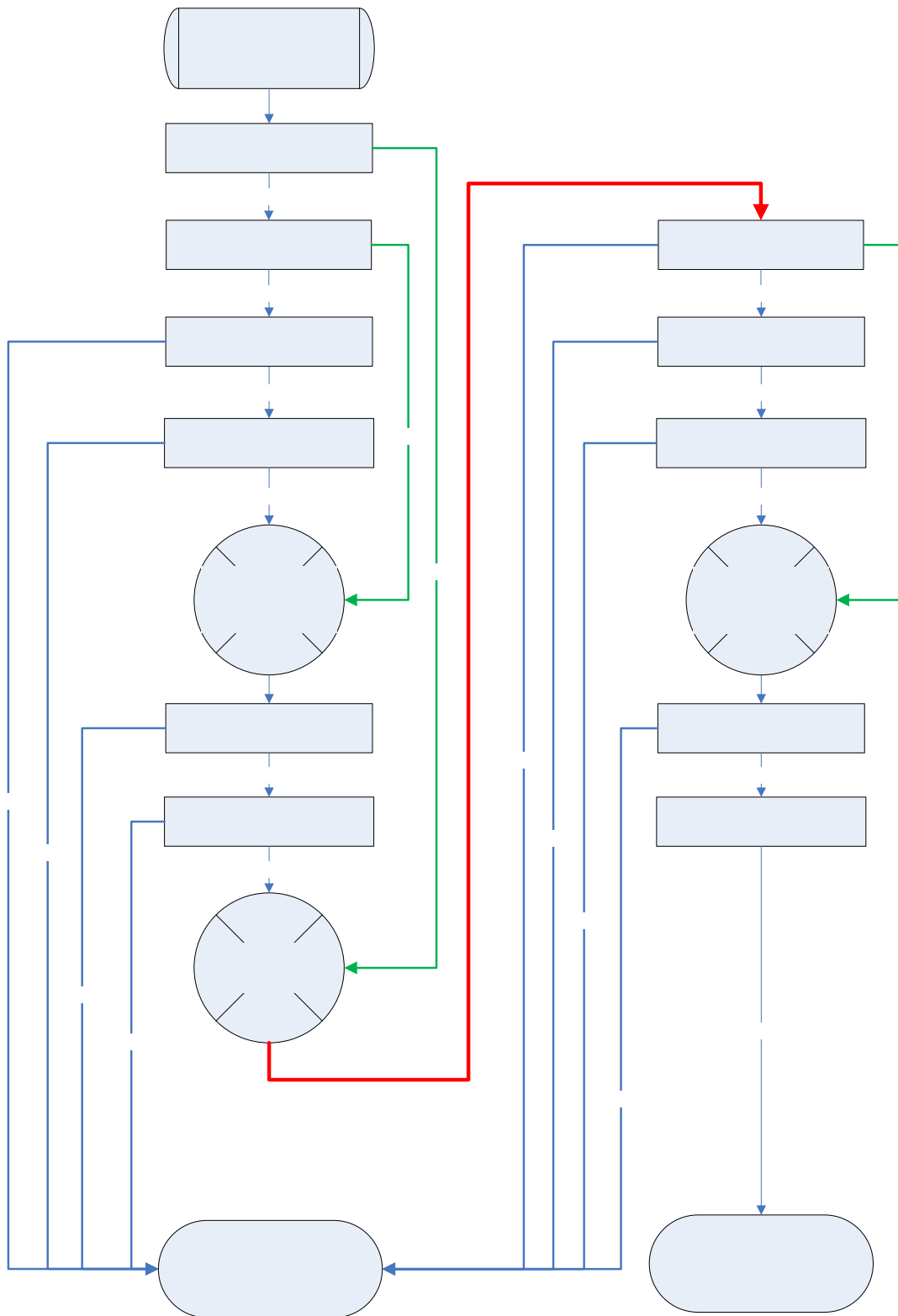

**TFC DP Program – Batch Management – Import Blend Recipes**

The rules of importing blend recipes from Datacolor match or ITM process is the same as G1 software. Take an example below

```

500M171506.1/1  10F  569080 Flick 239.E.1/1  00.000  84.000YN0120070404
DN
500C171506.1/1  1G  DST120  0.50000
500C171506.1/1  2G  C502    2.00000
500C171506.1/1  3G  DST102  1.93640
500C171506.1/1  4F  DST103  0.44592
500C171506.1/1  5   C621    10.00000
500C171506.1/1  6F  18-D/103 0.00000
500C171506.1/1  7F  19-D/301 0.04096
500C171506.1/1  8F  20-D/400 0.05200
500C171506.1/1  9F  18-D/504 0.13000
500C171506.1/1  10W  WATER   15.10528
500M171506.1/2  4G  569080 Flick 239.E.1/2      84.000YN0120070404  DN
500C171506.1/2  1G  DST118  3.00000
500C171506.1/2  2G  C131    3.00000
500C171506.1/2  3G  DST115  1.00000
500C171506.1/2  4W  WATER   7.00000

```

Recipe “171506” is a 2 sides blend recipe.


When the import parameter set “Merge recipes”, there will be only one recipe, which is recipe “171506”, imported. The recipe code is “171506” and the products of 2 sides will be put in the same recipe content and there is an empty line between products of 2 sides for identification.

When the import parameter set “Don’t merge recipes”, there will be 2 recipes imported into G2 recipe database. One is recipe code “171506/1” and it contains only the products of side 1. Another one is recipe code “171506/2” and it contains only the products of side 2. G2 software will treat these 2 sides as 2 different recipes.

The industry of the recipe will be identified automatically.

## 2-6 TF DP Program – Solution Stock

The screen below is “Solution stock” screen.

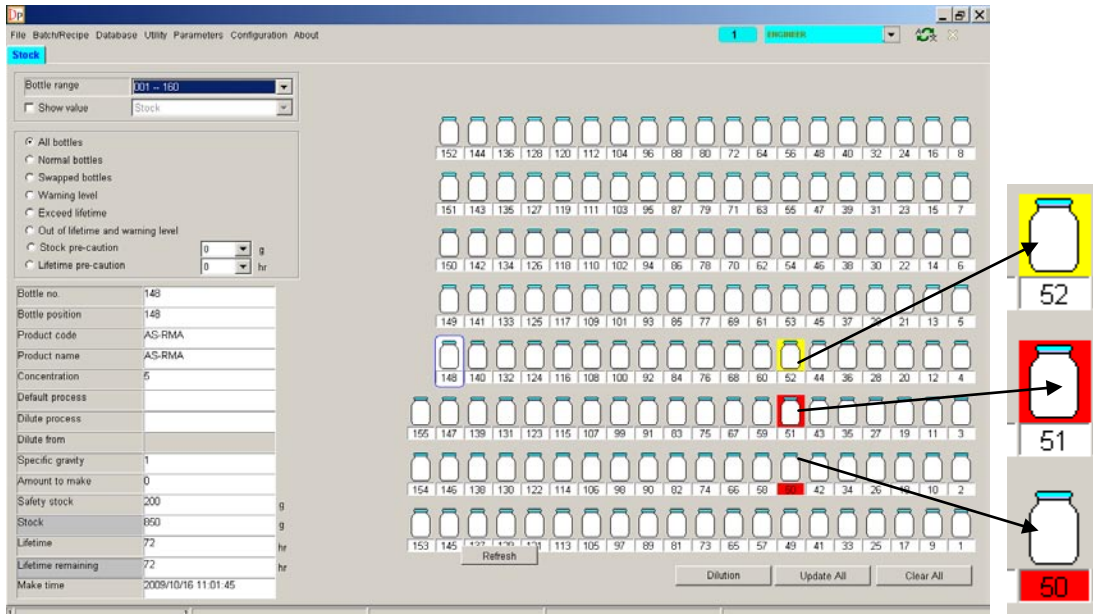

After moving mouse cursor on a bottle, the information of the solution in the bottle will be showed on the left side of screen.

There are filters can be chosen on the left side. By different filter, the bottles will show different information.

When there is solution, which the stock is in warning level, yellow shade will surround the bottle, such as the bottle 52 in the picture above.


When there is solution, which the stock is in low level, red shade will surround the bottle, such as the bottle 51 in the picture above.

When there is solution is out of life time, the background of bottle number will be shade in red color, such as the bottle 50 in the picture above.

Select a bottle and right click mouse, there are 3 functions can be selected.


- **Set Stock**

It is an easy way for operator to see all the information of the selected solution, and update the stock individually. Input the weigh of the new solution, and click ok to save the stock.

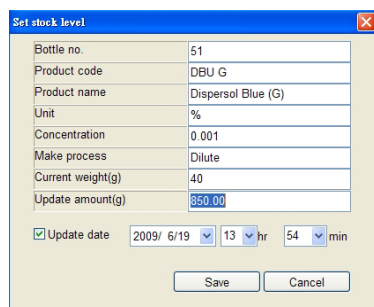

- **Swap Bottle**

This is the function of swap extended bottle to dispensable position. The extended bottles are from 241 to 9999.

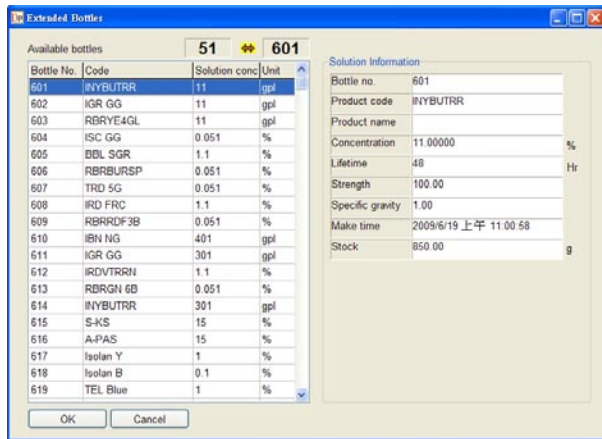

Select an extended bottle in the list of left, click “OK”. The bottle is swapped into the dispensable position, and original bottle number will be changed to the number of the extended bottle, and the color of bottle is change to black. See the picture below.

- **Swap back**

It is for reset a swapped bottle back to original bottle.

- **Make Dilution in TF**

Using special bottle trays, holding 2 empty solution bottles at a time, the system can automatically make dilutions, by loading the bottle tray, dispensing the required amount of stronger solutions from the relevant solution bottle on the system using the injector, then dispense cold water from the water tank to make the solution up to the required final solution strength. Once completed, the solutions can then be taken off the bottle trays and placed on the dispenser system in the correct bottle position for that solution strength. There are 2 ways to enter the “Dilution” screen in DP program, for making dilutions. The first way is to first enter the stock status screen, and click on the “Make Dilution on TF” button. The second way is click on the “Make Dilution on TF” button in solution stock and life time alarm list screen, when transferring recipes from batch list to dispense queue.

Product code	Target bottle no.	dilute target conc.	Unit	Dilute from	dilute source conc.	amount to make
<input checked="" type="checkbox"/> TYE 4G	1	0.05000	%	6	1.00000	850.00
<input type="checkbox"/> TOR 2RL	2	0.05000	%	7	1.00000	850.00
<input checked="" type="checkbox"/> TRD R	3	0.05000	%	8	1.00000	850.00
<input type="checkbox"/> TVT BL	4	0.05000	%	9	1.00000	850.00
<input type="checkbox"/> RYE R	12	10.00000	gpl	17	100.00000	850.00
<input type="checkbox"/> TBRBLJRL	13	0.05000	%	18	1.00000	850.00
<input checked="" type="checkbox"/> DRD C4G	14	0.05000	%	19	1.00000	850.00
<input type="checkbox"/> RYE GR	15	10.00000	gpl	20	20.00000	850.00
<input checked="" type="checkbox"/> BRD S3B	16	0.05000	%	19	1.00000	850.00
<input type="checkbox"/> BYE S8C	23	0.05000	%	24	1.00000	850.00
<input type="checkbox"/> ROYE RNL	27	15.00000	gpl	28	500.00000	850.00
<input type="checkbox"/> LBRUS EB	29	0.05000	%	5	3.00000	850.00
<input type="checkbox"/> BRU SGW	32	0.05000	%	31	0.50000	850.00
<input type="checkbox"/> BGN S4B	34	0.05000	%	35	2.00000	850.00

Product code	Target bottle no.	dilute target conc.	Unit	Dilute from	dilute source conc.	dilute source weight	water weight	amount to make
TYE 4G	1	0.05000	%	6	1.00000	42.50	807.50	850.00
TRD R	3	0.05000	%	8	1.00000	42.50	807.50	850.00
DRD C4G	14	0.05000	%	19	1.00000	42.50	807.50	850.00
BRD S3B	16	0.05000	%	10	1.00000	42.50	807.50	850.00

Lists of solutions that can be made up by diluting from a stronger solution with then appear.

To change the default amount, double in the column “amount to Make” and you can sett the desired amount.

To further filter the solutions in the list, you can enable the check box of “Low level” and “Out of life time” to list only those solution which fulfil these criteria.

Select the bottles which are required to be made up from the list, and the solution selected will appear on the “selected bottle” list on the right side of the screen. If you want to change the amount of the solution to make up, double click on the amount box of the bottle in available bottle list.

After selecting all the solutions required to be made up, click on “OK”, and this will then open the “Dilution batch generator screen”. This dilution batch generator will automatically create a dilution batch code for every solution selected, DiXXXX-YYYZZZ, where XXXX is a dilution batch sequential number. YYY is first solution number to make in this batch. ZZZ is the second solution number in this batch (each batch consists of 2 solutions to make, since each bottle tray holds 2 bottles). In this screen, you can also see the detail of dilution bottle and decide whether put the batch as first the batch in dispense work batch. Click “OK” to move the batches to the work batch. The dilution batches can not be transferred back to the pre-batch box, like normal dispense batches can be.



## 2-7 TF DP Program – History Report

The screen below is “History Report” screen.

Operators can use this function to check all dispensed reports, by batch or by recipe

Batch | Recipe | Dilute

Date Range: 2008/10/1 – 2009/10/20 Search

Select User: User Code SUPERVISOR

Select Result: ☐ All ☒ Successful ☐ Failed

Batch code	Date	Recipe code	Quality code	Shot no	Error
SUPERVISOR_0900001	2009/10/16	1		1	
SUPERVISOR_0900002	2009/10/16	2		1	
SUPERVISOR_0900003	2009/10/16	3		1	
SUPERVISOR_0900004	2009/10/16	4		1	
SUPERVISOR_0900005	2009/10/16	5		1	
SUPERVISOR_0900006	2009/10/16	6		1	
SUPERVISOR_0900007	2009/10/16				
SUPERVISOR_0900008	2009/10/20				
SUPERVISOR_0900009	2009/10/20				
SUPERVISOR_0900010	2009/10/20				
SUPERVISOR_0900011	2009/10/20				
SUPERVISOR_0900012	2009/10/20				
SUPERVISOR_0900013	2009/10/20				
SUPERVISOR_0900014	2009/10/20				
SUPERVISOR_0900015	2009/10/20				
SUPERVISOR_0900016	2009/10/20				
SUPERVISOR_0900017	2009/10/20				
SUPERVISOR_0900018	2009/10/20				
SUPERVISOR_0900019	2009/10/20				
SUPERVISOR_0900020	2009/10/20				
SUPERVISOR_0900021	2009/10/20				

Product code	Recipe ts	Target Wt.	Actual Wt.
A-IW	1	5	5
Water	0	45	45

Total: 26

Preview Print

### Batch Report

In the batch report screen, users can filter the batches by selecting “User”, “Dispensing Date Range” and “Dispense Result”.

Once chosen the “User”, “Date Range” and “Dispensed Result”, click on the “Search” button. The program will follow the condition you choose and find out the batches by the set condition.

In the screen, the box on the left will show you the batches which dispensed by the selected user in the date range. The box will show you the information of “Batch Code”, “Dispensed Date” and “Dispensed time”.

Select a batch and the recipes in the batch will appear in the box on the middle. The box on the middle will show you the “Recipe Code” and “Shot No.” information.

Select a recipe in the middle box and the recipe detail will appear in the box on the right.

Click “Print” button, the program will print a batch report.

## **Recipe Report**

In recipe report, you can filter the recipes by “Dispensing Date Range” and “Dispensed Result”. Or if you need find a specific recipe, you can key in the recipe code in “Select Recipe Code” box. The recipe you want to find will appear.

Once chosen the “Date Range” and “Dispensed Result”, click on the “Search” button. The program will follow the condition you choose and find out the recipes by the set condition.

In the screen, the box on the left will show you the recipes which dispensed in the selected date range. The box will show you the “recipe code”, “Shot No.” and “Dispensed Date” information.

Select a recipe and the batch which content the recipe will appear in the box on the middle. The box on the middle will show you the “Batch Code” If the recipe with the shot number has been dispensed in more than one different batch, all the batches will appear in the box on the middle.


Select a batch in the middle box and the recipe detail of the specific dispensing will appear in the box on the right.

Click “Print” button, the program will print a recipe report.

## 2-8 TF DP Program – File

---

The screen below is “File” screen.

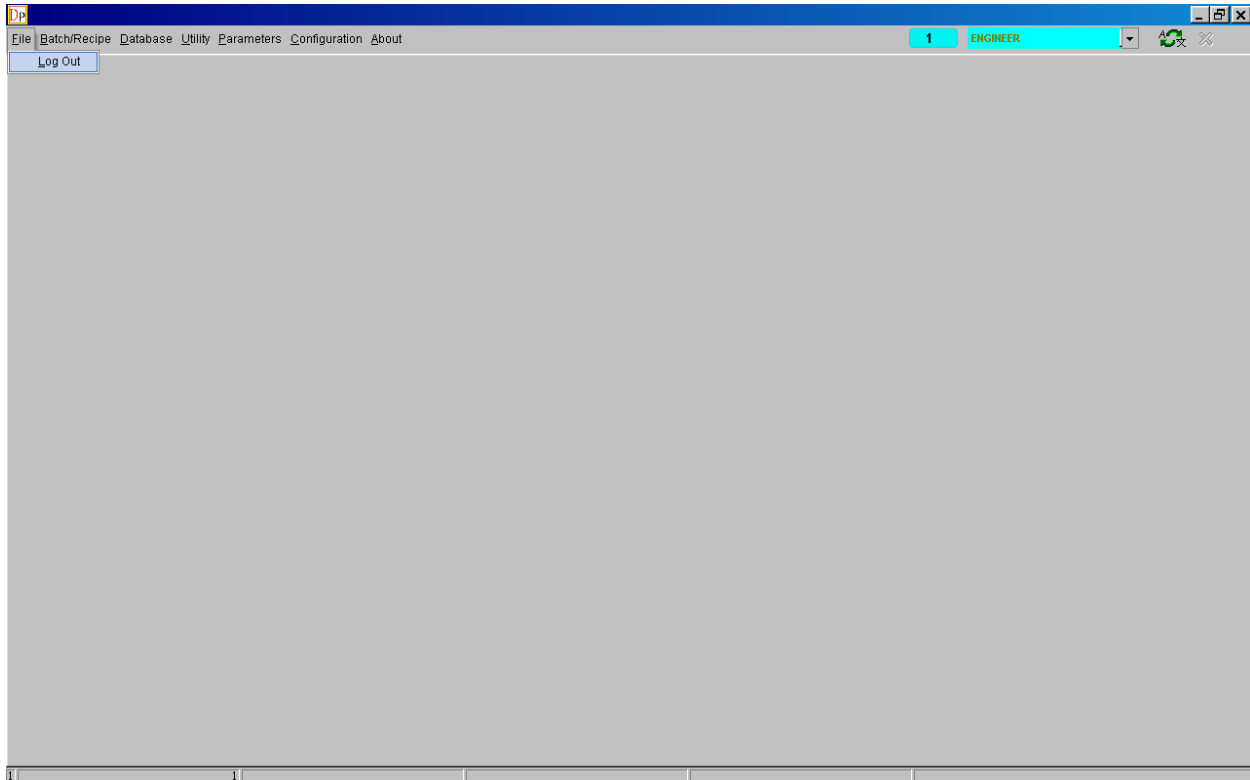


There is only one function “Log out” in File menu.

Note after selecting log out function will log out all users.

## CHAPTER 3: TF/TFC Ctrl program

### 3-1 TF/TFC Ctrl program – Overview

The Control Program is the main program for controlling and maintaining the Dispenser. The main screen looks like this

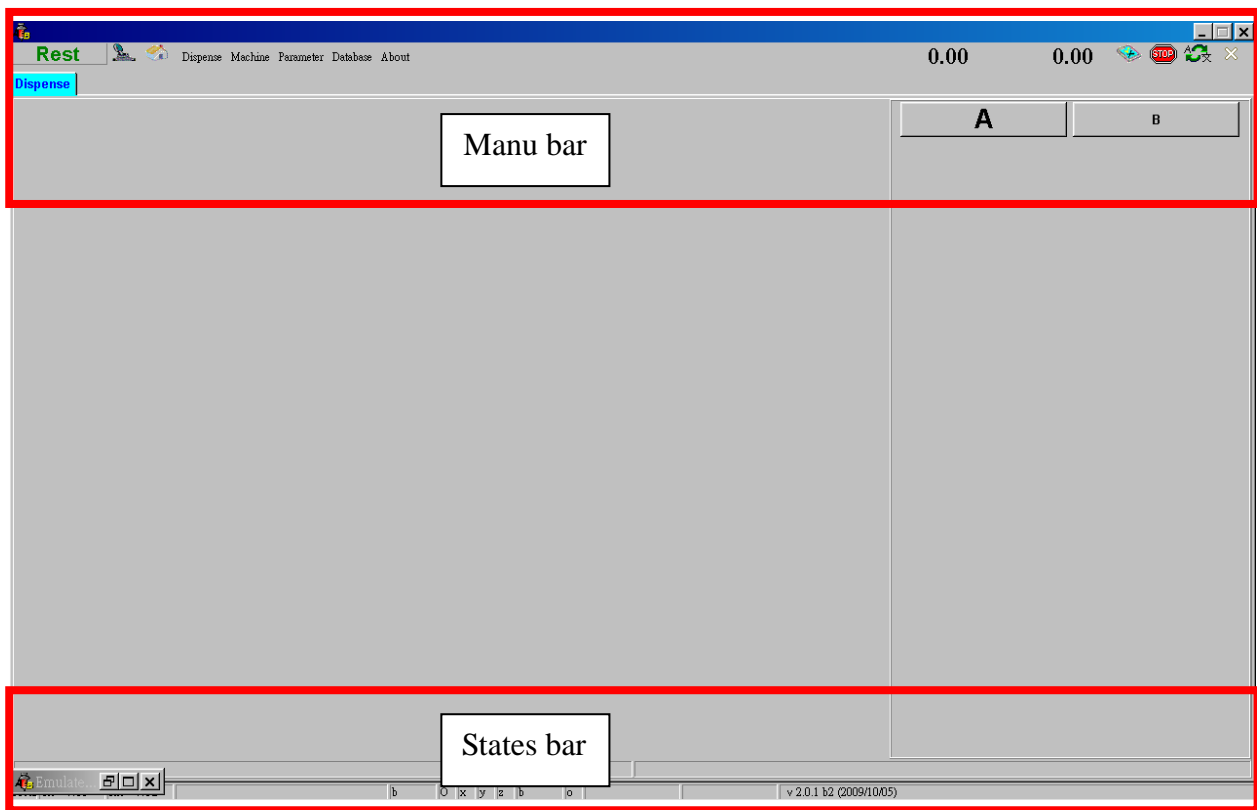

## TF/TFCCtrl program – Overview – Manu Bar

---

### Status :



The system is in running mode.



The system is in stand by mode.

### Home:



Reset the x, y and z back to home positions. Operator has to click the button every time after starting the Control program.

**Dispense:** Show the existing condition of butch dispensing.

**Machine:** Diagnostic the system, and calibrate the x, y positions of injectors.

**Parameter:** For setting parameters.

**Database:** For checking the dispensing histories.

### Scale:



For switching the view of Module A and B. Big letter means the current view of which Module. (This is only for TF systems, and will not show in TFC systems)

### Emergency Stop :



Click this button for stopping all the action of TF machine.

### Language Setting :



Select Language, start the language selection form.

### Exit :



Exit, close the Control program.

## Warning :



Warning, if there is a problem with the valve coefficient and they need rechecking, the warning button will appear. Click on the button for more information of which valve has problem.

## Dispensed Log :



Click the button to view the dispensed log.



Delete the selected dispensing record. First select the batch/batches that are required to be deleted from the list by clicking on the relevant batch, using “Ctrl” or “Shift” key to multi-select batches. Then press the delete button to delete these batches from the dispense history.



Print the dispensing record. Only one batch can be selected at a time.




Show the successful dispensed records.



Show the unsuccessful dispensed records



Sort batches by batch code.

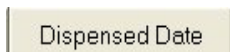

Sort batches by dispensed date.



Show recipe detail. When printing out recipe detail, the recipes in selected batch will be printed out individually. Below is recipe detail screen shot.



Shows dilution solution batches record. This option shows all the solutions made by dilution method on TF and will also show the actual concentrations of solutions made.

**Dispensing Monitoring Screen :** It shows you the process of dispensing.

## **TF/TFC Ctrl program – Overview – Status Bar**

---

The bar on the bottom only shows communication status between computer and system.

**Local :**

It means that TF is operated by host PC.

**Tx / Rx:**

Showing the communications between PC and TF machine

Tx: transmitted counts, Rx: received counts.

If the system is still online, the number will keep changing rapidly.


**Version:**

It shows you the version of the software

## 3-2 TF/TFC Ctrl program – Dispense

### The procedures before dispensing:

At DP program, moving a batch from **Batch list** (in **Current Batch Management**) to (**Current Work Batch**), a \*.A file will be created in “C:\ALAB\Exchange\ALTF\PreBatch” directory. Control Program will scan if there is \*.A file present in the “PreBatch” for every 5 seconds. The control program will start the dispensing automatically once find the \*.A file. The screen below is the Control is dispensing recipes:

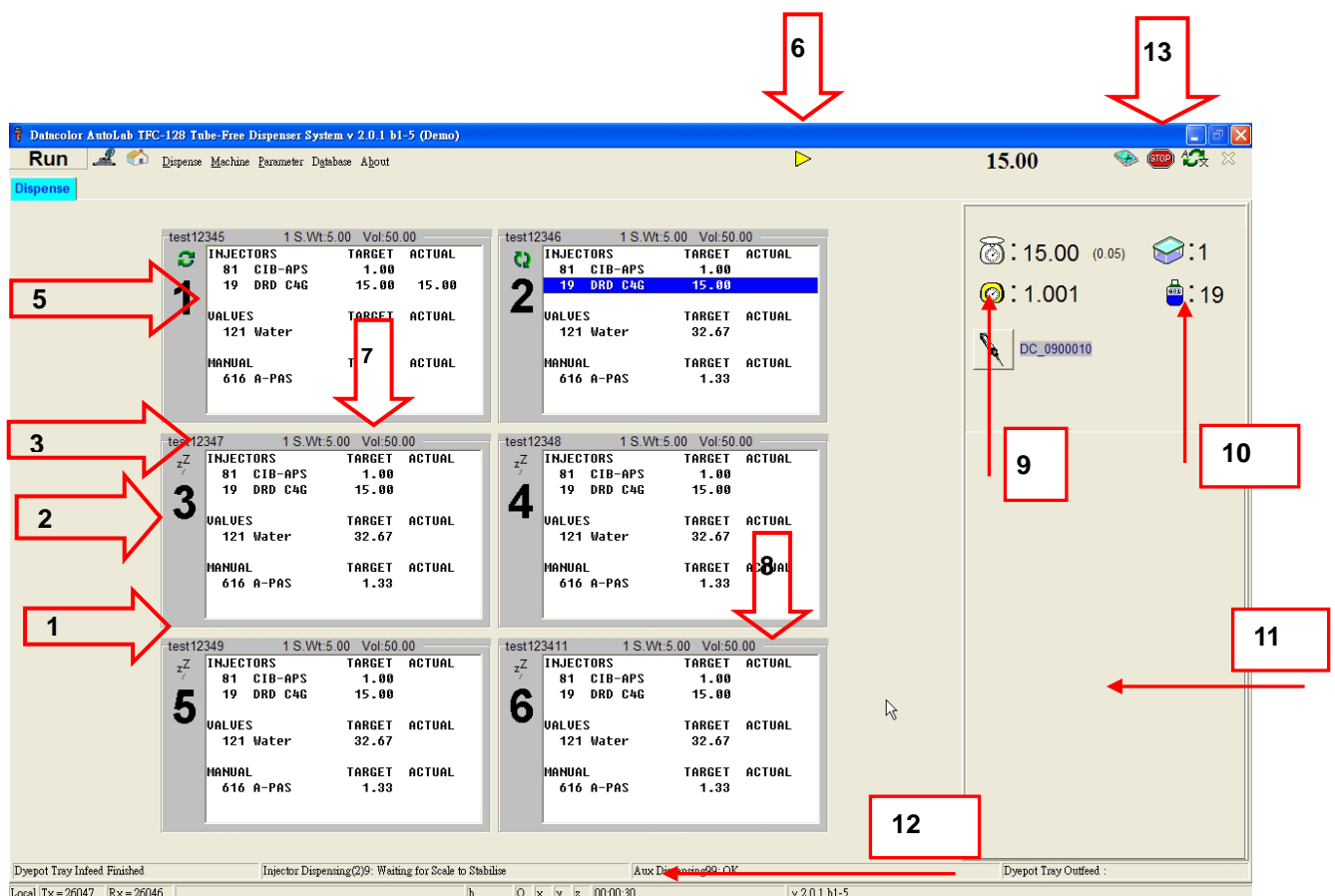


1. **Recipe Code** : The recipe code of the recipe is being dispensed in this dye pot. The number beside on the right side is Shot No.
2. **Dye Pot Number** : The dye pot position on the dye pot tray.
3. **Dispensing by Injectors** : The products dispensing by injector.
4. **Error Dispensing** : If the dispensing result is over or under the tolerance, there will be a plus or minus mark showed beside the injector/valve number. The plus mark means over dispensing and the minus mark means under dispensing. The system will warn you if the recipe need re-dispense. Follow the instruction and change the dye pot to re-dispense the recipe.
5. **Under Dispensing Solution** : The blue blank will cover the dye solution when it is



under dispensing status.


6. **Pause Button** : Click this button will stop the action of dispensing temporarily and the button will change to “Play” button. Click “Play” and machine will continue dispensing.
7. **Target Weight** : the target dispensing weight of the dye solution.
8. **Actual Weight** : the actual dispensing weight of the dye solution.
9. **Specific Weight** : The specific weight of the current solution being dispensed. The system will calibrate the specific weight automatically by the weight it dispensed.
10. **Bottle Number** : The solution bottle number which is being dispensed.
11. **Waiting List** : The batches here will be dispensed in sequence.
12. **Dispensing Time** : The time has been used for dispensing the batch.
13. **Emergency Stop** : If the dispensing is stopped in the middle by clicking on the button, all actions of TF will be stopped and the batch will be returned to DP program. You need to re-arrange the batch. A “confirmed window” will be show for put the syringe back to solution bottle.
14. **Empty bottle warning**: When dispensing solutions by injector, after the first dispense shot, if the actual dispensed weight is at least 1g less than it should be (the target weight), then the control program will show a warning message to inform the user. Typically an undershot so large on the first shot indicates the solution bottle may be empty, or there is a problem with the injector. The warning message states this, and the user can then check the solution quantity and injector condition. The warning message also gives the option to abandon dispensing, or continue.

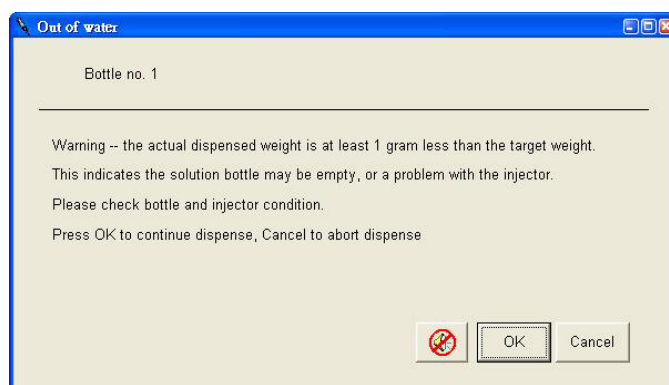

### 3-3 TF/TFC Ctrl program – Machine

#### 3-3-1 TF/TFC Ctrl program – Diagnostic

This is for engineers to diagnostic the system only. If you do not so familiar with TF, please don't use this function avoiding causing damages. The system will ask you to key in a password for opening it.

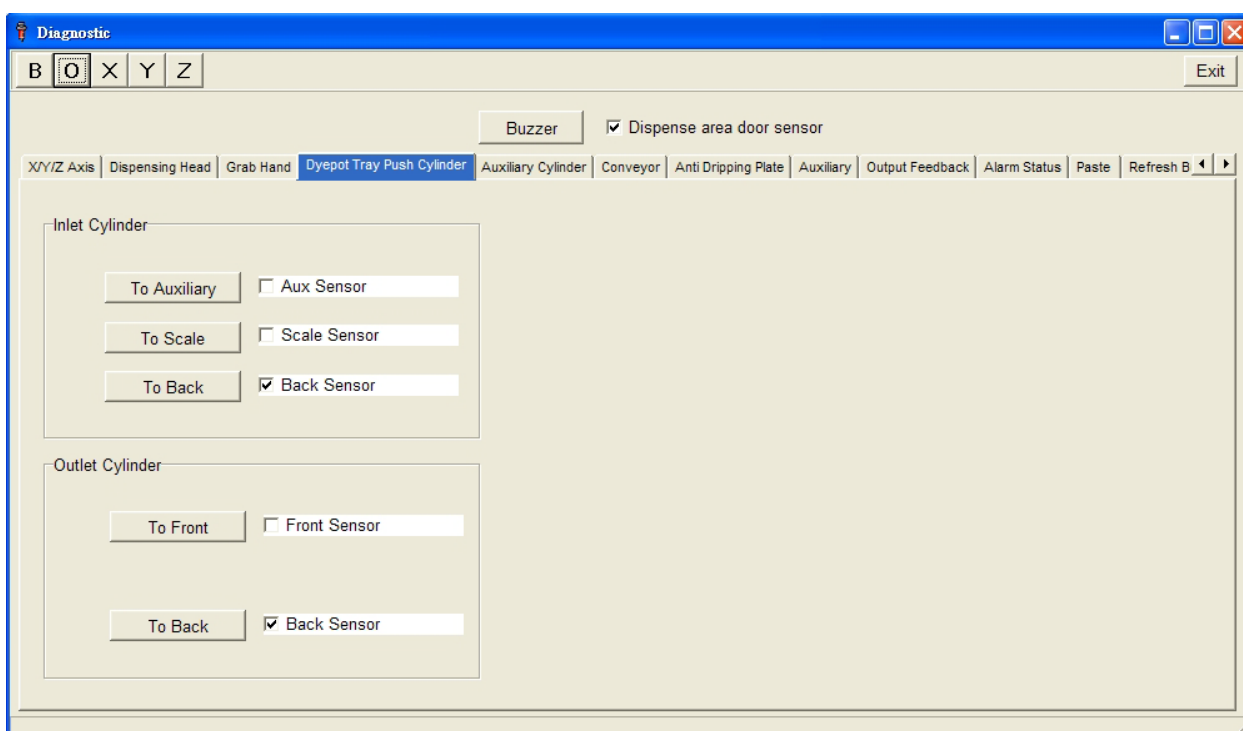

#### Online Buttons Bar:



Click on the buttons for connecting the branch control system.

“b”: balance

“o”: outputs

“x, y and z”: the three axis of robot arm

Once the system is connected, the letters on the bar will be changed to capital letters. See the picture below:



#### Warning Message :

The warning message here for warning that damage may be caused if you send a command to TF.

#### Exit :

Click the button to exit the diagnostic.

**Dispense area sensor :**


The tick box will be checked if all the safety doors are closed.

**Buzzer :**

Click the button to test the buzzer.

## TF/TFC Ctrl program – Diagnostic – Output feedback


The picture below is the “Output Feedback” screen:

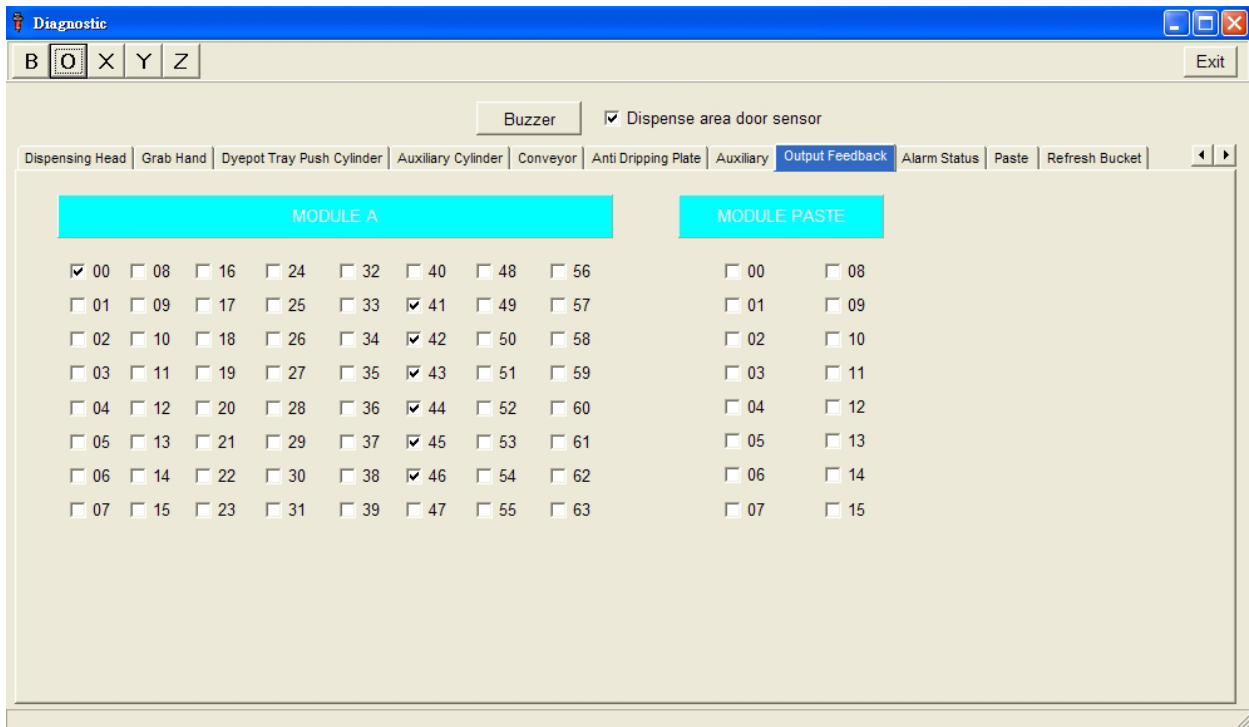

It is just the information about the current status of I/O.

## TF/TFCCtrl program – Diagnostic – Alarm Status


The picture below is the “Alarm Status” screen:

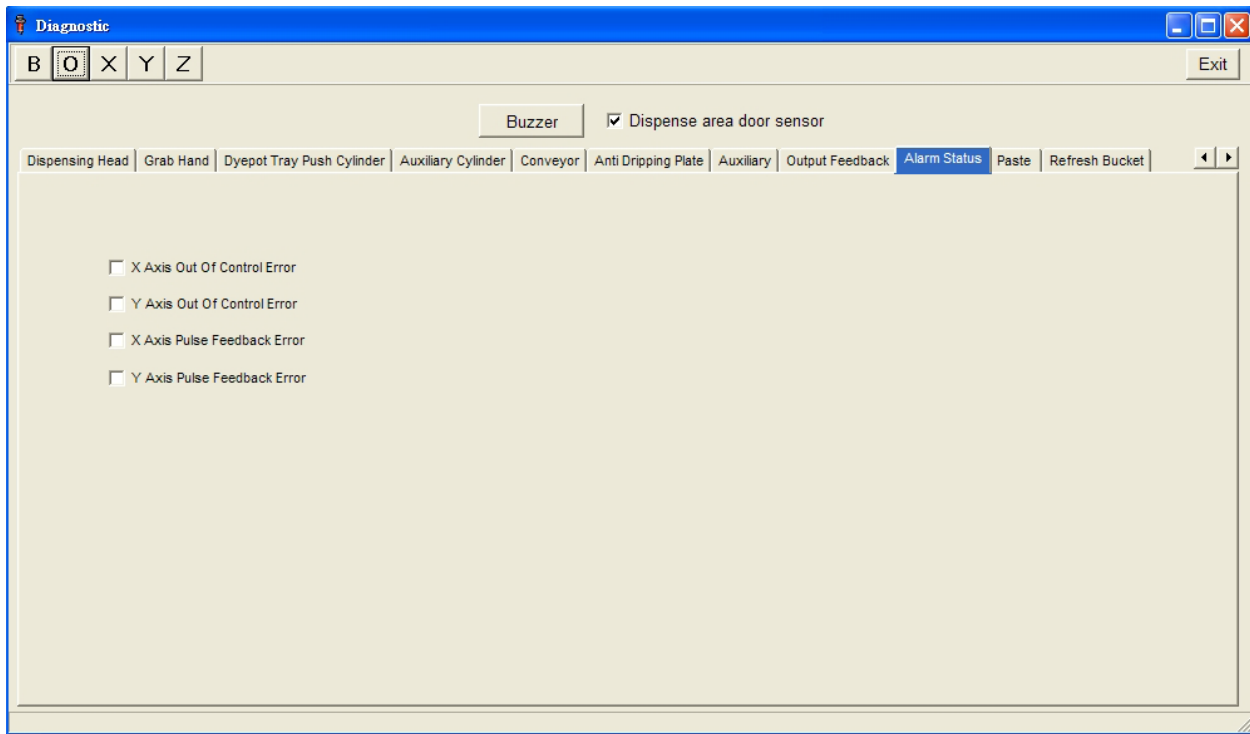

This function shows abnormal movement of the robot arm. When there is abnormal movement, the check box will be marked and the buzzer will alarm also. You have to reset the robot back to home positions. It is designed for security reason.

## TF/TFCCtrl program – Diagnostic – Ack button


The picture below is the “Ack Button” screen:

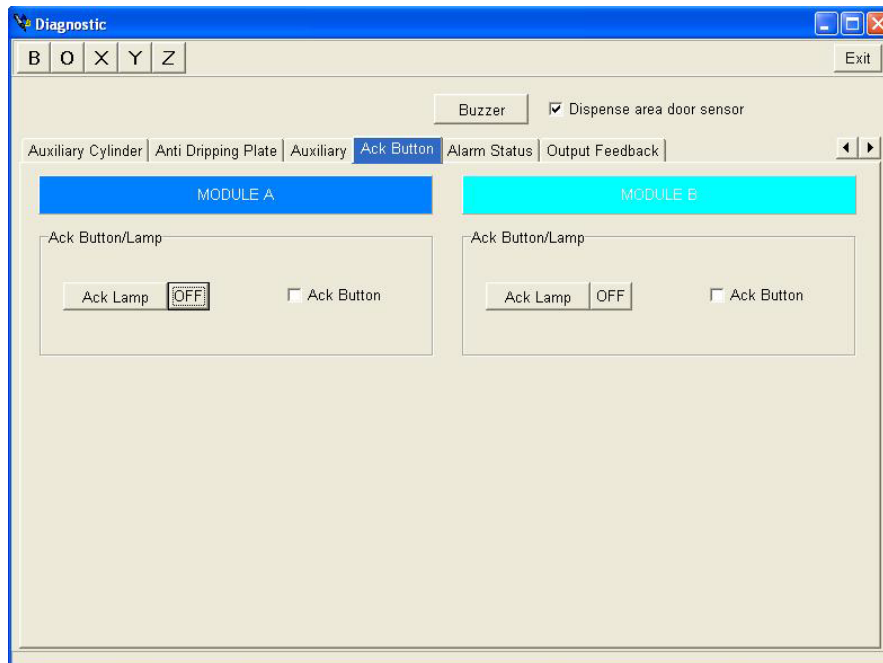

This is just for TF 2 scales systems. For checking if the “Confirm” buttons work fine or not.

## TF/TFCCtrl program – Diagnostic – Anti Dripping Plate

The picture below is the “Anti Dripping Plate” screen:

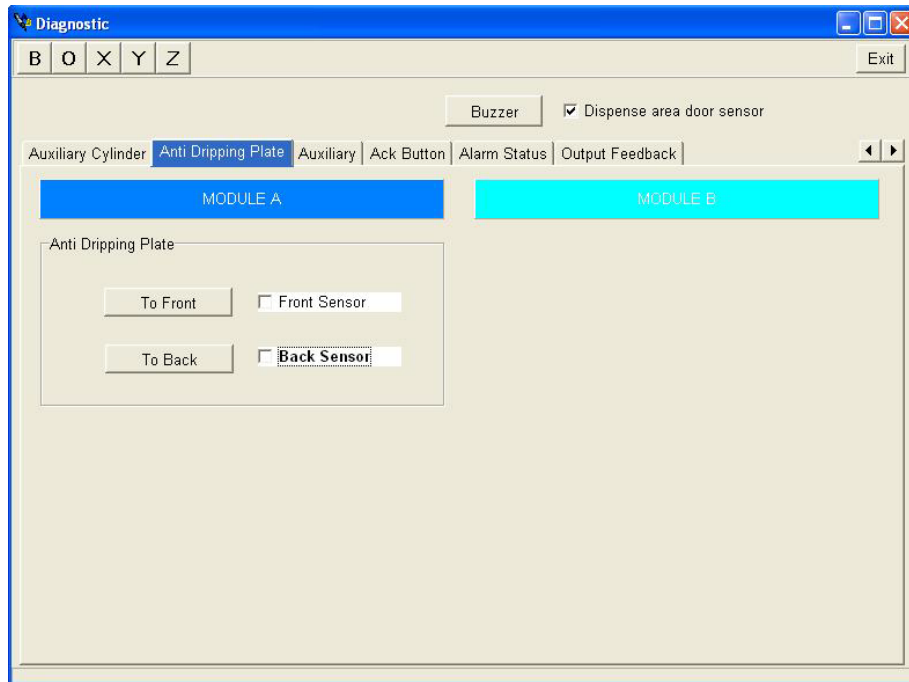


**To Front:** Click on the button, the anti-dripping plate will be moved outwards, and the check box of front sensor will be marked.

**To Back:** Click on the button, the anti-dripping plate will be moved backwards, and the check box of back sensor will be marked.

## TF/TFCCtrl program – Diagnostic – Auxiliary Cylinder

The picture below is the “Auxiliary Cylinder” screen:

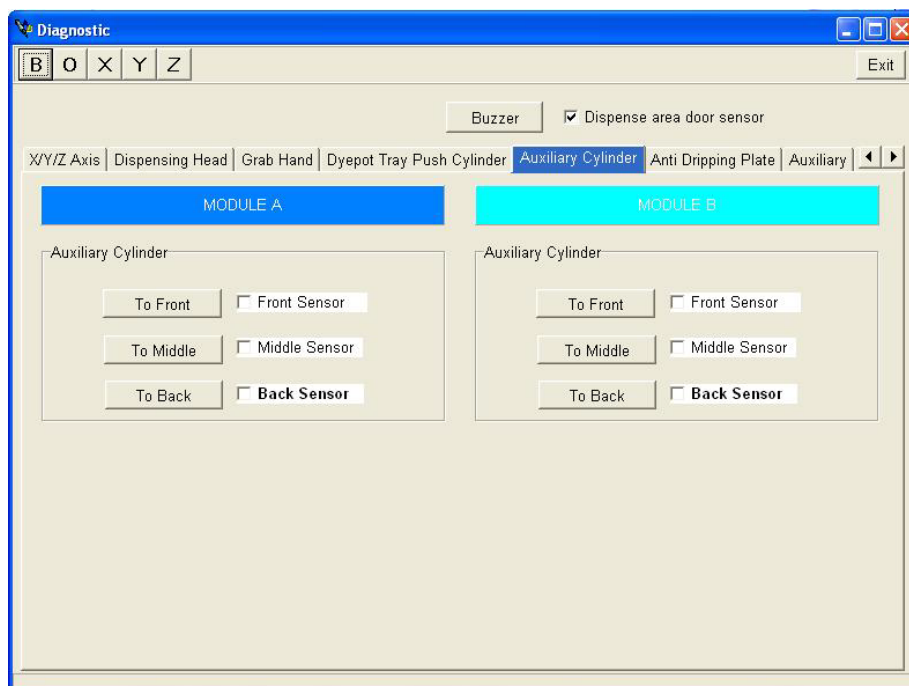

- To Front:** Click on the button, the auxiliary dispensing head will be moved to front position, and the check box of front sensor will be marked.
- To Middle:** Click on the button, the auxiliary dispensing head will be moved to middle position, and the check box of middle sensor will be marked.
- To Back:** Click on the button, the auxiliary dispensing head will be moved back, and the check box of back sensor will be marked.


**MODULE A:** The scale A (It is just for 2 scales system),

**MODULE B:** The scale B (It is just for 2 scales system).



## TF/TFCCtrl program – Diagnostic – Dyepot tray push cylinder

The picture below is the “Dye Pot Tray Push Cylinder” screen:

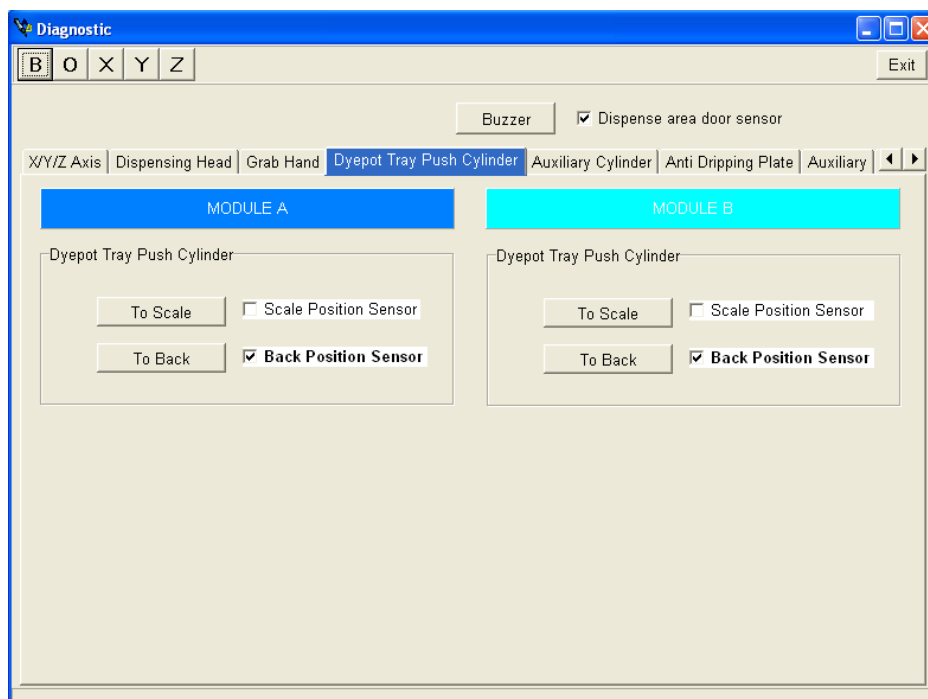

**To Scale:** Click on the button, the dye pot tray will be moved out onto the scale, and the check box of Scale Position Sensor will be marked.


**To Back:** Click on the button, the dye pot tray will be moved back, and the check box of Back Position Sensor will be marked.

**MODULE A:** The scale A (It is just for 2 scales system),

**MODULE B:** The scale B (It is just for 2 scales system).

## TF/TFCCtrl program – Diagnostic – Grab Hand


The picture below is the “Grab Hand” screen:

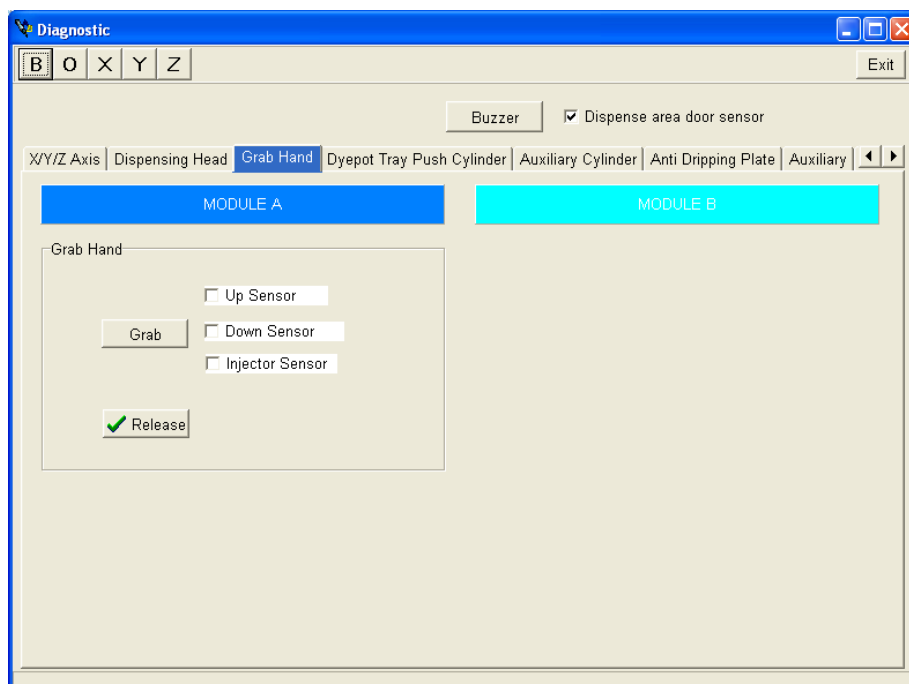

- Grab:** Click on the button to close the grab hand, and a “√” will appear on the button.
- Release:** Click on the button to release the grab hand, and a “√” will appear on the button.
- Up Sensor:** If the grab hand is on up position, the check box of up sensor will be marked.
- Down Sensor:** If the grab hand is on down position, the check box of down sensor will be marked.
- Injector Sensor:** In presentation of injector inside the grab hand, the check box of injector sensor will be marked.

## TF/TFCCtrl program – Diagnostic – Dispensing Head

The picture below is the “Dispensing Head” screen:

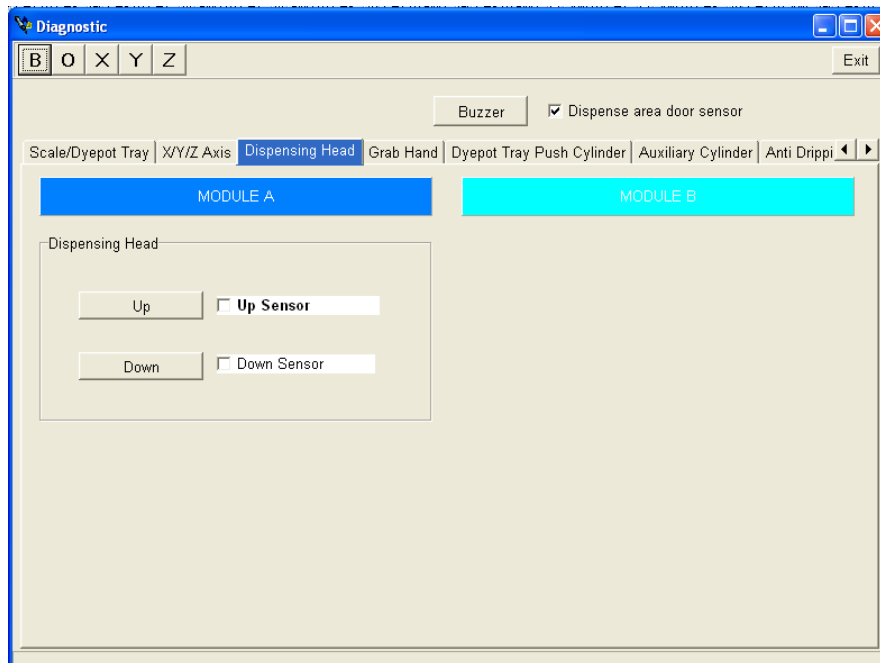


**Up:** Click on the button, the dispensing head will be moved up, and the check box of Up Sensor will be marked.

**Down:** Click on the button, the dispensing head will be moved down, and the check box of Down Sensor will be marked.

## TF/TFCCtrl program – Diagnostic – X/Y/Z axis

The picture below is the “x, y and z” screen:

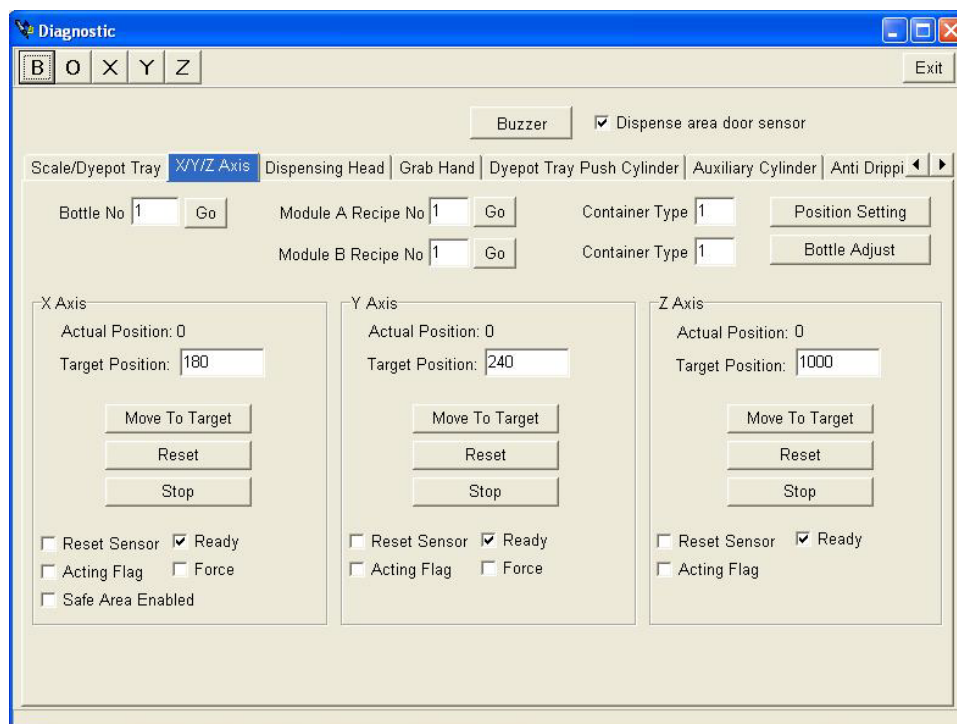

This function is for testing and calibrating the x, y and z coordinates of all the solution bottles, dye pot trays and movement of the XYZ motors. The coordinates are all calibrated for each machine prior to the system shipping. The x-y-z calibration data for each machine are shipped with each system (on the software CD )

### 1. Bottle No. :

Key in the bottle No. then click the button “Go”. The robot arm will move to the position of the bottle.

### 2. Module A Recipe No. :

Key in the dye pot number of the dye pot then click the button “Go”. The robot arm will move to the position of the pot on A scale.

### 3. Module B Recipe No. :

Key in the dye pot number of the dye pot then click the button “Go”. The robot arm will move to the position of the pot on B scale.

### 4. Container Type:


For set the type of dye pot tray.

## 5. Position Setting (Engineer password login only) :

This is engineer setting, for checking/setting the x, y and z axis of solution bottles and dye pot trays.


### 5.1. Whole Section :

#### 5.1.1. Section (1 to 4)

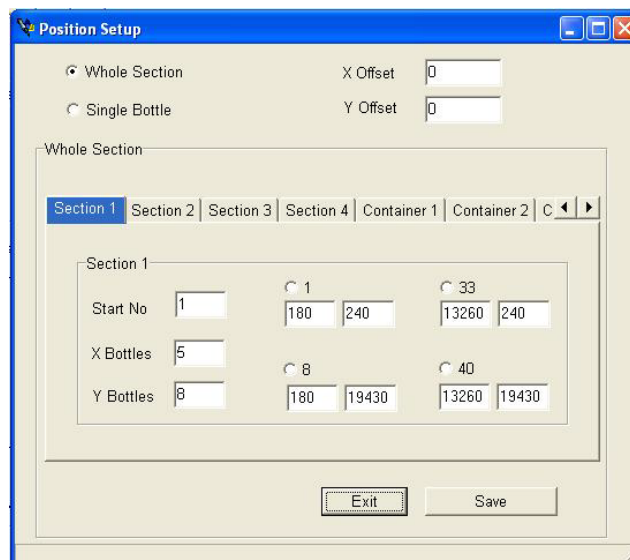


- **[Section\_1]:** coordinate of 4 corners of the 1<sup>st</sup> bottle tray (bottles 1 to 40).
- **[Section\_2]:** coordinate of 4 corners of the 2<sup>nd</sup> bottle tray (bottles 41 to 80).
- **[Section\_3]:** coordinate of 4 corners of the 3<sup>rd</sup> bottle tray (bottles 81 to 120).
- **[Section\_4]:** coordinate of 4 corners of the 4<sup>th</sup> bottle tray (bottles 121 to 160).

### 5.1.2. Container (1 to 4)

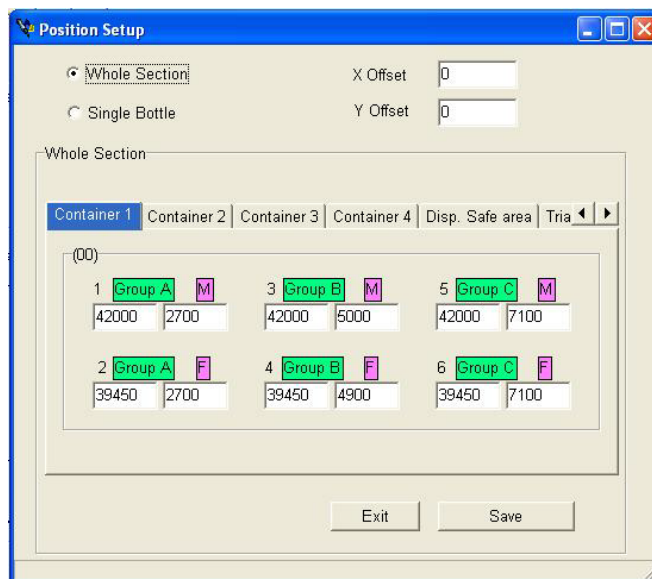


- **[Container 1]:** The X, Y position of 1<sup>st</sup> dye pot adapter.
- **[Container 2]:** The X, Y position of 2<sup>nd</sup> dye pot adapter.
- **[Container 3]:** The X, Y position of 3<sup>rd</sup> dye pot adapter.
- **[Container 4]:** The X, Y position of 4<sup>th</sup> dye pot adapter.

### 5.1.3. Disp. Safe area

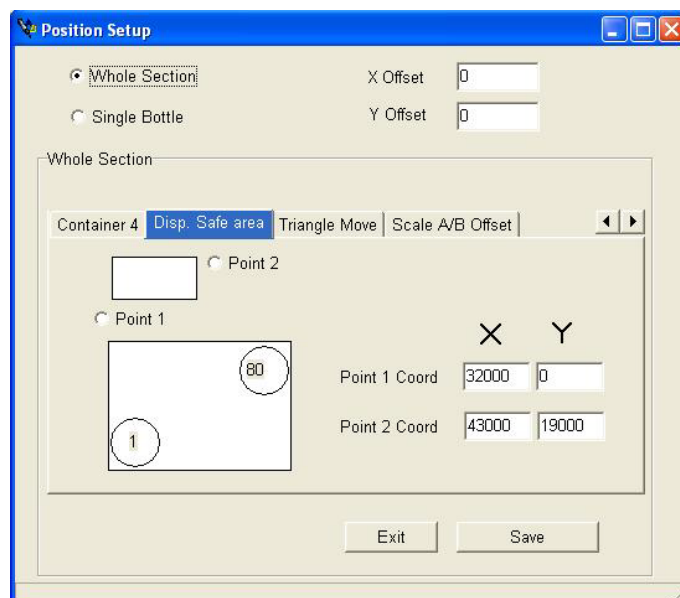


For set the area robot arm can only moving in that range. This is for safety issue. The default setting will be included in the calibration data in the software CD, you should not change it.

#### 5.1.4. Triangle Move

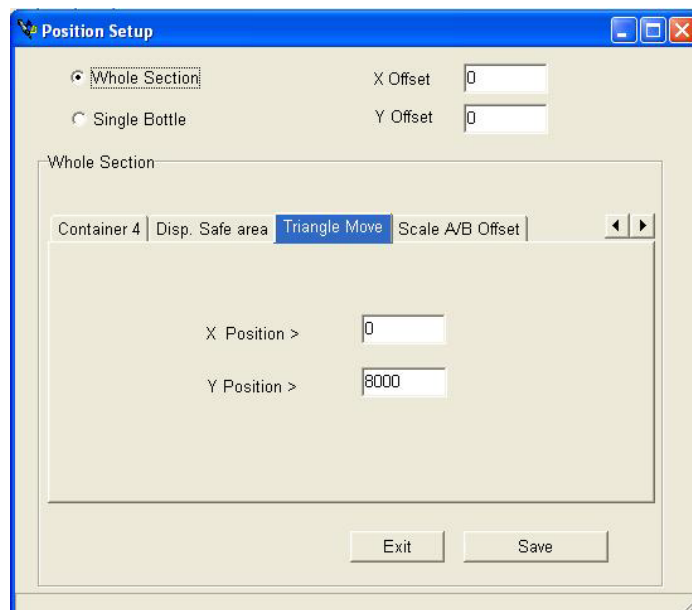

For set the area robot arm moving curve. The default setting will be included in the calibration data in the software CD, you should not change it

#### 5.1.5. Scale A/B Offset

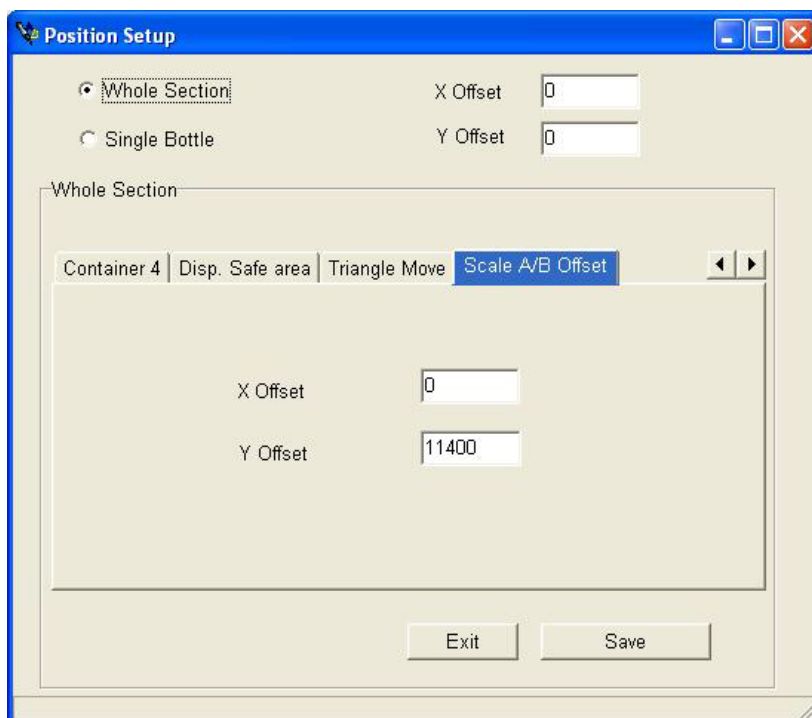

For set the offset between scale A and B. The default setting will be included in the calibration data in the software CD, you should not change it

## 5.2. Single Bottle :

The **Position Setup** dialog box is shown with the **Single Bottle** option selected. The **Position Test** section includes checkboxes for **Head Down**, **Grab**, and **Retry**, along with a **Test** button. The **Single Bottle** section features a list of bottle ranges (1-8, 9-16, 17-24, 25-32, 33-40, 41-48, 49-56, 57-64, 65-72, 73-80) with **1 -- 8** selected. An **Auto Change Row** checkbox is also present. A table displays X and Y coordinates for rows 1 through 8, with a **No Change** checkbox for each row.

	X	Y	No Change
1	180	240	<input checked="" type="checkbox"/>
2	180	2960	<input checked="" type="checkbox"/>
3	150	5770	<input checked="" type="checkbox"/>
4	130	8460	<input type="checkbox"/>
5	120	11220	<input type="checkbox"/>
6	140	13960	<input type="checkbox"/>
7	130	16730	<input type="checkbox"/>
8	130	19440	<input type="checkbox"/>



## 5.2.1. Position Test

This function is for testing if all x, y coordinates of solution bottles are correct.

- Test: Start the testing.
- Head Down: Moving the dispensing up/down
- Grab: Close/open grab hand
- Retry: Continue the testing with ONE injector only.

## 5.2.2. Bottle numbers: Select the bottle range (row) to start the testing.


## 5.2.3. Auto Change Row: Change to another row automatically.

## 5.2.4. X, Y: For manually change the coordinates.

## 5.2.5. No Change: If the check box is marked, the x, y coordinates will not be changed. So, if you want to change the x, y coordinates, you have to unmarked the check box, change the coordinates, and remarked the check box.

## 6. Bottle Adjust: (Power user password login only)

This is for calibrating the x, y axis of solution bottles. Click on the “Bottle Adjust”, the screen will show as below:

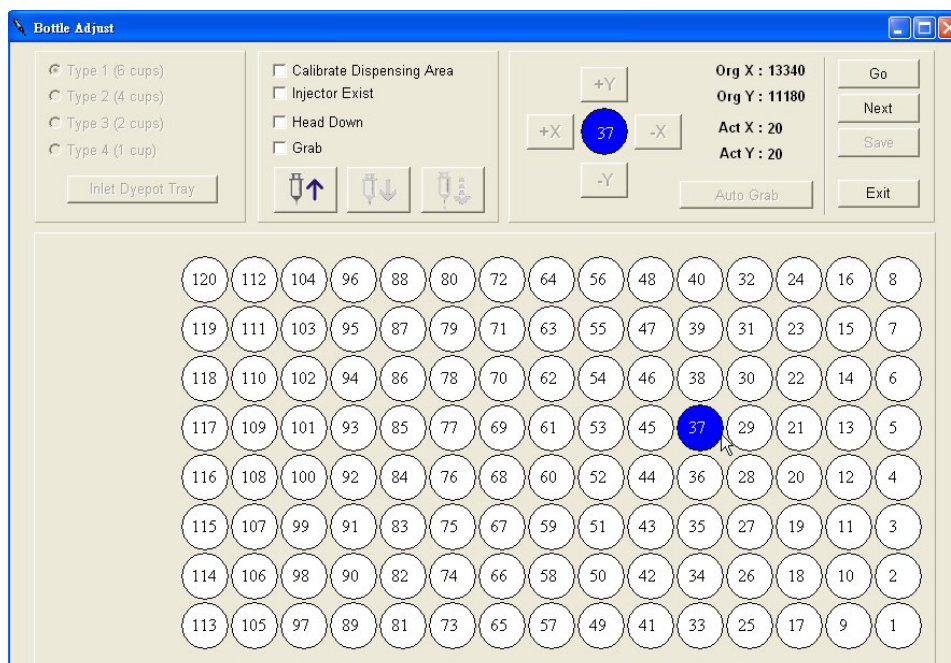


- 6.1. Calibrate Dispensing Area:** Shows the status of calibration, if engineer is calibrating
- 6.2. Injector Exist:** For sensing the injector, the check box will be marked if there is injector.
- 6.3. Head Down:** Enable the option, the dispensing head will move up/down in doing calibration.
- 6.4. Grab:** Enable the option; the grab hand will close/open in doing calibration.
- 6.5. +X, +Y, -X and -Y:** Click on the buttons to move the robot arm.
- 6.6. Act X:** Showing the actual position of X
- 6.7. Act Y:** Showing the actual position of Y
- 6.8. "Next":** Move to next position.
- 6.9. "Go":** Depends on the enabled options of "Head down" and "Grab", the system start the calibration.
- 6.10. "Exit":** Exit the function.
- 6.11. Bottle numbers:** For selecting bottle for doing calibration.

## 7. Dyepot Dispense Position Adjust: (engineer login only)


This is for calibrating the x, y axis of positions for dispensing solution into the dyepots in the dispensing area.

Enter “Bottle Adjust” screen and pick a bottle which you are going to take solution


from and the button “Get Injector”  will be enabled.

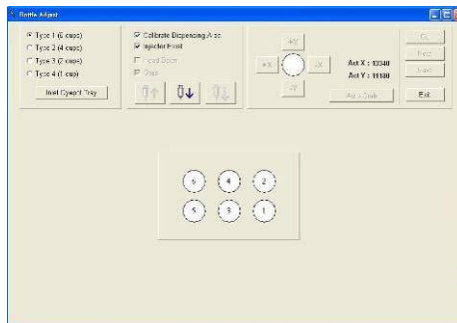



After clicking “Get Injector” button, the robot arm will physically go pick up the injector chosen and take 60cc solution. Meanwhile, the screen will automatically change to dyepot position calibration screen.

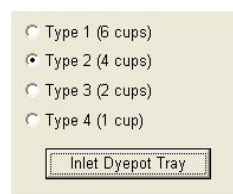


Choose the dyepot tray type you are going to calibrate positioning for, and click on the button “Inlet Dyepot Tray” in TFC system, or physically press the confirm button on the TF system, to feed dyepot tray into dispensing area.

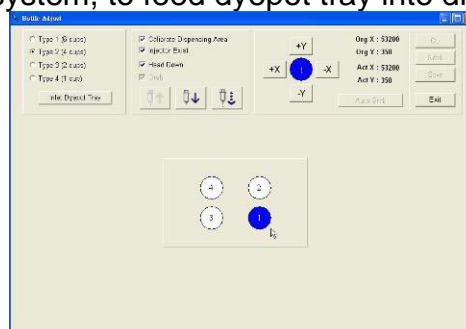

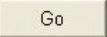
Select the dyepot position you are going to calibrate. Click button “Go”  and the robot arm holding the injector will go to the dyepot position chosen.

When the robot arrive the target position, the button “Dispense”  will immediately become enabled on a TF system. For TFC systems, you have to first move the dispensing head to the down position, then the “Dispense” button will become enabled.

Every time the “Dispense” button is pressed, the Z axis will move 1000 pulses to dispense a small quantity of solution. Observe the dispense position when it dispenses solution into the dyepot. Using the same way as calibrate solution bottle position, adjust the robot arm X/Y coordinate to make sure the position is in line with the middle of the dyepot.

After calibrating every dyepot position, click on the button “Put Injector” . The robot arm will then replace the injector back into the bottle.

## 8. X, Y and Z axis (Engineer password login only):

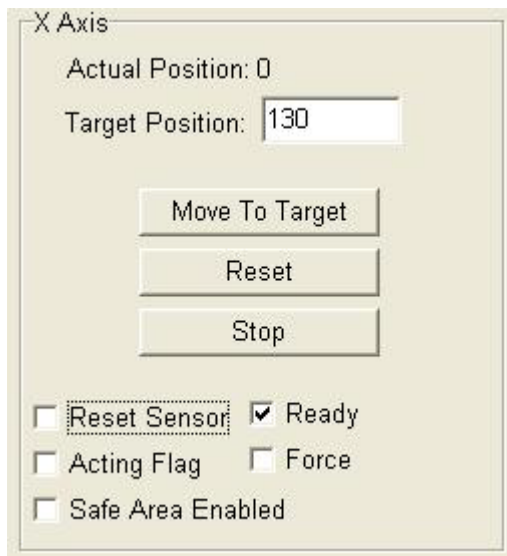


The screenshot shows a control panel titled "X Axis". It contains the following elements:

- Actual Position:** 0
- Target Position:** A text input field containing the value "130".
- Buttons:** Three buttons are stacked vertically: "Move To Target", "Reset", and "Stop".
- Checkboxes:**
  - ☐ Reset Sensor
  - ☒ Ready
  - ☐ Acting Flag
  - ☐ Force
  - ☐ Safe Area Enabled

- 8.1.1. "Actual Position": For showing the actual position (pulse).
- 8.1.2. "Target Position": Inputting target position (pulse), and click on the button "Move to target", the robot arm will move to the position.
- 8.1.3. "Reset": Reset the robot arm back to home position (pulse 0).
- 8.1.4. "Stop": Stop the movement of robot arm.
- 8.1.5. "Reset Sensor": The check box will be marked when the robot arm is at home position.
- 8.1.6. "Acting Flag" The safety doors are not closed properly.
- 8.1.7. "Safe Area Enabled": The safety area is enabled; the motion of robot arm is limited inside the safe area.
- 8.1.8. "Read": The system is ready.
- 8.1.9. "Force": Force the robot arm to move to the target position, and ignore all other safety issues.

## TF/TFCCtrl program – Diagnostic – Scale/Dyepot tray

The picture below is the “Scale/Dyepot Tray” screen:

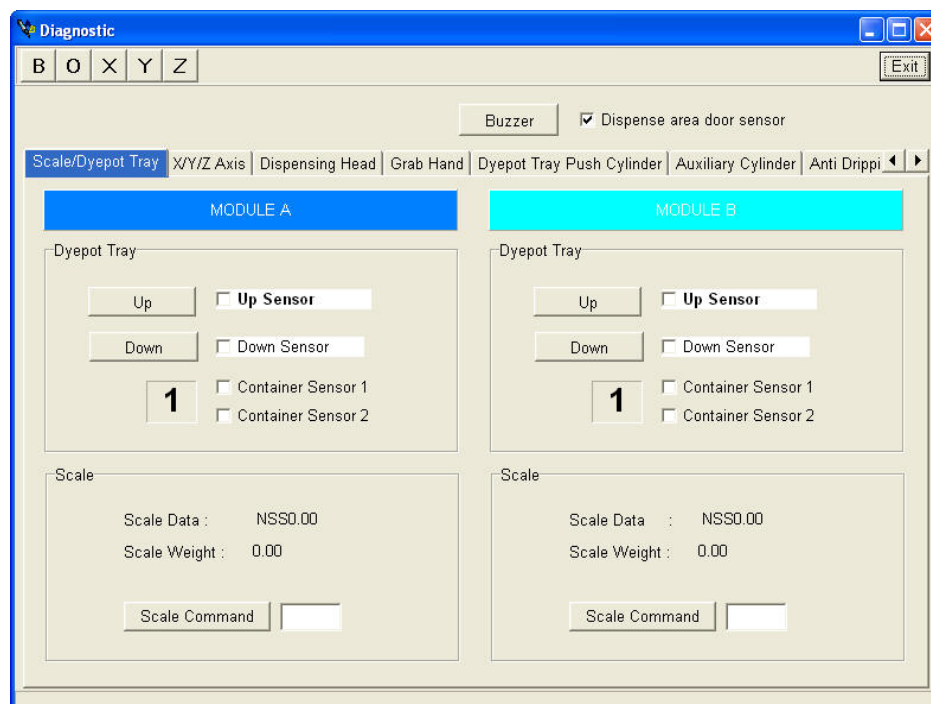

### 8.2. Dye pot Tray :


- 8.2.1. “UP”: Click the button to move the scale UP, and the check box for “Up sensor” will be marked.
- 8.2.2. “Down” Click the button to move the scale DOWN, and the check box for “Down sensor” will be marked.
- 8.2.3. “Container Sensor 1/2”: It is for testing the container type.

### 8.3. Scale :

- 8.3.1. “Scale Data”: Showing the reading on scale
- 8.3.2. “Scale Weight”:
- 8.3.3. “Scale Command”: For testing the scale. Inputting “sir” or “z” into the text box, and click the button, the scale will be reset.

### 3-3-2 TF/TFC Ctrl program – Specific wt. calibration

This is for calibrating the specific weight of dye solutions. Once all the stock solutions have been ready, engineer need to do the 1<sup>st</sup> specific weight calibration. After that, software will automatically do the calibration every time in dispensing dye solutions.

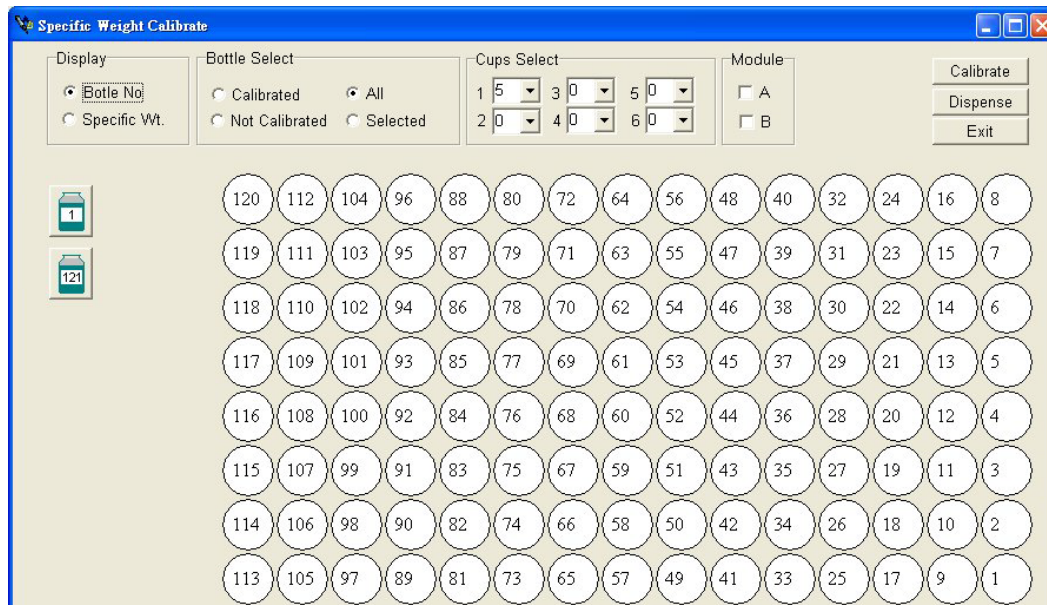

#### Display:

- Bottle No.: For showing bottle numbers.
- Specific Wt.: For showing the specific weight of every bottle.

#### Bottle Select:

- Calibrated: For showing calibrated bottles.
- All: Showing all calibrated bottles.
- Not Calibrated: For showing not calibrated bottles.
- Selected: Showing selected bottles only.

#### Cup Select:

For set the dispensing weight of on each dye pot.

#### Module:

Select module A or B for doing the calibration.

#### “Calibrate”:

Start the calibration.

#### “Dispense”:

#### “Exit”:

Exit the option

### 3-3-3 TF/TFCCtrl program – Cylinder Calibration


Entering “Cylinder Calibration” function, you will see the screen like this.

Sensor name	Standard time	Actual time	Result
<input checked="" type="checkbox"/> Dispensing head down	2000	1900	OK
<input checked="" type="checkbox"/> Dispensing head up	2000	1600	Too fast
<input checked="" type="checkbox"/> Grab hand release	800	900	OK
<input checked="" type="checkbox"/> Grab hand grab	800	900	OK
<input checked="" type="checkbox"/> Anti dripping plate front	700	50	Too fast
<input checked="" type="checkbox"/> Anti dripping plate back	1000	900	OK
<input checked="" type="checkbox"/> MA Dyept tray front	5000	4100	Too fast
<input checked="" type="checkbox"/> MA Dyept tray down	3000	4100	Too slow
<input checked="" type="checkbox"/> MA Dyept tray up	2000	2300	OK
<input checked="" type="checkbox"/> MA Dyept tray back	3000	2600	Too fast
<input checked="" type="checkbox"/> MA Aux cylinder front	3000	3100	OK
<input checked="" type="checkbox"/> MA Aux cylinder back	3000	3100	OK
<input type="checkbox"/> MB Dyept tray front	5000	-----	-----
<input type="checkbox"/> MB Dyept tray down	3000	-----	-----
<input type="checkbox"/> MB Dyept tray up	2000	-----	-----
<input type="checkbox"/> MB Dyept tray back	3000	-----	-----
<input type="checkbox"/> MB Aux cylinder front	3000	-----	-----
<input type="checkbox"/> MB Aux cylinder back	3000	-----	-----


1. Mark the checkbox to enable testing of this air cylinder.
2. On the m/c currently being tested there are no this kind of physical cylinders, so there is no checkbox.
3. Current cylinder being tested will be highlighted in blue.
4. “OK”: This means the cylinder speed is no problem.
5. “Too Slow”: It means the cylinder speed in that direction is too slow, please adjust it faster.
6. “Too Fast”: It means the cylinder speed in the direction is too fast, please adjust it slower.
7. “Select All”: Click on this button to select all air cylinders for testing.
8. “Repeat”: Mrk the checkbox to repeat the testing.
9. “Start/Stop”: To start/stop cylinder speed testing.
10. “Print”: Click this button to print out the test results



## How to adjust cylinder speed

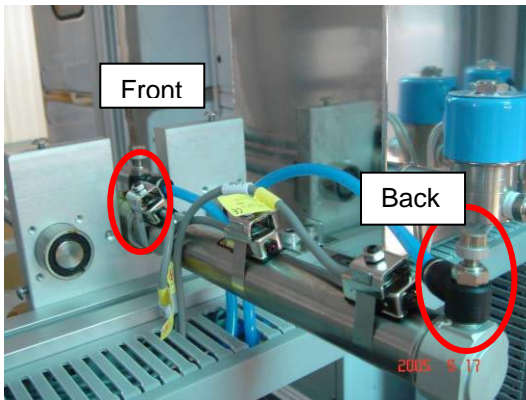


On most TF cylinders, you can find 2 speed adjustable quick-fit connectors. Turn the knob on the top of the connectors counter-clockwise to turn up the air volume so the cylinder speed will be faster. To change the cylinder speed to the front, adjust the knob on the front position. On the contrary, to

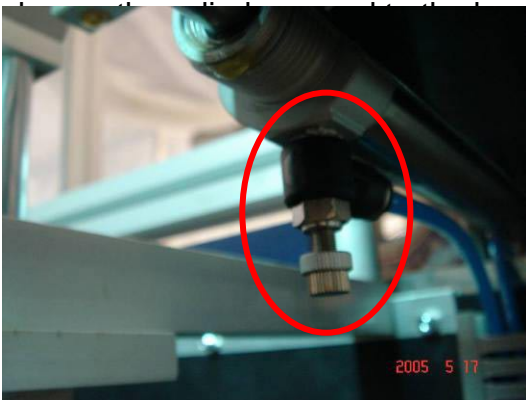

There is one special cylinder in TF system, the dyepot tray in/out cylinder. Due to the physical position of the cylinder, the front speed adjustable quick-fit connectors are installed on the solenoid valves in back cabinet so the users can easily adjust the speed. The speed adjust knob to back direction is the same as other cylinder which is installed on the cylinder body.

## Things you need to aware when using “Cylinder calibration” function

1. When checking dispensing head up/down speed, make sure you take the injector out from underneath the dispensing head to avoid damage to the injector or grab arm.

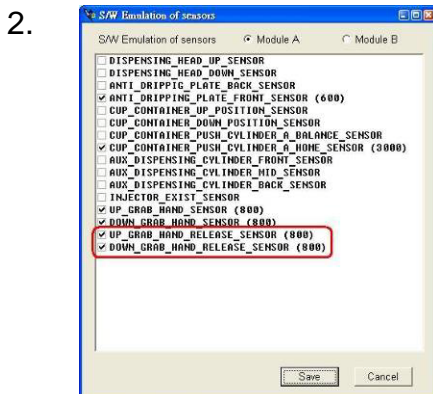

In software emulation function, the up grab hand and down grab hands are separated, so users can test each cylinders



speed.

3. If you find the cylinder speed varies, check the main air pressure. Make sure it is stable, then adjust the individual cylinder speeds again.
4. If you find the cylinder speed is getting slower, it might be caused by dirty stuff in the cylinder. Please check the air quality and water collect cup of main air regulator. If the air is not so clean, improve air quality. If the water is too full in the water collect cup, get rid of the water. Used correctly in the correct conditions, the air cylinders will have a longer life time.

### 3-4 TF/TFC Ctrl program – Parameter

#### 3-4-1 TF/TFC Ctrl program – Parameter – Setup

Input the password, and the setup screen will be show as below:

#### Basic Setting:

- “Batch Input Dir”: The directory for storing pre-batch data. Please leave the path settings as defined
- “Batch Output Dir”: The directory for storing completed batch (dispensed batch) data. Please leave the path settings as defined
- “Information Dir”: The directory for exchanging the machine status between DP & Control program.
- “Scale Type”: For selecting scale type. Currently, three types of scales can be used with TFC software, Mettler, Precisa and Sartorius.
- “User Password”
- “Power user password”: Power user is authorized to control all normal user functions and additionally to also use the “Bottle position adjust” function and “Cylinder speed calibration” function. The default password for power user is “DC”. This

password will only appear and only changeable under Power User or Engineer log in.


**MISC:**

- Handle First Component Error of Recipe
  - Enabled this option, software will stop the dispensing if there is problem with first recipe.
- Stop Dispense Rest Component Error.
  - Enabled this option, software will stop the dispensing if there is problem in dispensing.
- Wait on Batch Failure
  - Enable the option, if there is any error in dispensing, the Alarm will make sound last for time interval set to re-dispense the recipe.
- X axis low speed.
  - Enable the option, the robot will move in slower speed.
- Dispense paste then water for saving time (TF-P system only)
  - Enable the option, the system will dispense water immediately after completing paste dispensing. Default setting is function enabled.

**V5 Injector setup**

Once you click on this you will get option to setup V5 injectors. You need to select from which positions you would like to use these injectors and system will displayed on graph as shown below :

Example V5 injectors put from 1 to 10 :

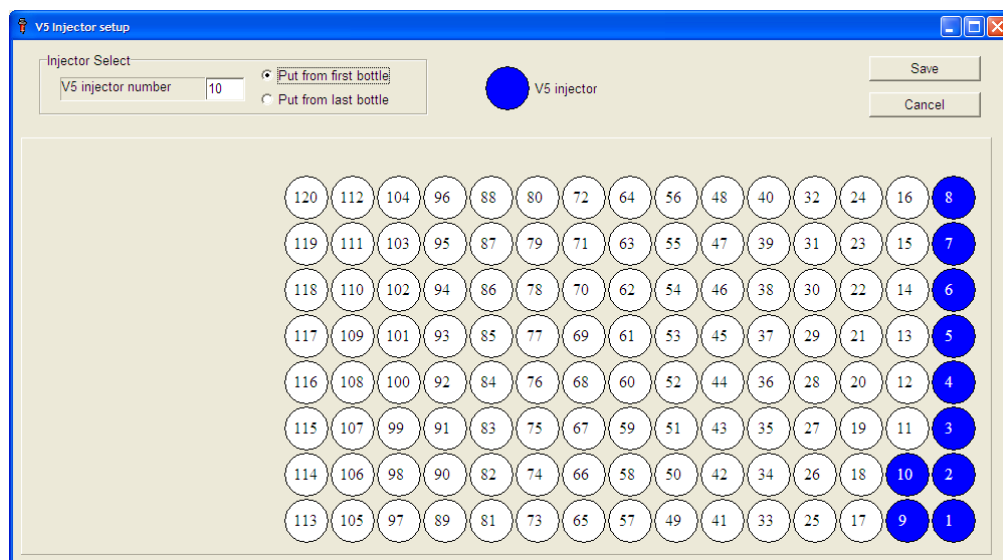

## TF/TFC G2 Operation Manual

If you put them from last it would show as follows:

V5 Injector setup

Injector Select  
V5 injector number: 10

☐ Put from first bottle  
☒ Put from last bottle

V5 injector

Save  
Cancel

120	112	104	96	88	80	72	64	56	48	40	32	24	16	8
119	111	103	95	87	79	71	63	55	47	39	31	23	15	7
118	110	102	94	86	78	70	62	54	46	38	30	22	14	6
117	109	101	93	85	77	69	61	53	45	37	29	21	13	5
116	108	100	92	84	76	68	60	52	44	36	28	20	12	4
115	107	99	91	83	75	67	59	51	43	35	27	19	11	3
114	106	98	90	82	74	66	58	50	42	34	26	18	10	2
113	105	97	89	81	73	65	57	49	41	33	25	17	9	1

**TF/TFC Ctrl program – Parameter – Setup – Engineering Setup**

Click on the “Engineer” button, the engineer setup will show. This is only available with engineer login.

The screenshot shows the 'Engineer Setup' window with the following sections and values:

Parameter	Value
Number of Injector Bottles	120
Grab Injector Z Position	60000
Leave injector bottom speed	1000
Leave injector bottom pulse	800
Middle Conveyor Take Time	10000
Balance stable time (ms)	1200
Z axis cw/ccw gap pulse	100
Valve minimum open time (ms)	20
Unable learning assume spec. wt.	1.000
Robot auto reset period (hour)	12

**Paste Setup**

☐ Including paste device

**RS232 Setting**

COM Port	1
Baudrate	57600
Timeout (ms)	300

**Mettler Scale Weighing Process Adaptor Setting**

☐ Medium Drop ☒ Fine Drop

**Scale Setting**

☒ Define Tare Command

**Other Scale Type Setting**

☐ Scale Cmd Pre Code

Scale Str Ignore Bytes

Scale Start Command

**Weighing Settings**

☐ Ignore unstable 0.01g floating scale

☒ Enable Multi Shots Function

**Injector Dispensing Speed**

☒ Normal ☐ Slower ☐ Slowest

☒ Demo Mode Available

Normally, for running the software at customer site, all the settings should only modified by engineers or under engineer's direction.


**Unable learning assume specific weight**

The control s/w will detect the solution specific weight if the solution's specific weight not yet acquired, but to detect solution specific weight need at least 4 grams target weight for dispensing. So if the target weight less than 4 grams the s/w will assume the solution's specific weight as defined until once the solution's target dispensing weight large than 4 grams then the s/w will automatically detect the specific weight of the solution.

While using V5 injectors for high concentration solutions you should set the value which is maximum Specific gravity obtained using calibration menu. Calibrate few of the bottles (no need to calibrate each bottle) having high concentration solution and then checked the specific weight in this menu and whatever maximum value you get among these bottles would be value for this parameter.

## TF G2 Operation Manual

The example is as below:

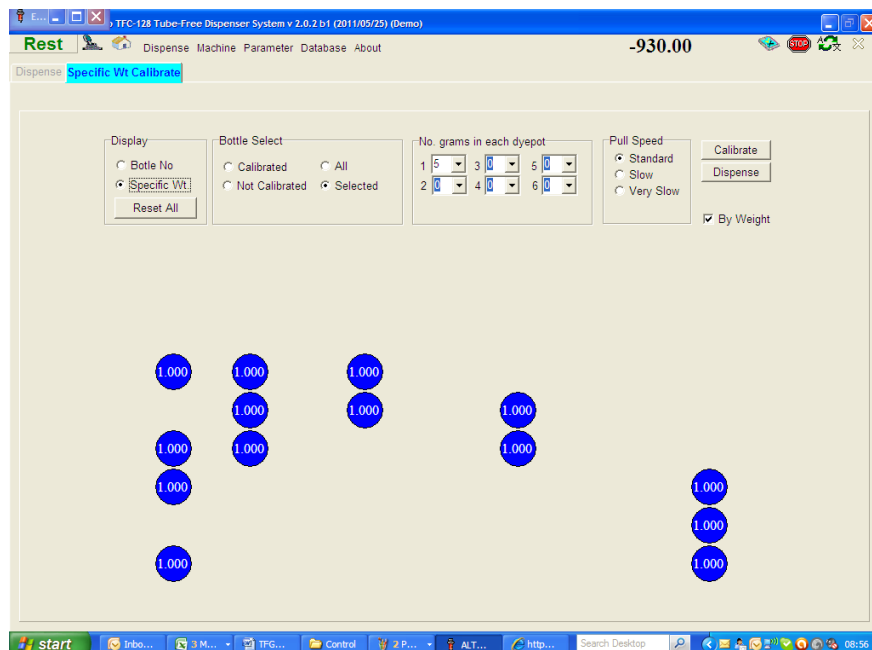

As all values for Specific weight is 1, so our maximum would be 1.000

### “RS232 Setting”:

- **COM Port:** It is the PC COM Port the system is connected.
- **Baud rate:** It must be 57600
- **Timeout (ms):** it must be 200

### “Use drip tray when moving between dyepots”:

Enable this function, in dispensing, the drip tray will be moved out in moving to different dyepots. This function is for avoiding any pollution from occurred when dispensing head move from dyepot to dyepot but will make dispensing speed slow down.

### “Ignore unstable 0.01g floating scale reading”:

Enable this function to eliminate the affection of scale by slightly vibration floor. If the vibration is severe, suggest put professional anti-vibration material under scale platform feet.

### “Enable multi-shots function”:

When the function multi-shots is enabled, the system will keep dispensing a solution until it achieves the target weight required, up to a maximum of 10 shots. Typically it takes 2 shots to reach target weight, but if this is not achieved in the first 2 shots, it will continue to dispense. When disabled, the system will stop dispensing after the first 2 shots, and move onto the next solution, even if the target weight has not been reached. It is recommended that Multi shots is always enabled, unless accuracy of dispense is not required and fast throughput is more important, in this case multi shot can be disabled.

### “Demo Mode Available”:

If you want to run the program in demo mode (without system connected), tick the option “Demo Mode Available”, and restart the program.

**“Mettler scale weighing process adaptor setting”:**

Set up Mettler scale weighing process to middle drop or fine drop during dispensing. It is recommended to always use fine drop for dispensing. Middle drop should only be used when specifically instructed by Datacolor.

**“Paste Setup”:**


If you choose to install TF-P model in the installation program, the parameters “Including paste device” will be automatically enabled.

**“Injector Dispensing Speed”:**


This setting can change the speed of moving Z axis when dispensing solutions to reduce the reflection force of dispensed solution and ease the situation of possible splashing. There are 3 options, “Normal”, “Slower” and “Slowest”. “Normal” is the standard speed which is used in all previous versions. The splashing problem is various with the diameter and height of dye pot. Shorter and larger diameter dye pot have more serious splashing problem. Slow dispensing speed can reduce the possibility of splashing, but it takes more time to finish the dispensing of the recipe. If splashing is occurring it is recommended to try the “slower” setting first before switching to “slowest”.

**S/W Emulation of sensors**

This is for setting the time of software emulating sensors.

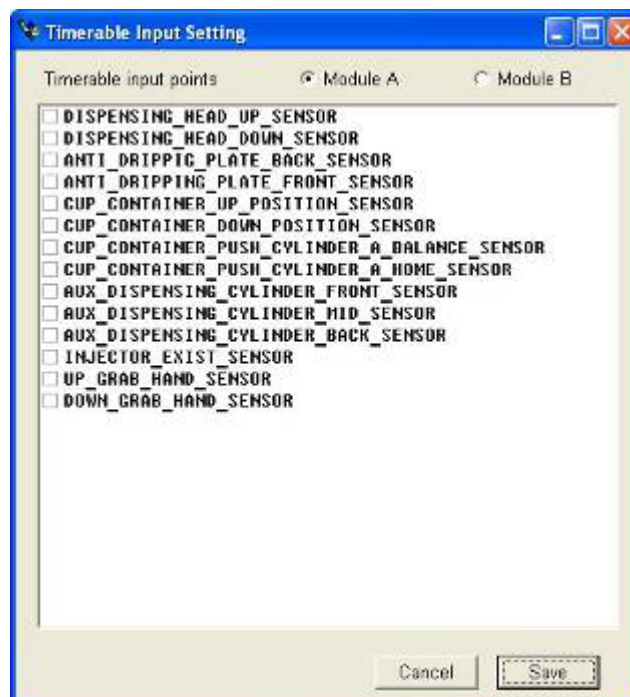

In case there is a faulty sensor as in the list below, to make the system temporally workable, you need to use the function of S/W emulation sensor. Click on the items, a dialogue will pop up with default time setting.

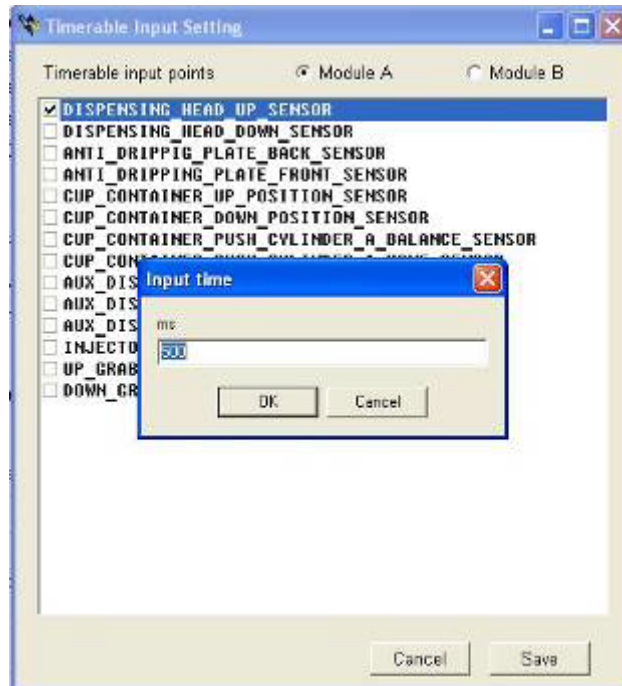

Click on “OK” to close the dialogue, and then click “Save” to save the setting.



### 3-4-2 TF/TFCtrl program – Parameter – Aux Setup

---

#### Emergency Stop :




: Click this button for stopping all the action of TFC.

#### Exit :



: Exit, close the “Tool” function.

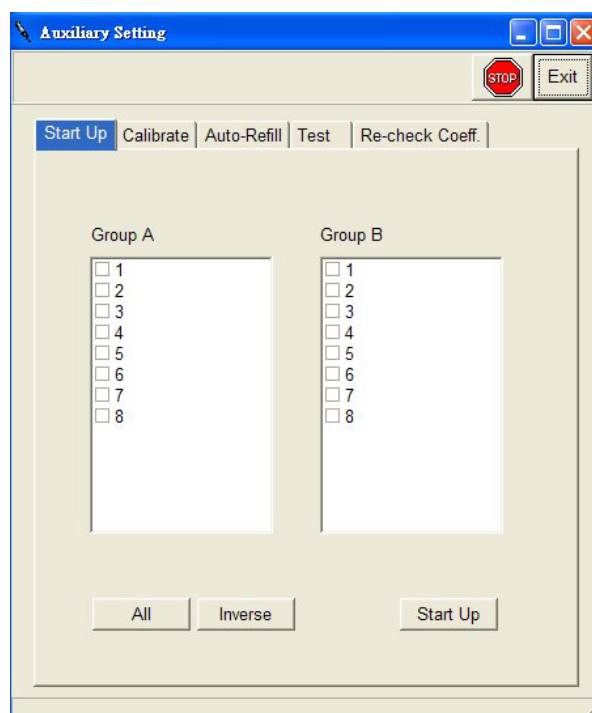

#### Start Up:

Select the valve needed to be started up and click “Start Up” button. The system will open the selected valve for starting up.

In TFC Ctrl program, the calibration tray will be automatically loaded in chemical valve “Start up” function. Users need only click “Start” button in the “Start Up” screen and the tray will be auto loaded.


After the tray is loaded and read by scale, the valves which the user has chosen to start-up will open in sequence. If the scale reads the dispensed chemical as being over 15g, or the valve dispense flow rate is over 0.5g/sec, the status of the valve will show “OK”. The completed valve will then close automatically after start up, and the system will then move onto the next valve selected.

If the flow rate of the valve is under 0.5/sec, the status of the valve will show as “Low”.

If the dispensed weight cannot achieve 15g in one minute, the status of the valve will show “Time out”.

If there is any “Low flow rate” and “Time out” status when all chosen valves are completed start up, the system will pop up with a dialogue to ask users if they want to redo the start up of valves which have failed.

## Calibrate:

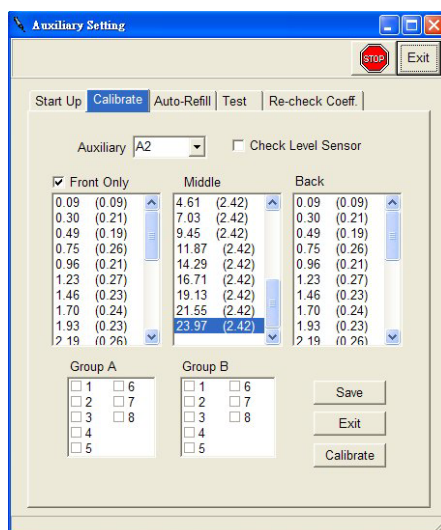

This is for calibrating the auxiliaries and water. For fast and accuracy dispensing of water and auxiliaries, suggest doing the calibration once a week.

Select A1, B1 or C1 in Module A or B for showing the calibrated result:

Module A	Module B
A1	B1
0.53 (0.53)	0.35 (0.35)
1.29 (0.76)	0.99 (0.64)
2.03 (0.74)	1.67 (0.68)
2.85 (0.82)	2.43 (0.76)
3.72 (0.87)	3.19 (0.76)
4.52 (0.80)	3.98 (0.79)
5.36 (0.84)	4.75 (0.77)
6.17 (0.81)	5.49 (0.74)
6.97 (0.80)	6.24 (0.75)
7.81 (0.84)	7.04 (0.80)
=====	=====
7.81 (7.81)	7.04 (7.04)
16.07 (8.26)	14.86 (7.82)
24.32 (8.26)	22.68 (7.82)
32.58 (8.26)	30.50 (7.82)
40.84 (8.26)	38.32 (7.82)
49.09 (8.26)	46.14 (7.82)
57.35 (8.26)	53.96 (7.82)
65.61 (8.26)	61.78 (7.82)
73.86 (8.26)	69.60 (7.82)
82.12 (8.26)	77.42 (7.82)

**“Check Level Sensor”:** Check the float sensor for refilling auxiliaries and water.

**“Module A/B”:** Selecting dye pots for doing calibration.

**“Cancel”:** Click the button to close the current screen.

**“Save”:** Click the button to save the settings.

**“Calibrate”:** Click the button to start the calibration.

## Test:

For test the accuracy of dosing auxiliaries.

The screenshot shows the 'Auxiliary Setting' window with the 'Test' tab selected. The window has a title bar with standard Windows controls. Below the title bar are 'Stop' and 'Exit' buttons. A tabbed interface at the top includes 'Start Up', 'Calibrate', 'Auto-Refill', 'Test' (active), and 'Re-check Coeff.'. Below the tabs, there are two dropdown menus: 'No. Recipes' set to '1' and 'Dyepot Type' set to '1'. The main area contains a table for valve settings:

Valve No	Target Weight(g)
<input type="checkbox"/> 1	5
<input type="checkbox"/> 2	5
<input type="checkbox"/> 3	5
<input type="checkbox"/> 4	5
<input type="checkbox"/> 5	5
<input type="checkbox"/> 6	5
<input type="checkbox"/> 7	5
<input type="checkbox"/> 8	5

To the right of the table, there is a checkbox for 'Gravimetric' and a text field for 'Tolerance(g)' set to '0.03'. A 'Dispense' button is located at the bottom right of the table area.

**“Recipes”:** Select how many recipes to be tested

**“Cup Container Type”:** For selecting the dye pot tray type.

**“Valve No.”:** For selecting valve to test.

**“Target Weight”:** For setting the dispensing weight.

**“By Weight”:** For setting the dispensing weight.


**“Tolerance”:** For setting tolerance in dispensing.

**“Module”:** For setting which module for dispensing.

**“Dispense”:** Click the button to start the testing.

### Re-check Coefficient:

This function is similar with “Calibrate” but it is faster. After doing auto refresh, the system will get a factor of valves to make water dispensing more accurate. We suggest users do it everyday before working.

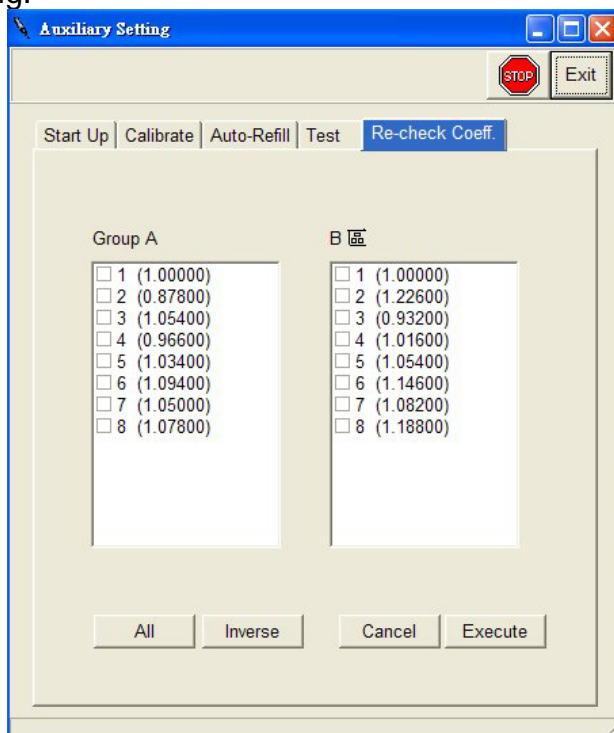

“Test 5/10 g”: For setting the dispensing weight.

“A1, B1 and C1 of Module A and B”: For selecting which valve for doing refresh.

“All”: Selecting all valves.

“None”: De-select all valves

“Exit”: Exit the program.

“Refresh”: Start the refresh.


### 3-4-3 TF/TFC Ctrl program – Parameter – Paste Setup

---


There are 3 functions for paste, which are “Clean”, “Test” and “Calibration” for paste module.

#### **Clean function**


“Clean” function is for cleaning the paste dispense pipes for regular cleaning maintenance. The paste tank should be filled with water and the tank and pipes can then be cleaned using this option, by dispensing water through the pipes. Alternatively the paste pipes can be connected to the mains water, instead of filling the tank with water, and clean in that way.

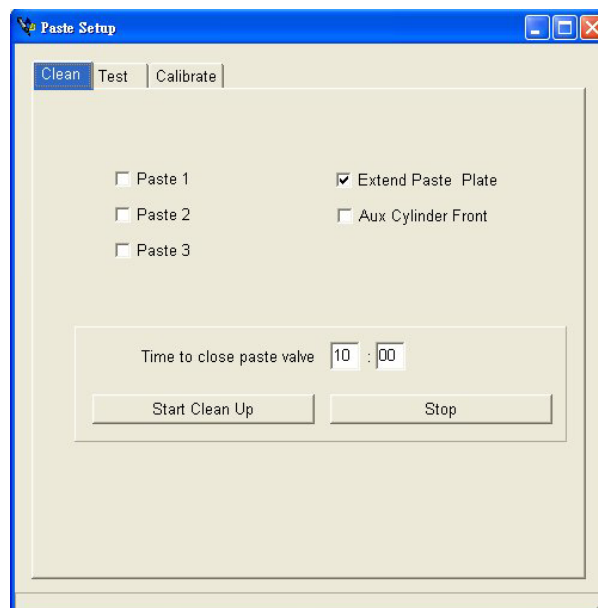

You can set the timer to automatically stop the cleaning paste automatically. The default setting is 10 minutes.

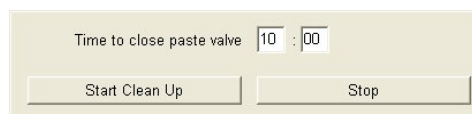


Click “Start clean up” button. The system will show a reconfirmation message to ensure cleaning tube has been fixed on the paste dispensing valve.

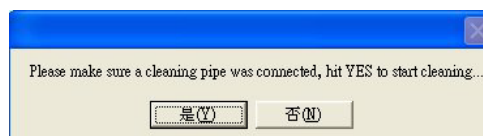

Make sure the paste cleaning tube has been fixed correctly. Click “Yes” and the paste dispensing valve will open to start cleaning paste pipe.


For more hardware detail, please refer to TF “Guidelines on correct operation of TFC dispensers to ensure best dyeing results” manual revision 7.0.

## **Test function**

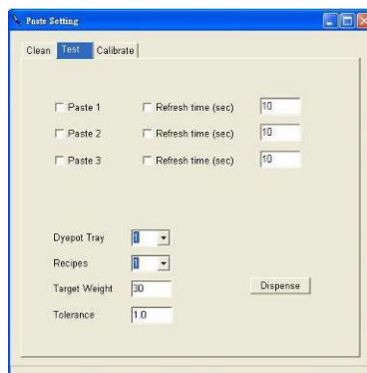

The “Test” function is for testing paste dispensing accuracy. The way to use this function is the same as chemical and water test function. Select dyepot tray type and recipe quantities then filling target weight and tolerance. Choose the paste you are going to test. Click “Dispense” button to start dispensing paste.

## **Calibrate function**

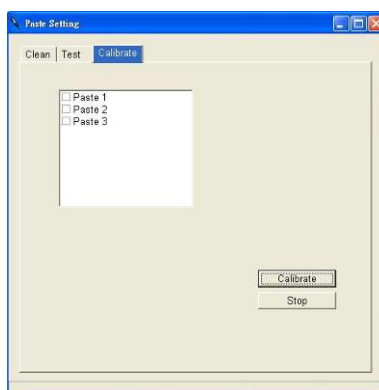

The “Calibrate” function is for calibrating paste dispensing coefficient. The way to use this function is the same as chemical and water calibrate function. Choose the paste you are going to calibrate then click “Calibrate” button to start calibrating.

It is recommended to carry out a calibration after changing/refilling paste.

### 3-4-4 TF/TFC Ctrl program – Parameter – Print Setup

---

This is for setting the printer parameters.

**Printer Option**

☒ A4 size printer  
☐ Label printer

☒ Weight ☐ Concentration

**Print Option**

☐ Print dispense reports    ☐ Template TF\_SixBlock.prt

☐ Print Recipe Detail    ☐ Template TF\_Detail.prt

☐ Customer Report Program

☐ Print target weight less than 4.00 g

☒ Print all dispensed batches  
☐ Print complete batches with recipe failure  
☐ Print only failed recipes

Save

#### Printer Option

There are 2 types of printer can be select, A4 or Label printer. When the Label printer is selected, the weight or concentration needs to select when print out due to the width limit of label printer.

**Prints dispense reports:** Enable the option; it will print out a six-recipe report in a page.

**Print Recipe Detail:** Enable the option; it will print out a recipe detail report, which each recipe takes one paper.

**Customer report program:** When the customer need a special report, it need a special report program to print out, a program name need to be define here.

**Print target weight less than X g:** it will automatically print a report after a recipe is finished if the target weigh of any product is less than the predefined target weight. Then the supervisor will be aware that the recipe was created wrongly.

**Options for selecting when to print the report:**

1. Print all dispensed batches
2. Print complete batches with recipe failure
3. Print only failed recipes



### 3-4-5 TF/TFC Ctrl program – Parameter – Agitation Setup

---

This is for setting the agitation parameters:

**Mixer Setting**

Section 1

☒ Enable

CW Time  Sec

STOP Time  Sec

CCW Time  Sec

Section 2

☒ Enable

CW Time  Sec

STOP Time  Sec

CCW Time  Sec

Section 3

☐ Enable

CW Time  Sec

STOP Time  Sec

CCW Time  Sec

Section 4

☐ Enable

CW Time  Sec

STOP Time  Sec

CCW Time  Sec

☒ PC Control Mixer

Cancel Save

**“Section 1, 2, 3 and 4”:** Maximum 4 sections agitation module in TFC systems, depends on the model of the system.

**TFC 40:** 1 section

**TFC 80:** 2 sections

**TFC 120:** 3 sections

**TFC 128:** 3 sections

**TFC 168:** 4 sections

**“CW Time”:** Time for clockwise agitation

**“STOP Time”:** Time for stopping the agitation when switching between CW and CCW.

**“CCW Time”:** Time for counter clockwise agitation

**“PC Control Mixer”:** Enable the function to control the agitation by PC.

**“Cancel”:** Close the program

**“Save”:** Save the settings.

### 3-5 Summary of features of TFC-P system

- Paste dispensing function available for new TFC-P systems.

#### TF-P Dispensing Progress Program (DP)

In DP program Database->Product page, the paste data must be define first.

Bottle no.	Product code	Solution conc.	Solution unit	Default process	Specific gravity	Strength	Lifetime (hr)	Warning level	Color
231	Paste1	5.00000	%	P1	1.00	100.00	72	200.00	

Define the paste bottle no in solution page, the bottle number of paste is from 231 to 233.

Product code	Product name	Product type	Exhaust	Min. solution co.	Max. solution co.	Price	Supplier	Color
Paste 2		Paste		100.00000	100.00000	0.000		
Paste1		Paste		1.00000	10.00000	0.000		
Water	Water	Chemical	%	1.00000	10.00000	0.000		

**Dispensing Control Program (Ctrl)**

The screenshot shows the 'Engineer Setup' window with the following settings:

Number of Modules	1
Number of Injector Bottles	40
Grab Injector Z Position	60000
Leave Injector Bottom Speed	1000
Leave Injector Bottom Pulse	800
Scale Stable Time (ms)	1200
Z axis cw/ccw Gap Pulse	100
Valve Minimum Open Time (ms)	10
Unable learning assume spec. wt.	1.0
Robot auto reset period (hour)	12

**Paste Setup**

☒ including paste device

**RS232 Setting**

COM Port	1
Baudrate	57600
Timeout (ms)	300

**Mettler Scale Weighing Process Adaptor Setting**

☐ Medium Drop ☒ Fine Drop

**Injector Select**

V5 injector number: 0 ☒ Put from first bottle ☐ Put from last bottle

☒ Demo Mode Available

**Module A Scale**

☒ Define Tare Command: Z

**Other Scale Type Setting**

☐ Scale Cmd Pre Code: 0

Scale Str Ignore Bytes: 3

Scale Start Command: SIR

**Module B Scale**

☒ Define Tare Command: Z

**Other Scale Type Setting**

☐ Scale Cmd Pre Code: 0

Scale Str Ignore Bytes: 3

Scale Start Command: SIR

**Other Weighing Settings**

☐ Use drip tray when moving between dyepots

☒ Ignore unstable 0.01g floating scale reading

☒ Enable Multi Shots Function

**Injector Dispensing Speed**

☒ Normal ☐ Slower ☐ Slowest

Buttons: Save, S/W Emulation of sensors, Exit/Cancel

In “Engineer Setup” of Ctrl program, there is a new setting “Paste Setup” which includes the options “enable paste device”.

## PART 4: Autolab SPS Program

### 4-1 Autolab SPS Program – Overview

TFC SPS is for preparing stock solutions. The main screen is as below:

**In G2 software, all solution database and making solution process have to be created through AutolabDP software and link SPS software to the DP database.**

The SPS has 2 types of user interface: menu and the Visual Display area.

#### Menu bar:

**Auto**

: **Auto Mode**

It means that the system is computer controlled.


**12.00**

: **Displays the weight on the scale**

It is just beside the “Auto” Mode, for display the reading of scale.

: **Mixing Areas**

There are 12 mixing bottles data in the mixing area. Each mixing bottle data includes machine number, bottle number, and remaining time of mixing.

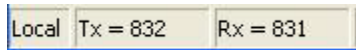

**: Status bar**

On the bottom of main screen displays information such as current status of the machine, software edition, and so on.



**: Emergency Stop**

Press it to stop making solution immediately.



**: Change Languages**

Change languages as your request. The default language is English.

## 4-2 Autolab SPS Program – Dispense

### 4-2-1 Autolab SPS Program – Dispense – Making Solution

You can choose the option “Make solution” in Dispense function of menu bar or double click the bottle in solution browse screen to start the solution making view as below:

**Make Solution**

Machine No.	1	Product Code	
Bottle No.		Product Name	the material is designed by logcart aux dye machine
Total	800 g	Specific Gravity	1.1
Concentration		Life Time	48 hr
		Dilute From	

Please Input Bottle number

0.00  
TARE

A 2 4 6 1 3 5 1 3 5 B

- “Machine”:** For Selecting machine number.
- “Bottle Number”:** This shows the currently selected bottle position. To change, type in the bottle number
- “Target”:** Show the target weight. For TF system, it should be under 850cc.
- “Product Code”:** Show the selected bottle product code.
- “Concentration”:** The target solution concentration of the make.
- “Specific Gravity”:** Show the specific gravity of the selected solution.
- “Life Time”:** Show the life time of the current selected solution.
- “Dilute from”:** If there is one or more same solution with higher concentration, the filed will be enabled for selecting a bottle of high concentration solution for making solution by diluting from the high concentration bottle.

Please enter Bottle number


- “Message bar”** Show the message of current step.



**“Process Detail”:** Show the detail of the current solution making process. If the process has been done, the check box will be marked.



**“The Agitation bar”** Show the current status of solution mixing.

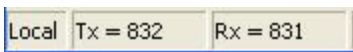

**“The status bar”** Show the online current status PC and system.  
**Procedures**

- Step 1: Select the bottle to make up
- Step 2: Specify whether to make from pure dyestuff, or dilute from a stronger solution.
- Step 3: Put the correct bottle on the scale with a stirrer bar already inside it.
- Step 4: Follow the steps of the process for that solution, which may be a combination of the following:
  - Add Dyestuff manually
  - Add Chemical manually
  - Dispense Cold/Warm/Hot Water
  - Dispense Chemical Automatically (Optional specification)
  - Mix
- Automatically Make up solution to required concentration.



#### 4-2-2 Autolab SPS Program – Dispense – Manual dispense water

---

You can choose the option “Immediately dispense water” in Dispense function of menu bar or double click the bottle in solution browse screen to start the solution making view as below:

Immediately Dispense Water

Target Weight 100 g

Target Temp. 30 °C

Start

Stop

0.00

TARE

A 2 4 6 2 4 6 B

1 3 5 1 3 5

This option can be used to dispense water, which may be useful for various operations within the Laboratory. To do this, select the option, and then specify the Temperature and Target Weight required, and then click “Start” to dispense.

## 4-2-3 Autolab SPS Program – Dispense – Solution List

You can choose the option “All bottle list” in Dispense function of menu bar.

Move cursor onto the bottle, the detail of solution in the bottle will be showed on the left.

### Solution Pre-caution:

**“Low Stock Pre-Caution Weight”:** Set the warning quantity of solution.

**“Life Time Pre-Caution Hour”:** Set the warning life time of solution.

## 4-3 Autolab SPS Program – Database

### 4-3-1 Autolab SPS Program – Database – Solution Data

You can choose the option “Solution Data” in Database function of menu bar.

The solution here is only viewable. To modify the data, operators have to use DP software to modify.

This is for defining solution data.

- “**Machine**”:
  - “**Bottle No.**”:
  - “**Product Code**”:
  - “**Product Name**”:
  - “**Concentration**”:
  - “**Unit**”:
  - “**Dilute From**”:
  - “**Process Code**”:
  - “**Specific Gravity**”:
  - “**Amount to Make**”:
  - “**Use Hot Water**”:
  - “**Life Time**”:
  - “**Print**”:
- For Selecting machine number.
- For key in bottle number.
- For show the product code, operator can change the data also.
- For show the product code.
- For show the concentration of the solution.
- For show the unit of the solution.
- For show the default bottle from diluting.
- For show the process code.
- For show the specific gravity of the solution.
- For show the default capacity of solution bottle, generally, it is 850cc.
- If value is 1, it will open hot water pipe while fill water up.
- For defining the life time of the product.
- Click the button for printing the solution data.

## TF G2 Operation Manual

The data about how and when the solution was made.

Raw Weight	0.40	<b>g</b>
Dilute Source Solution Wt.	---	<b>g</b>
Life Time Remaining	48.0	<b>hr</b>
<b>Made :</b>	26/7/2004 21:35:04	
<b>Use :</b>	Raw Dye	

The solution making process detail


Process Steps
Add Dye
Add Hot Water 30%
Mix Time 1 Min.
Add Chemical 10.0000%
Mix Time 2 Min.

### 4-3-2 Autolab SPS Program – Database – Process Data

---

You can choose the option “Process Data” in Database function of menu bar.

The solution here is only viewable. To modify the data, operators have to use DP software to modify.

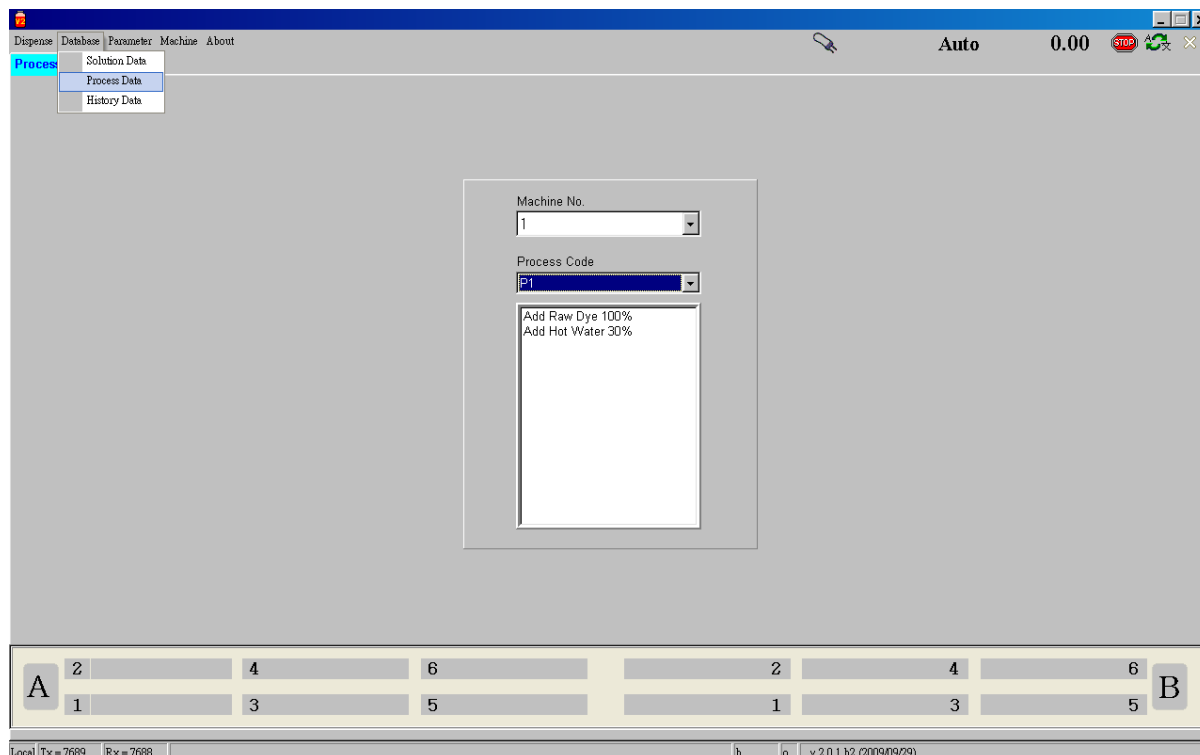

**“Process Code”:** For selecting a process code.

### 4-3-3 Autolab SPS Program – Database – History Data

You can choose the option “History Data” in Database function of menu bar

This function shows the solution making history.


The screenshot shows the 'History Data' window of the TF G2 Autolab SPS Program. The window has a menu bar with 'Dispense', 'Database', 'Parameter', 'Machine', and 'About'. The 'Database' menu is open, showing 'Solution Data', 'Process Data', and 'History Data'. The 'History Data' option is selected. The window displays a table with columns: Time, Machine, Bottle No, From, Product code, Process Code, Amount to make, Concentration, Actual Conc., Unit, and Accuracy. The table is currently empty. At the bottom of the window, there are two rows of buttons labeled 'A' and 'B'. The 'A' row has buttons for 2, 4, 6, 2, 4, 6. The 'B' row has buttons for 1, 3, 5, 1, 3, 5. The status bar at the bottom shows 'Local Tx = 8224', 'Rx = 8223', and 'v 2.0.1 b2 (2009/09/29)'.

Time	Machine	Bottle No	From	Product code	Process Code	Amount to make	Concentration	Actual Conc.	Unit	Accuracy
------	---------	-----------	------	--------------	--------------	----------------	---------------	--------------	------	----------


Local Tx = 8224 Rx = 8223 v 2.0.1 b2 (2009/09/29)

## 4-4 Autolab SPS Program – Parameter Setup

You can choose the option “Setup” in Parameter function of menu bar.

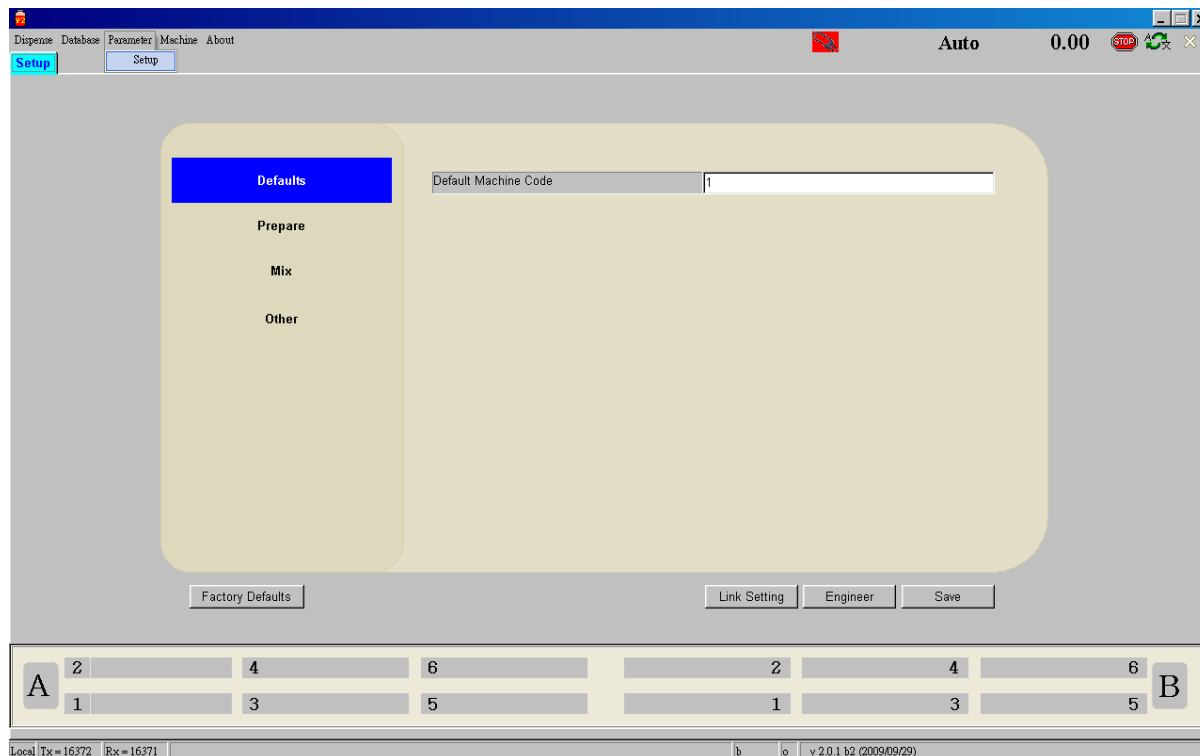

### Default:

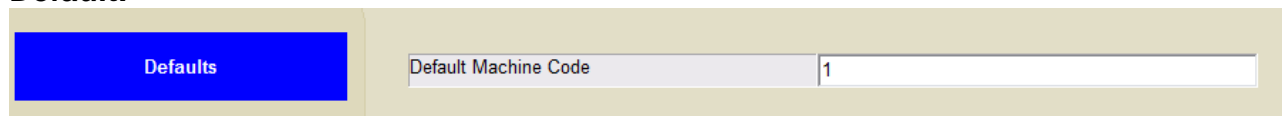

**“Default Machine Code”:** Machine number of dispenser the SPS is linked.

**Prepare:**

Defaults	Hot Water Temperature	95	°C
Prepare	Cold Water Temperature	25	°C
	Water Dispense Tolerance	0.20	g
	Min. Empty Bottle Weight	200	g
	Max. Empty Bottle Weight	500	g
	Add Dye Tolerance (+)	5	%
	Add Dye Tolerance (-)	10	%
	Manual Add Chem. Tolerance (+/-)	5	%
	Min. allowed amount of Dye to Add	0.10	g
	Press CONFIRM Between Steps	Yes	
	Allow change of amount to make	Yes	
	Enable Sound Function	No	
	Mix		
Other			

- “Hot water Temperature”:** The temperature of hot water for this SPS system.  
**“Cold Water Temperature”:** The temperature of cold water.  
**“Water Dispense Tolerance”:** Dispensing tolerance.  
**“Min. Empty Bottle Weight”:** Minimum empty solution bottle weight.  
**“Max. Empty Bottle Weight”:** Maximum empty solution bottle weight.  
**“Add Dye + Tolerance”:** The plus range of manual add dyestuff.  
**“Add Dye – Tolerance”:** The minus range of manual add dyestuff.  
**“Manual Add Chem. Tolerance”:** The tolerance of adding Aux.  
**“Min. Allowed amount of dye to add”:** The minimum weight of dyestuff to be added for making solution.  
**“Press CONFIRM between steps”:** When making solution, enable this function press Confirm button on control panel will be necessary.  
**“Allow change of amount to make”:** Allow to change the quantity of solution in solution making screen.  
**“Enable sound function”:** Use sound to assist adding dyestuff.

**Mix:**

Defaults	Delta Tolerance after Mix (+/-)	1.00	g
Prepare	Confirm Mix Result	No	
	Auto Select Mix Position	Yes	
	Secondary Mix Enable	No	
Mix			

- “Delta Tolerance after mix”:** The tolerance of weight loose after mixing.  
**“Confirm Mix Result”:** Confirm the weight after mixing.  
**“Auto Select Mix Station”:** Allow system select mixer position automatically.  
**“Secondary Mix Enable”:** Allow use second mixer if there are 2 mixers.



**Other:**

<p>Defaults</p> <p>Prepare</p> <p>Mix</p> <p><b>Other</b></p>	Sound Wave Path	c:\ALAB\ALSPS\Wav\	...
	Scale Type	METTLER ▼	
	User Password		

**“Sound Wave Path”:**

The sound files folder.

**“Scale type”:**

Select the scale in the SPS system.

**“User password”:**

User password.

## “Link Setting”

Link SPS Solution database to DP solution database

Machine/Database Link Test Connect

	Link Setting	Server Name	Authentication Mode	Login Name	Password	Server Link Status	Database Link Status
1	YES	.\SQLEXPRESS	Windows Authentication	-----	-----		
2	NO		Windows Authentication				
3	NO		Windows Authentication	-----	-----		
4	NO		Windows Authentication	-----	-----		
5	NO		Windows Authentication	-----	-----		
6	NO		Windows Authentication	-----	-----		
7	NO		Windows Authentication	-----	-----		
8	NO		Windows Authentication	-----	-----		
9	NO		Windows Authentication	-----	-----		

Toggle Link
Server Name
Authen. Mode
User Name
Password
Cancel
Save

This function is to setup the database linking with DP software. The detail of linking setup can refer to AutolabTF S/W Install manual.

Only the system linked with the database, the software can making solution.

## “Engineer”

Click the “Engineer” button to get into the engineer set up :

**Engineer Setup**

**RS232 Setting**

Com Port: 5  
 Baud Rate: 57600  
 Timeout (ms): 1000

**Water Tank Refill Control Parameters**

☒ Enable Refill Control

Enable/Disable    Timeout to alert (ms)

☒ Cold Water    120000  
☒ Hot Water    120000

**Scale Parameters**

Scale Digits: 2  
☒ Define Tare Command: Z  
 Scale Stable Time (ms): 800

**Other Scale Type Setting**

☐ Scale Cmd Pre Code: 0  
 Scale Str Ignore Bytes: 3  
 Scale Start Command: SIR

**Valve Parameters**

Water Sample Time(ms): 10000

Use Valve	Sample Wt.	Valve Response (ms)
<input checked="" type="checkbox"/> Large Cold	55.25	10
<input checked="" type="checkbox"/> Large Cold	41.00	10
<input checked="" type="checkbox"/> Small Cold	14.97	20
<input checked="" type="checkbox"/> Large Hot	61.58	10

☒ Demo Mode Available

Save    Cancel

Normally, for running the software at customer site, all the settings should same as above. And the only setting you need to check is RS232 setting.

### RS232 Setting:

- **COM Port:** It is the PC COM Port the system is connected.
- **Baud rate:** It must be 57600
- **Timeout (ms):** It must be 1000

### “Demo Mode Available”


If you want to run the program in demo mode (without system connected), tick the option “Demo Mode Available”, and restart the program.

## 4-5 Autolab SPS Program – Machine

### 4-5-1 Autolab SPS Program – Machine – Diagnostic

You can choose the option “Diagnostic” in Machine function of menu bar.

The default password for customer configuration the software is “dci”, and there is also a higher priority password for Datacolor engineer use only.

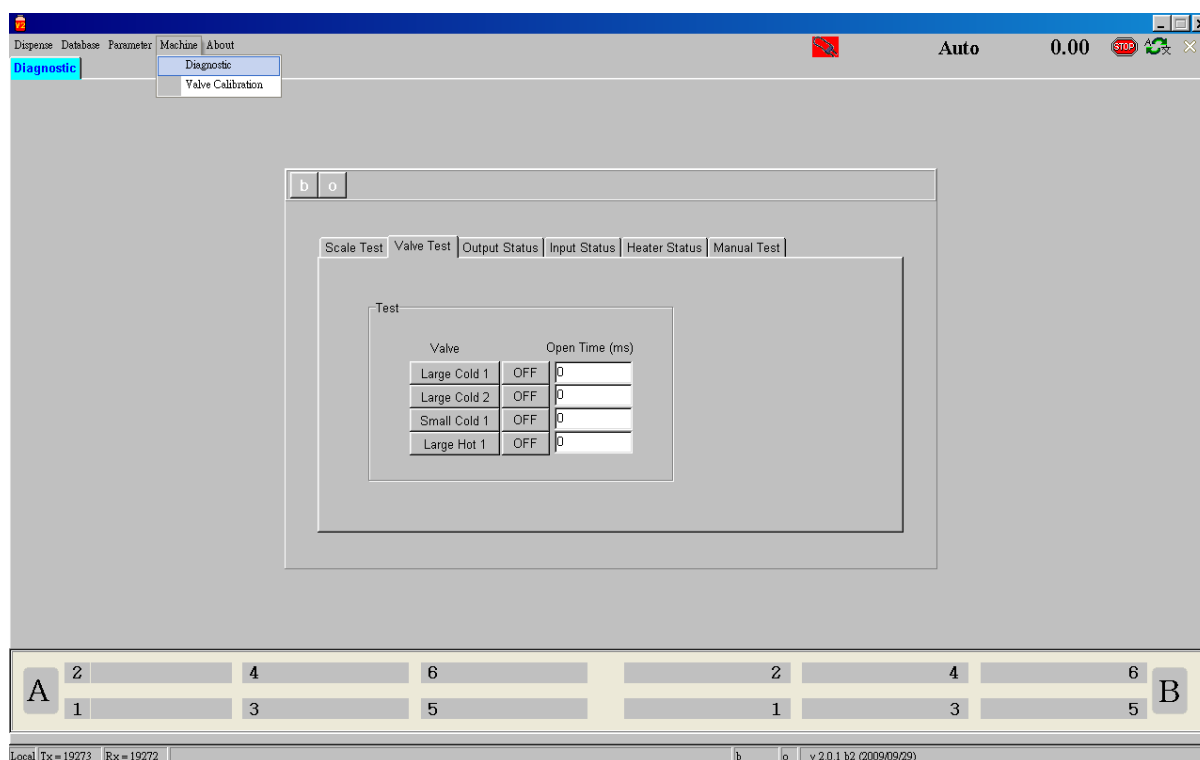

First of all, click on the buttons “b” and “o” to enable the connection between PC and system.

## Autolab SPS Diagnostic - Scale Test

Scale Test | Valve Test | Output Status | Input Status | Heater Status | Manual Test

Scale

Scale String : NSS0.00

Scale Weight : 0.00

Scale Command

This is for testing the scale. Key in “Z” or “SIR” inside the input box, click on the “Balance Command” to see if the scale is reset to zero.

## Autolab SPS Diagnostic - Valve Test

Scale Test	Valve Test	Output Status	Input Status	Heater Status	Manual Test															
<div>Test</div> <table> <thead> <tr> <th>Valve</th> <th></th> <th>Open Time (ms)</th> </tr> </thead> <tbody> <tr> <td>Large Cold 1</td> <td>OFF</td> <td><input type="text" value="0"/></td> </tr> <tr> <td>Large Cold 2</td> <td>OFF</td> <td><input type="text" value="0"/></td> </tr> <tr> <td>Small Cold 1</td> <td>OFF</td> <td><input type="text" value="0"/></td> </tr> <tr> <td>Large Hot 1</td> <td>OFF</td> <td><input type="text" value="0"/></td> </tr> </tbody> </table>						Valve		Open Time (ms)	Large Cold 1	OFF	<input type="text" value="0"/>	Large Cold 2	OFF	<input type="text" value="0"/>	Small Cold 1	OFF	<input type="text" value="0"/>	Large Hot 1	OFF	<input type="text" value="0"/>
Valve		Open Time (ms)																		
Large Cold 1	OFF	<input type="text" value="0"/>																		
Large Cold 2	OFF	<input type="text" value="0"/>																		
Small Cold 1	OFF	<input type="text" value="0"/>																		
Large Hot 1	OFF	<input type="text" value="0"/>																		

Click the button “Calibrate” to calibrate the valves. You need to do it for new installation. Key in the time interval in the “Open Time” input boxes, click on the buttons “Large Cold 1”, “Large Cold 2”, “Small Cold 1” and “Large Hot 1”, the valve should open as the time defined.

## Autolab SPS Diagnostic - Output Status

Scale Test	Valve Test	Output Status	Input Status	Heater Status	Manual Test
<input type="checkbox"/> 00	Large cold 1	<input type="checkbox"/> 08	ACK light		
<input type="checkbox"/> 01	Large cold 2	<input type="checkbox"/> 09	Auto/Manual light		
<input type="checkbox"/> 02	Small cold 1	<input type="checkbox"/> 10	Reserved		
<input type="checkbox"/> 03	Large hot 1	<input type="checkbox"/> 11	Reserved		
<input type="checkbox"/> 04	Cold refill	<input type="checkbox"/> 12	Reserved		
<input type="checkbox"/> 05	Hot refill	<input type="checkbox"/> 13	Reserved		
<input type="checkbox"/> 06	Heater	<input type="checkbox"/> 14	Reserved		
<input type="checkbox"/> 07	Alarm	<input type="checkbox"/> 15	Reserved		

This is just for showing the status of all the OUTPUT.

## Autolab SPS Diagnostic - Input Status

Scale Test	Valve Test	Output Status	Input Status	Heater Status	Manual Test
<input checked="" type="checkbox"/> 00	Auto/Manual input		<input type="checkbox"/> 08	Hot water level sensing	
<input type="checkbox"/> 01	Large cold manual disp		<input type="checkbox"/> 09	Heater enable	
<input type="checkbox"/> 02	Small cold manual disp		<input type="checkbox"/> 10	Overheat	
<input type="checkbox"/> 03	Large hot manual disp		<input type="checkbox"/> 11	Heater Out	
<input type="checkbox"/> 04	Cold water refill		<input type="checkbox"/> 12	Reserved	
<input type="checkbox"/> 05	Ack button		<input type="checkbox"/> 13	Reserved	
<input type="checkbox"/> 06	Hot water refill		<input type="checkbox"/> 14	Reserved	
<input type="checkbox"/> 07	Cold water level sensing		<input type="checkbox"/> 15	Reserved	

This is just for showing the status of all the INPUT.



## Autolab SPS Diagnostic - Heater Status

Scale Test	Valve Test	Output Status	Input Status	Heater Status	Manual Test
------------	------------	---------------	--------------	---------------	-------------

Check Points

Hot Water Level : High Level

Heater Switch : **Off**

Heater Alarm : Off

Heater Out : Off

Heating In Progress : No

This is just for showing the status of the Heater.

## Autolab SPS Diagnostic - Manual Test

Scale Test	Valve Test	Output Status	Input Status	Heater Status	Manual Test
------------	------------	---------------	--------------	---------------	-------------

Refill

Cold Refill On

Cold Refill Off

Hot Refill On

Hot Refill Off

Ack Light

Ack Light On

Ack Light Off

Buzzer

Buzzer On

Buzzer Off


- “Cold Feeding On”:** Click on the button, for opening the cold water refill.
- “Cold Feeding Off”:** Click on the button, for closing the cold water refill.
- “Hot Feeding On”:** Click on the button, for opening the Hot water refill.
- “Hot Feeding Off”:** Click on the button, for closing the Hot water refill.
- 
- “Ack Light On”:** Click on the button, for switching on the Ack light.
- “Ack Light Off”:** Click on the button, for switching off the Ack light.
- “Buzzer On”:** Click on the button, for switching on the buzzer.
- “Buzzer Off”:** Click on the button, for switching off the buzzer.

## 4-5-2 Autolab SPS Program – Machine – Valve calibration

---

You can choose the option “Valve Calibration” in Machine function of menu bar.

The function is for calibrating valve.

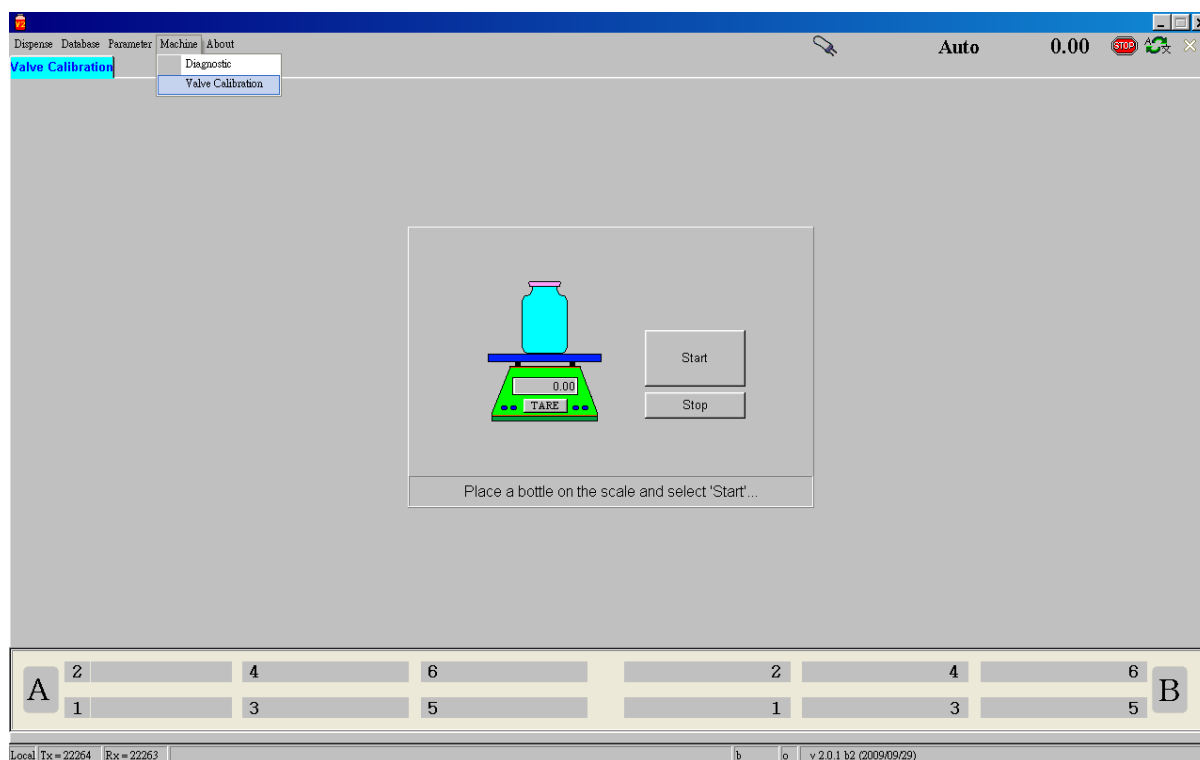

### Trial: Low Liquor Ratio (1:4) Dyeing in Ahiba IR

Machine: Ahiba IR Pro  
Beaker: 300cc beaker with Dosing Lids  
Material: 5 gm Cotton Single Jersey knitted fabric  
Dyes: Reactive Dyes  
Chemicals: Glauber's Salt, Soda Ash, Caustic Soda.  
(Caustic Soda if required as per Shade was taken in liquid form in dosing lid.)

#### Process:

Three recipes were selected (Red, Green and Blue). Recipes were dispensed using Dispenser with maintaining liquor ratio 1:4. As Caustic was taken in liquid form, the quantity to be added later was reduced from total dispensing volume. Salt in solid form was added to the beaker. 5gm fabric was wetted with water and squeezed to maximum using padder.

Fabric was added to the beaker. Soda Ash & Caustic Soda were kept in solid form in dosing lids and loaded to machine.

Machine RPM keeping at 45, temperature was raised to 60°C at gradient of 1.5 °C/min. The machine was run with dyes and salt at this temperature for 45 mins. Temperature was cooled to 45 °C and hold for dosing.

Dosing was done through dosing lids. Both Soda Ash and Caustic were added together. After Dosing Complete, Temperature was raised to 60 °C at 1.5 °C/min. Temperature was maintained at 60 °C and machine was run for 60 mins. After this the program ends. Samples were done a cold wash out side. Neutralization with acetic acid. Then washed with soap at 80°C for 10 mins.

#### Observations:

1. As we could simulate bulk conditions for low liquor, samples were compared with bulk dyed fabrics and results were acceptable.
2. Dosing time was reduced to large extent. This also help as temperature does not falls during dosing time.
3. Found that dosing in liquid form is better than solid form as some particle may remain undissolved during dosing that may affect shade. Particularly in light shades.